DECLARACIÓN DE CONFORMIDAD CE
CE DECLARATION OF CONFORMITY
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita magnetic stirrer with heating series 691. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita magnetic stirrer with heating series 691, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the equipment avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the equipment to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the magnetic stirrer.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to info@auxilab.es or by telephone: +34 807 117 040 (1.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

 **ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

INDEX OF LANGUAGES

Spanish 2-8
 English 9-15
 French..... 16-23

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT 10
 2. DESCRIPTION 10
 3. TECHNICAL SPECIFICATIONS 11
 4. INSTALLATION / SETTING UP 12
 5. MAINTENANCE AND CLEANING 15
 ANNEX I: CE CERTIFICATE 24

1. USES OF THE INSTRUMENT

Nahita magnetic stirrers with heating adapt perfectly to daily lab work allowing to obtain homogeneous solutions and speeding up assays. Thus, they result appropriated to any type of analysis in biology, chemistry or medicine.

2. DESCRIPTION

1.1 Heating plate	1.7 Button Set/Meas: visualization of selected or real temperature
1.2 Display	1.8 Stirring function indicator
1.3 Speed adjustment knob	1.9 Heating function indicator
1.4 Temperature adjustment knob	1.10 Bosshead (only models 691/3 and 691/4)
1.5 Button Temp/Speed: visualization of temperature or speed	1.11 Rod (only models 691/3 and 691/4)
1.6 Button Plate/Sensor (only model 691/4): selection of temperature control in plate or in solution	1.12 Temperature probe (only models 691/3 and 691/4)

5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of the magnetic stirrer it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The magnetic stirrer is supplied with a standard Schuko wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Change of fuse: unplugged the stirrer. The fuse compartment (2.2) is located at the base of the socket. With the help of a sharp object remove the fuse compartment and change the fuse.
- ◆ Never try to repair the magnetic stirrer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the magnetic stirrer (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the magnetic stirrer and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

model) will be displayed.

Note: Due to the resistance of material to thermal transmission, the volume to be heated and the heat lost due to radiation, the temperature inside the recipient will always be lower than the temperature of the plate.

ATTENTION!

There is danger of burns when the heating plate operates over 50 °C. Do not touch the heating plate, the beaker or the liquid.

For safety reasons, when heating flammable samples the selected plate temperature must be always lower than the ignition point of the sample.


Accessories

Other magnets: cylindrical with round ends, oval, orthogonal. Please, visit our web page: www.auxilab.com

Security


- ◆ The magnetic stirrer must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the magnetic stirrer on top of a horizontal, plane and stable table making a free space at least at 30 cm per side.
- ◆ When working at temperatures over 50 °C, do not touch the heating plate, the recipient or the liquid, since you could suffer burns.
- ◆ Maximize precautions when working with liquids that reach temperatures over 100 °C, such as oil.
- ◆ Do not put the magnetic stirrer near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ For safety reasons, when heating flammable samples the selected plate temperature must be always lower than the ignition point of the sample.
- ◆ When you are not using the magnetic stirrer for a long period of time please make sure it is unplugged in order to avoid possible accidents. It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the magnetic stirrer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the equipment (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

Code 50691040. Model 691/4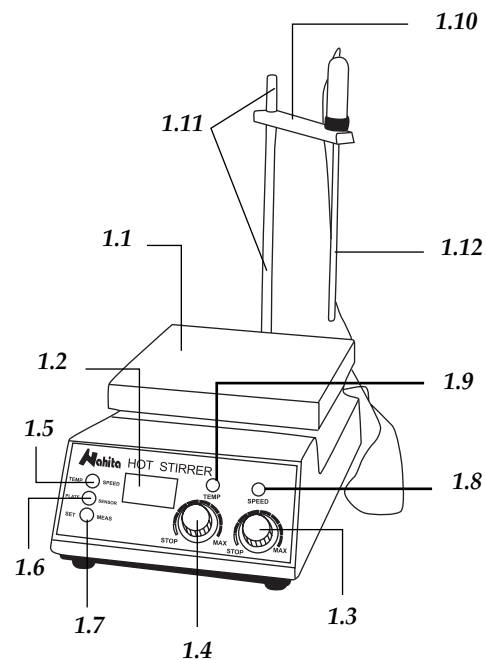

Parte trasera

- 2.1 Socket
- 2.2 Fuse
- 2.3 Rod stand (only models 691/3 and 691/4)
- 2.4 Plug for temperature probe (only models 691/3 and 691/4)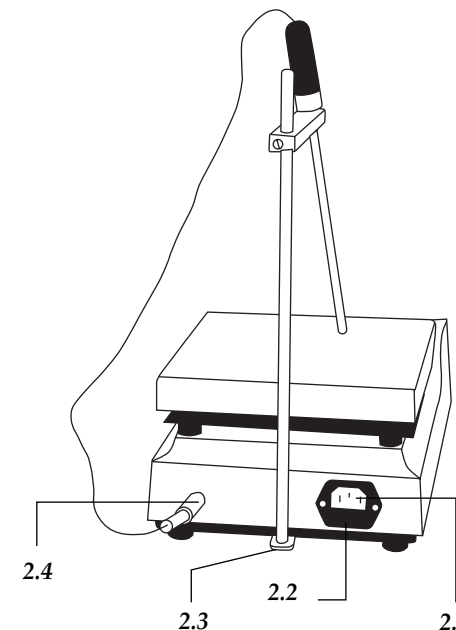

3. TECHNICAL SPECIFICATIONS

Code	50691020	50691030	50691040
Model	691/2	691/3	691/4
Speed range	100-2000 rpm		
Max. temperature	325 °C (placa)	150 °C (sol.)	325 °C (placa) 150 °C (sol.)
Temperature control	Plate	Solution	Solution and plate
Max. volume	5000 mL		
Plate dimensions	175 x 175 mm		
Power	220 V/ 50 Hz		
Dimensions (LxWxH)	300x190x110 mm		
Weight	4 Kg		

4. INSTALLATION / SETTING UP

Preliminary inspection

Unwrap the magnetic stirrer, take off the involving plastic and take off the polyspan protection in which it comes fitted. Take off all protections and without connecting the magnetic stirrer to the power supply, make sure that the equipment does not present any damage because of the shipment. In case it presents any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- Rod (only models 691/3 and 691/4)
- Bosshead (only models 691/3 and 691/4)
- Temperature probe (only models 691/3 and 691/4)
- Teflonated magnet
- Schuko wire
- User's manual
- Warranty certificate

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

Before using the magnetic stirrer, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

- ◆ Put the magnetic stirrer on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put the magnetic stirrer near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
- ◆ During operation, dangerous material such as flammable or infectious liquids must be out of the working area.
- ◆ Installation of the temperature probe (only model 691/3 and 691/4).
 - Screw the rod (1.11) in the stand (2.3) located at the back side of the equipment.
 - Insert the bosshead (1.10) in the rod and fix it at the desired height.
 - Connect the temperature probe (1.12) in its corresponding plug (2.4) and place it into the bosshead.
- ◆ The magnetic stirrer is supplied with a Standard Schuko wire. Make sure that the speed (1.3) and temperature (1.4) adjustment knobs are off and insert the Schuko wire that feeds the AC electric current in the base of current 220 V, 50 Hz \pm 10% provided with earth wire

and to the other end to the stirrer socket (2.1).

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz \pm 10%.

If you are not using the stirrer for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

Setting up

◆ Put the recipient (beaker, Erlenmeyer flask, etc.) with the solution to be stirred and the magnet, in the middle of the plate (1.1). Then, adjust the stirring speed and temperature.

Speed adjustment

Note: Stirring speed must be increased gradually to avoid the magnet spinning asynchronously; if this happens, decrease speed or turn the stirrer off if necessary, turn it on again and increase speed slowly.

- ◆ Put the Temp/Speed button (1.5) in Speed position (outwards).
- ◆ Put the Set/Meas button (1.7) in Set position (inwards).
- ◆ Turn the temperature adjustment knob (1.4) to the minimum.
- ◆ Turn the speed adjustment knob (1.3) clockwise; the stirring function indicator (1.8) will light up and the current speed will be displayed. Adjust the stirring speed according to your necessities.

Temperature adjustment

The temperature control mode depends on the model of stirrer:

- Model 691/2: temperature control in plate
- Model 691/3: temperature control in solution
- Model 691/4: temperature control in plate or in solution

To adjust the temperature, proceed as follows:

- ◆ Put the Temp/Speed button (1.5) in Temp position (inwards)
- ◆ Put the Set/Meas button (1.7) in Set position (outwards)
- ◆ Only for model 691/4: by using the Plate/Sensor button (1.6) select between temperature control in plate (position Plate; inwards) or in solution (position Sensor; outwards).
- ◆ Turn the temperature adjustment knob (1.4) clockwise and select the desired temperature.
- ◆ Put the Set/Meas button (1.7) in Meas position (outwards); the heating function indicator (1.9) will light up and the real temperature (in plate or in solution depending on the