DECLARACIÓN DE CONFORMIDAD CE  
CE DECLARATION OF CONFORMITY  
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

Declara que el equipo | Declare that the equipment | Declare que l'appareil:

CENTRÍFUGA | CENTRIFUGE | CENTRIFUGEUSES

Código | Code | Code: 52818000

Modelo | Model | Modèle: 2818

Cumple las siguientes directivas | Meet the following directives | Accomplit les directives suivantes:

73/23/CE   Directiva de seguridad eléctrica Directive for electrical safety Directive de sécurité électrique	89/336/CE   Directiva de Compatibilidad electromagnética (CEM) Directive for electromagnetic compatibility y (EMC) Directive de compatibilité électromagnétique (CEM)
--	---

Cumple las siguientes Normas: | Meet the following Standards | Accomplit les normes suivantes:

EN 61326 | Material eléctrico para medida control y uso en laboratorio  
Requisitos de compatibilidad electromagnética (CEM.)  
Electrical equipment for measurement, control and laboratory use  
EMC requirements.  
Matériel électriques de mesure, de commande et de laboratoire  
Prescriptions relatives à la CEM.

EN 61010-1 | Requisitos de seguridad de equipos eléctricos de medida, control y uso en laboratorio  
Parte 1: Requisitos generales  
Safety requirements for electrical equipments for measurement, control and laboratory use  
Part 1: General requeriments.  
Règles de sécurité pour appareils électriques de mesurage, de régulation et de laboratoire  
Partie 1: Prescriptions générales.

EN 61010-2-020 | Requisitos particulares para centrifugadoras de laboratorio.  
Particular requirements for laboratory centrifuges.  
Prescriptions particulières pour centrifugeuses de laboratoire.

Fdo: Alfonso Ainciburu Sanz  
DIRECTOR | GERENTE

BERIAIN a 18 de MARZO de 2009

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◆ The Nahita refrigerated centrifuge 2818 can reach 18000 rpm. If any tube breaks, the interior of the centrifuge must be cleaned and disinfected with a disinfectant agent and the broken parts must be removed; otherwise different components of the centrifuge will suffer important damages and due to the lack of cleaning mistakes in subsequent analysis will take place.

**¡ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

## 7. TROUBLESHOOTING

ERROR	CAUSE	SOLUTION
E1	Unbalance	Relocate the tubes in the rotor so as they are perfectly balanced
E2	Spin speed control loss	Contact the Nahita Technical Assistance Department
E3	Starting up with lid open	Close the lid properly
E4	Lid opening during operation	Close the lid and restart
E5	Rotor missing or placed wrongly	Put the rotor properly according to point Placing the rotor
E6	FCR or rpm too high for the rotor used	Select the adequate speed depending on rotor used (see Rotor chart)
E7	Impossibility in detecting rotor (magnet loss)	Check the bottom of the rotor. The rotor can not work without magnets

### INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita refrigerated centrifuge 2818. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita refrigerated centrifuge 2818, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the centrifuge avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the centrifuge to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the equipment.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to [asistencia@auxilab.es](mailto:asistencia@auxilab.es) or by telephone: +34 807 117 040 (0.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the centrifuge warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

## INDEX OF LANGUAGES

Spanish .....	2-14
English .....	15 -26
French .....	27-39

## INDEX OF CONTENTS

1. USES OF THE INSTRUMENT .....	16
2. DESCRIPTION .....	16
3. TECHNICAL SPECIFICATIONS .....	18
4. THEORETICAL FOUNDATIONS .....	18
5. INSTALLATION / SETTING UP .....	19
6. MAINTENANCE AND CLEANING .....	25
7. TROUBLESHOOTING .....	27
ANNEX I: CE CERTIFICATE .....	40

## 1. USES OF THE INSTRUMENT

The Nahita refrigerated centrifuge model 2818 has been especially designed for using in Biochemistry, Cellular Biology, Medicine and research laboratories to carry out protocols of techniques such as PCR, RIA, ELISA, separation and purification of nucleic acids, etc.

## 2. DESCRIPTION

- 1.1 Lid
- 1.2 Rotor arm
- 1.3 Control panel
- 1.4 Switch on/off
- 1.5 Socket

## 6. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

*Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.*

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The centrifuge is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the centrifuge (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.
- ◆ Use the appropriate tubes resistant to the R.C.F. used.
- ◆ Balance the centrifuge by using tubes with the same weight, shape and size and put them in geometrically symmetrical positions. If it is necessary, use tubes filled with water. In a centrifuge it is very important the concept of weight so it is advisable to have a balance to compensate the weight of the different tubes.
- ◆ This is absolutely necessary because an upset of the rotor may produce vibrations that debilitate the centrifuge and rotor increasing the rate of broken tubes and producing a worse sedimentation.
- ◆ The centrifuge is equipped with a safety device that, in case of detecting severe vibrations of the rotor, will automatically stop centrifugation.
- ◆ Never force the stoppage of the centrifuge, it is a risk of accident and debilitation and it may produce as well a mixing of the separated parts.
- ◆ After every use, wipe the condensed water inside the centrifuge chamber. The centrifuge is supplied with a tray that must be placed at the end of the pipe where the water flows out. Empty this tray when needed.
- ◆ Change of fuse: unplugged the centrifuge. The fuse compartment is located at the base of the socket. With the help of a sharp object remove the fuse compartment and change the fuse.

### Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the centrifuge and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

**Visualization of errors**


In case an error appears the centrifuge automatically stops. The error number will be shown on display (2.13) (see Error chart).

**Error chart**

Number	Type of error
E1	Unbalance
E2	Spin speed control loss
E3	Starting up with lid open
E4	Lid opening during operation
E5	Rotor missing or placed wrongly
E6	FCR or rpm too high for the rotor used
E7	Impossibility in detecting rotor (magnet loss)


**Security**

- ◆ The centrifuge must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the centrifuge on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. The ventilation grilles must be always completely free of obstructions.
- ◆ Do not put the centrifuge near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ During operation, dangerous material such as flammable liquids or pathological material should be out of the working area.
- ◆ Nahita centrifuges are not suitable for treatment of flammable liquids, explosive samples or samples capable of forming explosive mixtures.
- ◆ The centrifuge is provided with a safety switch that avoids operation if the lid is open and that in case of opening the lid during operation automatically stops centrifugation. It is also provided with an electromagnetic lock that locks the lid when the rotor starts spinning. Never try to open the lid while centrifuge is operating.
- ◆ When you are not using the centrifuge for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the centrifuge (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.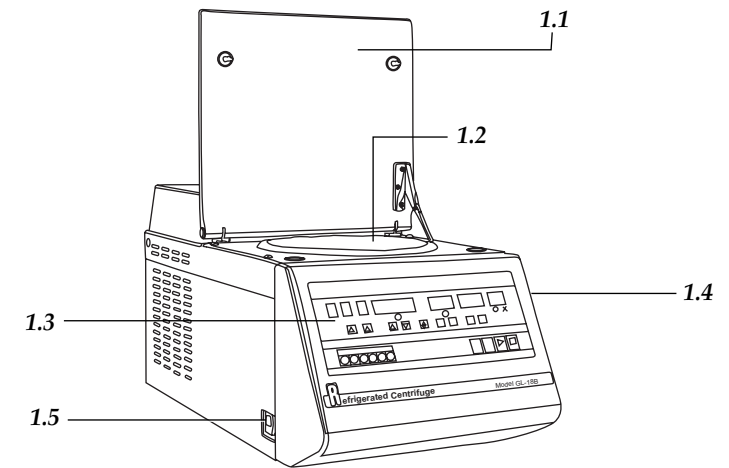

**Panel of controls**

- 2.1 State indicator
- 2.2 Acceleration rate display
- 2.3 Acceleration rate selection button
- 2.4 Braking rate display
- 2.5 Braking rate selection button
- 2.6 Speed display
- 2.7 Speed adjustment buttons
- 2.8 FCR/rpm selection button
- 2.9 Time display
- 2.10 Time adjustment buttons
- 2.11 Temperature display
- 2.12 Temperature adjustment buttons
- 2.13 Rotor/Error display
- 2.14 Program selection buttons
- 2.15 Lid lock indicator
- 2.16 Pre-cooling function button
- 2.17 Quick spin function button
- 2.18 Start
- 2.19 Stop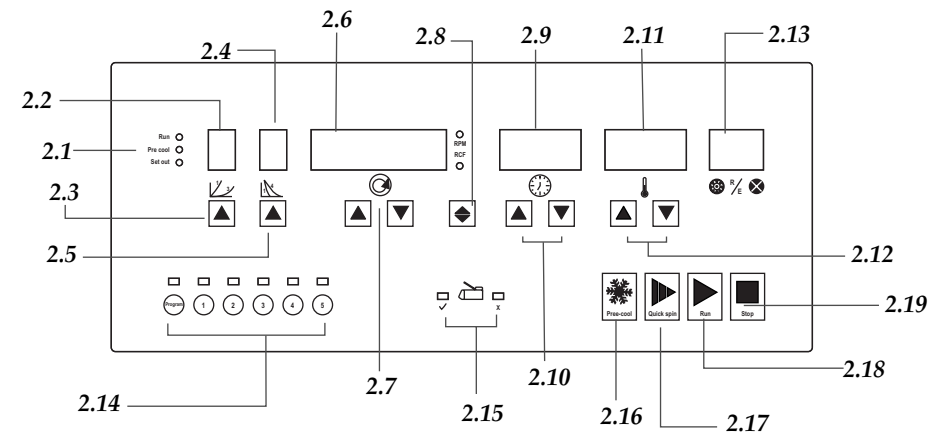

### 3. TECHNICAL SPECIFICATIONS

<b>Model</b>	2818
<b>Rotor</b>	Fixed angle (not included)
<b>Automatic detection of rotor</b>	Yes
<b>Temperature</b>	-10 °C to 40 °C (± 2 °C)
<b>Timer</b>	0-9 h 59 min
<b>Memory</b>	5 programs
<b>Acceleration rate</b>	3 levels
<b>Braking rate</b>	4 levels
<b>Other functions</b>	Quick spin, pre-cooling
<b>Motor</b>	Induction, maintenance-free
<b>Noise</b>	< 68 dB
<b>Power</b>	220-230 V, 50 Hz
<b>Dimensions</b>	74x38x38 cm
<b>Weight</b>	60 Kg

#### Maximum speed according to rotor

Code	92818000	92818001	92818002	92818003	92818004
Capacity	12x1.5 mL	24x1.5 mL	12x5 mL	12x10 mL	6x50 mL
Max. speed	18000 rpm	15000 rpm	15000 rpm	12000 rpm	10000 rpm
Max. RCF	21300 g	20600 g	18600 g	13000 g	10500 g

### 4. THEORETICAL FOUNDATIONS

Centrifugation is a separation technique based on the movement of the particles that are displaced to the distal end of the rotation axis depending on their shape and mass.

The Relative Centrifugal Force (R.C.F.) is the required force to separate particles. It is given as some value times that of gravity (g).

The R.C.F. is calculated with the following formula:


$$F.C.R. = 1.118 \cdot 10^{-5} \cdot r \cdot n^2$$

$1.118 \cdot 10^{-5}$  is a constant.

$r$  = radius, the horizontal distance in centimetres

from the rotation axis to the bottom of the sample tube.

$n$  = rotation speed expressed in revolutions per minute (r.p.m.).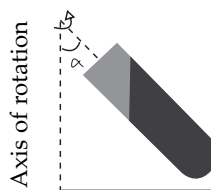

Radio of draft

#### Pre-cooling function

This function allows the centrifuge chamber to reach quickly the pre-selected temperature. For this purpose:

- Once temperature has been set press the pre-cooling function button (2.16).
- The rotor will spin at 3000 rpm and the timer will start counting up. The temperature inside the centrifuge will go down quickly until reaching the pre-selected temperature; once it has been reached, the rotor will go on spinning to keep it.

**Note:** To allow the rotor itself reach the pre-selected temperature pre-cooling time must be longer. The larger the rotor is, the longer the time must be.

**Note:** if temperature decreases slowly, check that the ventilation grilles of the equipment are completely free of obstructions

- ◆ To start with the centrifugation protocol, press stop (2.19) and then start (2.18).

#### Programs

The centrifuge 2818 presents a memory to store up to 5 different centrifugation programs. To save a program, proceed as follows:

- ◆ Select the parameters of the protocol to be saved.
- ◆ Press Prog. button of the program selection buttons (2.14)
- ◆ Then, press the number in which the program is going to be saved.
- ◆ To select a program preciously saved, press the corresponding button number and the different parameters will be displayed.

#### Rotor detection

Once it is operating, the centrifuge detects the rotor used and shows its corresponding number (Rn) on display (2.13). If pre-selected speed is not adequate for the rotor used, the centrifuge will stop automatically and will display the error.

#### Rotor chart

Number	Descripción	Capacity	Max. speed (rpm)	Max. RCF (g)
R0	Rotor waiting to be detected			
R1	Fixed angle rotor 45°	12x1.5 mL	18000	21300
R2	Fixed angle rotor 45°	24x1.5 mL	15000	20600
R3	Fixed angle rotor 32°	6x50 mL	10000	10500
R4	Fixed angle rotor 45°	12x5 mL	15000	18600
R5	Fixed angle rotor 30°	12x10 mL	12000	13000


Once the different parameters have been set (acceleration and braking rates, speed, temperature and time), proceed as follows:

- ◆ Put the tubes in the rotor making sure they are perfectly balanced.

*Note: an upset of the rotor caused by tubes not well balanced will produce vibrations that debilitate the centrifuge. The equipment presents a safety device that automatically stops centrifugation in case of detecting severe vibrations of the rotor.*

- ◆ Put the lid of the rotor and hold it well with the screw at the top (Fig. 3).
- ◆ Close the lid of the centrifuge.
- ◆ Press start (2.18) to begin operation. Automatically, the lid will be locked by means of an electromagnetic lock and the lid indicator (2.15) will show that the lid is blocked (orange light).
- ◆ The centrifuge will stop after the selected time has passed. When the rotor stops, the lid will be automatically unlocked and the lid indicator (2.15) will show that the lid can be open (green light). To stop the centrifuge before the operation time has passed, press stop (2.19).

*Note: If power cut occurs during operation, the lid must be unlocked manually. For this, introduce the needle supplied into the hole placed at the right side of the centrifuge (Fig. 4).*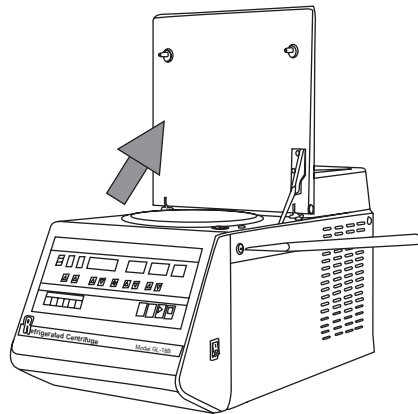

### Other functions

#### Quick spin function

To carry out quick spins, keep pressed the quick spin function button (2.17). The rotor will spin at 8000 rpm until loosing the button.

## 5. INSTALLATION / SETTING UP

### Preliminary inspection

- ◆ Unwrap the centrifuge, take off the involving plastic and take off the polyspan protection in which it comes fitted.
- ◆ Take off all protections and, without connecting the centrifuge to the net, check manually that the rotor spins freely, without any obstacle. Make sure that the equipment does not present any damage because of the shipment. In case it presents any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- Water tray
- Schuko wire
- Spanner
- Needle
- Fuse
- User's manual
- Warranty certificate

*We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.*



### Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

**PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.**

- ◆ Please put the centrifuge on top of a horizontal, plane and stable table making a free space around so as the ventilation grilles are completely free of obstructions.
- ◆ Do not put the equipment near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. During operation, people and dangerous material such as flammable or infectious liquids must be out of the working area.
- ◆ Please insert the Schuko wire that feeds the AC electric current in the base of current 220 V, 50 Hz  $\pm$  10% provided with earth wire and to the other end to the centrifuge socket (1.5).

*Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz  $\pm$  10%.*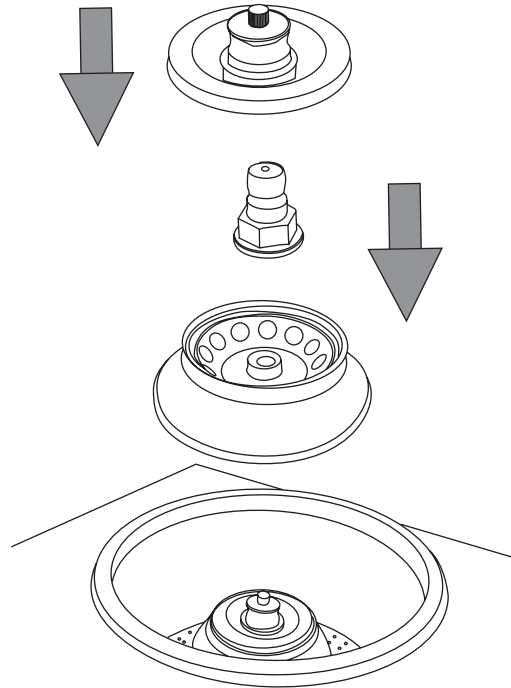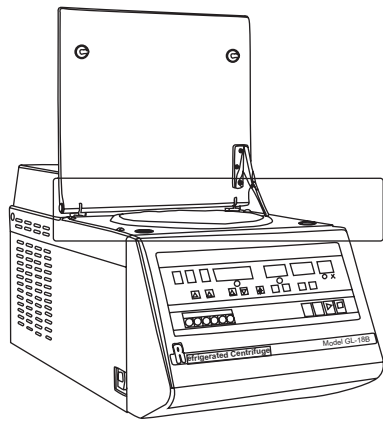

### Placing the rotor

- ◆ Carefully insert the rotor in the rotor arm.
- ◆ By using the spanner, tighten firmly the nut supplied with the rotor.
- ◆ Once tubes have been placed inside the rotor, put the lid so as the point of the rotor and the lid match, and tighten manually the screw on the top of the lid, until it is well fixed.

**Note:** Please be careful to avoid mixing the nuts and lids of the different rotors.

**Note:** The centrifuge automatically detects the installed rotor. Please, check the rotor bottom before installation to make sure the magnets that allow the centrifuge to detect the rotor are in perfect conditions

If you are not using the centrifuge for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

### Setting up

Press the on/off switch (1.4) to turn the equipment on. Then, proceed with the selection of the different parameters of the centrifugation protocol

- This parameter is related with the time that the rotor needs to reach the pre-selected speed. The centrifuge is provided with 3 acceleration rate levels, being the level 1 the fastest and the level 3 the slowest:

- Select the acceleration level by using the corresponding button (2.3)
- The selected level will be shown on display (2.2)

### Braking rate

This parameter is related with the time that the rotor needs to stop when spinning at the selected speed. The centrifuge is provided with 4 braking rate levels, being the level 1 the fastest and the level 4 the slowest:

- Select the braking level by using the corresponding button (2.5)
- The selected level will be shown on display (2.4)

### Speed adjustment

The maximum centrifugation speed depends on the rotor used (see Rotor chart). Once it is operating, the centrifuge automatically detects the rotor, and warns and stops if the selected speed is not adequate for that rotor.

- Select between rpm or FCR by using the corresponding button (2.8).
- Press the speed adjustment buttons (2.7) to select the desired centrifugation speed. Speed adjustment is performed in steps of 100 rpm or 100 g.

### Time adjustment

The centrifuge can operate in periods of time up to 9 h 59 min. For this purpose:

- Select the centrifugation time by using the time adjustment buttons (2.10)
- When rotor spins at the selected speed, the timer starts counting down; the remaining operation time will be shown on display (2.9).

### Temperature adjustment

Nahita refrigerated centrifuge 2818 allows centrifugation of samples at low temperatures down to -10 °C. For this purpose:

- Temperature display (2.11) shows the real temperature inside the centrifuge.
- Press the temperature adjustment buttons (2.12) to visualize the pre-selected temperature and to select the desired temperature. 5 seconds after setting temperature the real temperature inside the centrifuge will be displayed again.

**Note:** Please read the point of Pre-cooling function.