AUXILAB S.L.

Material de laboratorio  
Laboratory supplies

DECLARACIÓN DE CONFORMIDAD CE  
CE DECLARATION OF CONFORMITY  
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

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**ANNEX I**

**Refractive index correction chart with distilled water temperature**

Temperature °C	Refraction Index	n <sub>F</sub> -n <sub>C</sub> Mean dispersion
10	1.33369	0.00600
11	1.33364	0.00600
12	1.33358	0.00599
13	1.33352	0.00599
14	1.33346	0.00599
15	1.33339	0.00599
16	1.33331	0.00598
17	1.33324	0.00598
18	1.33316	0.00598
19	1.33307	0.00597
20	1.33299	0.00597
21	1.33290	0.00597
22	1.33280	0.00597
23	1.33271	0.00596
24	1.33261	0.00596
25	1.33250	0.00596
26	1.33240	0.00596
27	1.33229	0.00595
28	1.33217	0.00595
29	1.33206	0.00594
30	1.33194	0.00594
31	1.33182	0.00594
32	1.33170	0.00593
33	1.33157	0.00593
34	1.33144	0.00593
35	1.33131	0.00592
36	1.33117	0.00592
37	1.33104	0.00591
38	1.33090	0.00591
39	1.33075	0.00591
40	1.33061	0.00590

Thanks for choosing a Abbe Refractometer ZUZI. It is advisable to read carefully this manual and follow the processes of use.  
 ZUZI develops its products putting more emphasis on Ergonomics and Security for users. Its high mechanical and optical quality will allow you to have this equipment available for many years.

**BEFORE CONNECTING OR OPERATING WITH THIS INSTRUMENT, READ THESE INSTRUCTIONS THOROUGHLY.**

- In order to use this equipment, bear in mind the following points:
- ◆ As it is a precision instrument, it should be handled carefully avoiding sudden movements and knocks.
  - ◆ Never dismantle the different pieces of the balance to repair them yourself. In addition to loosing the warranty it could lead to a faulty performance of the equipment.
  - ◆ In order to prevent fire or electrical unloadings, avoid atmospheres excessive humidity.
  - ◆ To prevent fire or shock hazard, do not expose the balance to rain or moisture.
  - ◆ Any doubts can be clarified by contacting the ZUZI Technical Service, by e-mail asistencia@auxilab.es, by telephone to the 948310513 or by fax 948310500 - 948312071.
  - ◆ Do not forget to fill up and send it's certificate of warranty in 15 days. It has a year of warranty.

**ATTENTION!!  
 IF AN EQUIPMENT IS NOT PROPERLY CLEAN AND DISINFECTED IT  
 WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

**INDEX OF LANGUAGES**

Spanish ..... 2-10  
 English..... 11-19  
 French ..... 20-28

**INDEX OF CONTENTS**

1. APLICACIONES OF THE INSTRUMENT ..... 12  
 2. DESCRIPTION ..... 12  
 3. WORKING PRINCIPLES ..... 13  
 4. TECHNICAL SPECIFICATIONS ..... 14  
 5. OPERATING INSTRUCTIONS ..... 14  
 6. MAINTENANCE ..... 17  
 ANNEX I: REFRACTIVE INDICES WITH TEMPERATURE TABLE ..... 18  
 ANNEX II: WARRANTY INFORMATION ..... 19

## 1. APPLICATIONS OF THE INSTRUMENT

The refractometer is used to measure the refractive indice nD and mean dispersions nF- nC of transparent and translucent liquid or solid, mainly the transparent liquid. Zuzi 320 refractometer can be connected to a circulation thermostat. Working temperature range is 0- 100° C. See joint table to refraction index correction with temperature.


Refractive index is an important optical constant in substance and can be used to determine optical performance, purity and concentration, etc.

It is possible to get sugar percentage in saccharose solutions within a range of 0 - 95% Brix, equal to refractive indice 1.333 - 1.531.

This instrument can be widely used as indispensable tool in petrological, oil and fat, pharmaceutical, food, cosmetic or chemical industry, as well as universities and research institutions.

## 2. DESCRIPTION

Picture 1

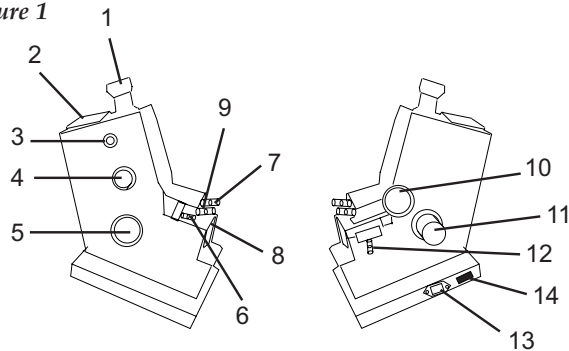- 1.1. Eyepiece
- 1.2. Reading scale
- 1.3. Calibration screw
- 1.4. Scale knob for dispersion graduation
- 1.5. Refraction index adjustment knob
- 1.6. Top prism circuitation thermostat connections
- 1.7. Bottom prism circuitation thermostat connections

- 1.8. Illumination window of solid samples
- 1.9. Illumination window of liquid samples
- 1.10. Close prism control
- 1.11. Lamp
- 1.12. Thermometer socket
- 1.13. Current intake
- 1.14. ON/ OFF switch

## 6. MAINTENANCE

- ◆ To ensure accuracy and to avoid instrument damage, please, pay attention to your refractometer maintenance.
- ◆ Follow the instructions and warnings relating to this manual.
- ◆ You should have this manual always by hand, so that any person can consult it.
- ◆ Preserve the balance from heavy shocks. Look for a dry place away from dust, direct sun beams or air flows. Since it is a precision instrument, it should be handled carefully.
- ◆ For security never disconnect it by pulling the cable.
- ◆ Never remove any of its mechanical components. In addition to losing the warranty, it could also lead to a faulty performance of the balance.
- ◆ It must be kept in a dry and well ventilated room to preserve optical items from humidity and mold.
- ◆ If you deal with corrosive liquids, you must clean all the instruments immediately to prevent them from rust damages you should include in this list optical and mechanical parts and also painted surfaces.
- ◆ After each use, you must clean the refractive prisms with a non fuzz gauze, use alcohol or water.
- ◆ Liquid samples must not contain solid impurities. Solid samples must have polished and flat faces to avoid scratching the prism.
- ◆ When not using the refractometer for long periods, keep it in a box with a draining substance (Silica Gel).
- ◆ Never take pieces apart to try repairing it by yourself because you may not only miss warranty but also make it does not work.
- ◆ If a lamp flux, replace it by an original ZUZI one of 6V 20W. Pull control 3.10 and take out the lamp from the lamp holder and replace it by the new one.
- ◆ In the rear you can find the fuse ("FUZE"). If the lamp does not light, check the fuse is in good condition. If it is fluxed, replace it by a new one.
- ◆ For any inquiry, please contact Zuzi official service.

**a) To test transparent and translucent liquids**

- ◆ Drop the tested liquid onto prism surface using a dropping tube. Cover the incident prism and lock with the knob (1.10). The sample must be homogeneous and spread prism surface wide without bubbles.
- ◆ Open the shield (1.9) and check the reflecting mirror (1.1) is closed, then adjust the ocular (1.1) based on the dioptic correction required until you see the crossed lines sharply.
- ◆ Turning the refraction index adjustment knob (1.5) look for light and dark field divisory line, then turn knob (1.4) to adjust the dispersion in a way this line is seen without any colour and as clear as possible. Once you have done this, turn the knob (1.5) again to move divisory line up to the middle of both crossed lines.
- ◆ The value in the scale tells us the accurate liquid refraction index.

Reading

- ◆ The range of the reading scale is in refractin index from 1.3000 to 1.7000 and from 0 to 95% in Brix degrees.

Example: In this scale we read 1.3355 nD and 12.5° Brix.


**b) To test transparent solids**

- ◆ Solid to be measured must be smooth and polished faced.
- ◆ Put one or two drops of naphthalene monobromure on the refraction prism surface. Clean all the tested object surface properly to get the best contact. Open the illumination window for solid samples (1.8) and read the value on the scale.
- ◆ When observing through the eyepiece, search for the divisory line between light and dark coloured fields, and do as the instructions in the previous case.

**c) To test translucent solid**

- ◆ The object to be measured must be smooth and polished faced.
- ◆ Put one or two drops of naphthalene monobromure on the solid and clean solid surface to place it on the refracting prism. Open the reflecting mirror (1.8) to solid samples illumination.

Picture 5

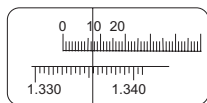**d) To measure sugar concentration in sugary liquids**

- ◆ The procedure is similar to liquid samples but the reading can be done on a Brix scale (larger scale 0- 95%) reading directly the sugar concentration in the liquid.

**e) Refractive index estimation at different temperatures**

- ◆ Insert the thermometer into its socket (1.11), connect the thermostat to the prepared connectors (1.6) for bottom prism and (1.7) for top prism, thanks to the thermostat temperature can be read in the ranges required. Before each reading, wait a couple of minutes to stable the temperature.

**3. WORKING PRINCIPLES**

Refractometry

- ◆ When a light beam arrives to the divisory surface between two different mediums, one part is reflected and the other is bended (it comes into the second medium), being spread with an angle and speed different to first medium. Both values are measured by Refractive index. Each substance has its own refractive index value.
- ◆ Refractive index is related to mass, charge and the number of particles in the substance through which the light beam is transmitted. There are different kinds of refractometers but they all are based on the same principle.
- ◆ The basic refractometer principle is based on the following refraction law:


$$n_1 \cdot x \cdot \sin \alpha' = n_2 \cdot x \cdot \sin \alpha''$$

Being  $n_1$  and  $n_2$  medium refractive index at both limit sides between different sides medium, where  $\alpha'$  is the incident angle and  $\alpha''$  refractive angle.


- ◆ When a light beam enters an optically lighter medium from a thicker one, the incident angle is smaller than refractive angle and if we change the first one until the second one reaches 90°, the incident angle is named critical angle.

- ◆ As shown in picture 3, when a light beam falls on surface AB from any angle, its refractive angle is always bigger than i, if we observe the emergent light beam through a microscope, it appears the view field divided in two halves, one dark and the other one light coloured. If we place the two halves separation line just in the middle, this is the critical angle position.

Picture 3 - Critical angle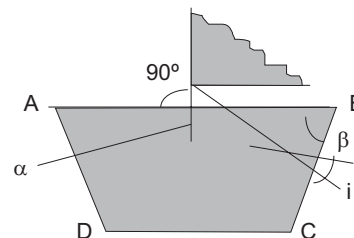

Picture 2

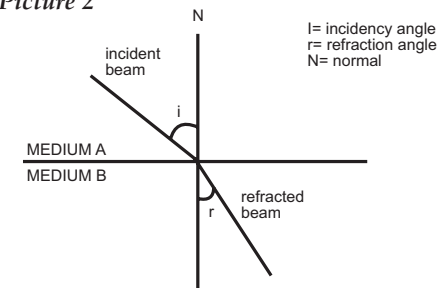- ◆  $ABDC$  is the refracting prism whose refractive index is  $n_2$ , and the sample to be measured (liquid or transparent solid) is placed on  $AB$  surface. Its refractive index is  $n_1$ .

- ◆ From refraction law, we obtain the following equation:

$$\begin{aligned} n_1 \cdot x \cdot \sin 90^\circ &= x \cdot \sin \alpha \\ n_2 \cdot x \cdot \sin \beta &= x \cdot \sin i \\ \phi &= \alpha + \beta \\ \alpha &= \phi - \beta \end{aligned}$$

Having this in the first equation:

$$n_1 = n_2 \times \sin(\alpha - \beta) = (\sin \phi \times \cos \beta - \cos \phi \times \sin \beta)$$

$$n_2^2 \times \sin^2 \beta = \sin^2 i$$

$$n_2^2 \times (1 - \cos^2 \beta) = \sin^2 i$$

$$n_2^2 - n_2^2 \times \cos^2 \beta = \sin^2 i$$

$$\cos \beta = \sqrt{\frac{n_2^2 - \sin^2 i}{n_2^2}}$$

$$n_1 = \sin \phi \times \sqrt{n_2^2 - \sin^2 i} - \cos \phi \times \sin i$$

Prism refractive angle  $\phi$  and prism refractive index  $n_2$ , are calculated when knowing the critical angle  $i$ . Sample refraction index can be known.

#### 4. TECHNICAL SPECIFICATIONS

##### ABBE Refractometer Model 320

<i>Measurement range</i>	
Refraction Index	1'3000 to 1'7000
° Brix	0 to 95%
<i>Measurement accuracy</i>	
Refraction Index	± 0'0005
° Brix	± 0'2
<i>Big sized prism</i>	6 cm <sup>2</sup>
<i>Thermometer range</i>	0 - 100° C
<i>Working temperature</i>	0 - 100° C
<i>External dimensions</i>	120 x 310 x 210 mm
<i>Weight</i>	2.6 Kg
<i>Power supply</i>	220V±10%.50Hz

#### 5. OPERATING INSTRUCTIONS

##### Preliminary examination

- ◆ Remove the Abbe Refractometer from the packing material and examine it carefully to make sure no damage has occurred during shipping. If there is any damage, notify to your dealer/forwarding agency.
- ◆ Keep all packing material until you are sure that the instrument works correctly. All defective items must be returned in the original packing together with any other supplied accessories.

- ◆ Verify you have received all the accessories that come with the Abbe Refractometer:
  - Thermometer up to 100° with steel socket included
  - Known refraction index (nD 1.5168) test part
  - Disolution of naphthalene monobromide
  - Allen Key
  - Instruction manual
  - Warranty

##### Installation

- ◆ Before starting this instrument, read carefully this manual.


**BEFORE CONNECTING OR OPERATING WITH THIS INSTRUMENT, READ THESE INSTRUCTIONS THOROUGHLY.**

- ◆ Place the refractometer in a flat horizontal table, avoiding humidity, heat, dust and vibrations. Keep it, if possible out of magnetic source, electric field or electric machine able to generate high frequency field.
- ◆ These microscopes have a Schuko plug. This plug must be connected to a plugging device connected to ground. The grounded device must be at hand to be able to disconnect the plug quickly in case of emergency.

*Neither the manufacturer nor the supplier will assume any responsibility for the damages caused to the microscope or for any other damage suffered by any person due to the neglect of the right process of electric connection. The alternating voltage must be 220-24 v, 50-60Hz.*

##### Calibration

- ◆ Before using the first time, you must check if the refractometer is well calibrated, that's why it is included piece to test the known refraction index.
- ◆ If you put two drops of naphthalene bromide on refraction prism and you place then the smooth test piece item over them, you can see though the eyefield the line that separates light and dark coloured parts (pic. 4). Once the dispersion is corrected, you agree inferior side scale with the amount in the test piece. If the line that divides light and dark parties do not meet in the center of the crossed lines seen in the top, you must to turn the calibration screw (3.3) until you have them met. This screw is covered by a lid, unbolt it and with the Allen key supplied with the equipment, turn the screw up to the point required.
- ◆ Once the refractometer is calibrated, the screw do not be manipulated while measuring. If you doubt about the refraction index estimation when testing, please, calibrate it again from the beginning.
- ◆ The ridged incident prism face, the smooth refraction prism face and test piece must be clean with the aid of an absorbent cotton gauze wet with a 1:1 absolute alcohol and ether mixture in order to avoid from the dirtiness that have an effect on image brightness and measurement accuracy.

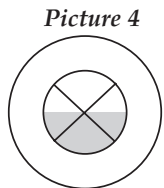##### Testing work

- ◆ Plug the refractometer and turn on the interruptor (1.13) to get the scale display lighted. Do as the instructions say depending on the kind of sample whose refraction index or Brix degrees you want to get.