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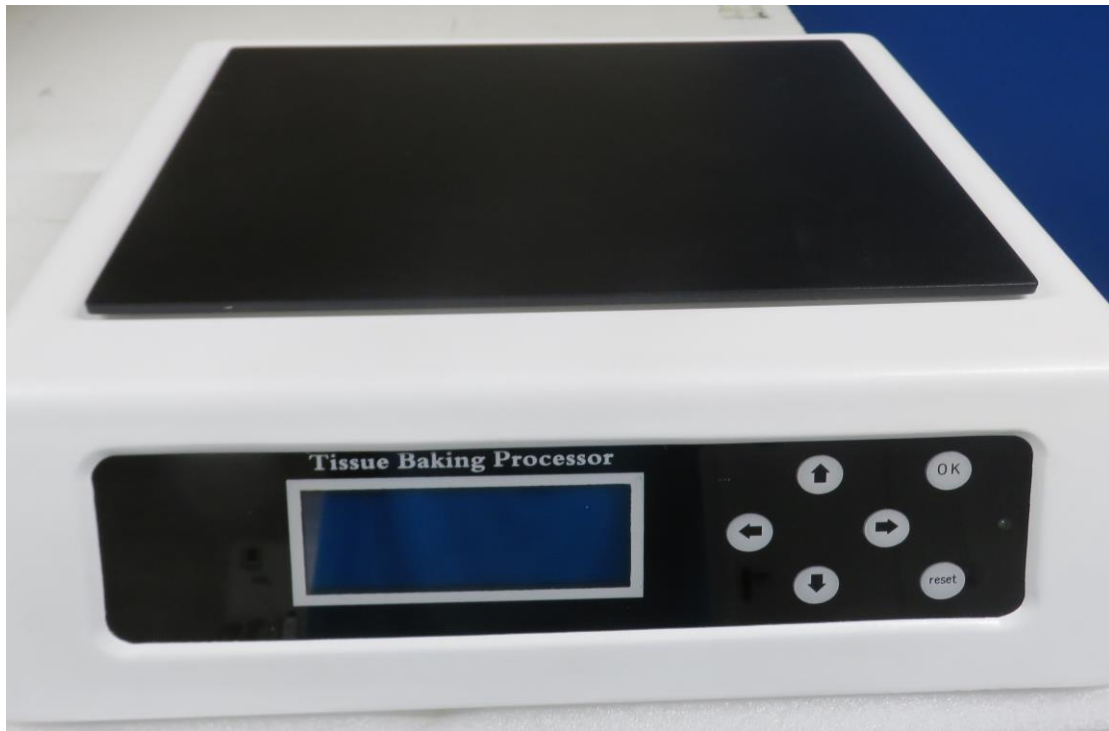
Please read this Manual before using the equipment and keep it for future use.

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Contents

1. Use of the equipment
2. Technical Parameters
3. Installation
4. Operation

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Use of the equipment

This equipment is used to dry the slides used in pathological studies carried out in

hospitals or research laboratories and in teaching in schools and universities.

It has a microprocessor control that allows the temperature to increase quickly and remain stable. The liquid crystal display and the control panel make it easy to configure. Automatically memorizes the last used temperature.

Technical parameters

1. Rated voltage: 220V
2. Rated current: 1.4A
3. Power: 300W
4. Temperature range: RT - 99°C.
5. Heating plate dimensions: 320 mm × 260 mm

Installation

Initial inspection

- Unpack the Tissue Baking Processor, remove the plastic that surrounds it and remove the poliespan protection.
- Without connecting the equipment to the power supply, make sure that it does not present any damage due to transport. If so, notify your distributor immediately so you can make the relevant claims within the deadline set by the transportation service.

Installation

- Place the Tissue Baking Processor on a horizontal and stable table, leaving the necessary work space around it.
- Comply with all the measures required by current legislation on occupational safety and health (electrical safety devices, flammable, toxic and / or pathological

products, etc.).

- Do not place the equipment in areas near heat sources, magnetic material or expose it directly to sunlight.
- The presence of flammable products (gases and liquids) in the workplace is prohibited.
- Insert the power cable into the 220V 50Hz socket provided with a ground connection and at the other end to the connector of the Tissue Baking Processor.
- When you are not going to use the equipment for long periods of time, make sure to disconnect it from the power.

Operation

Frontal panel

- Turn on the equipment by pressing the "ON-OFF" switch on the back.
- Press the OK key. The menu screen will appear with the options TIME, START, TEMPERATURE and AUTOMATIC. You can select one of them using the arrow keys and then pressing OK.
- By pressing the RESET key you can return to the menu screen.
- The Tissue Baking Processor has two modes of operation: normal and automatic.

- Normal operation: Select TEMPER. Press the key →. Adjust the heating temperature value by pressing ↑ or ↓. Then press OK.
- Select START to begin heating the plate. The set temperature (Setup) and the temperature of the plate (Baker) will appear on the screen.
- To end the plate heating, press RESET. Turn off the equipment by pressing the "ON-OFF" switch.
- Automatic operation: Select TIME. This screen shows SYS TIME, START TIME, and EXIT TIME.
- Press → to adjust the SYSTEM TIME. Press ↑ or ↓ to adjust the hours. Press →. Press ↑ or ↓ to adjust the minutes. Press OK.
- Select TIME again. Set the START TIME and the EXIT TIME. Press OK.
- Select AUTOMATIC. The display shows WAITING, the system time and the start time for heating.
- When the start time comes, the screen shows the set temperature (Setup) and the plate temperature (Baker).
- When the exit time comes, the screen shows again WAITING, the system time and start time for heating.
- If you are not going to use the equipment again, turn it off by pressing the "ON-OFF" switch.