
BALANZA DIGITAL NAHITA
NAHITA DIGITAL BALANCE
BALANCE DIGITAL NAHITA

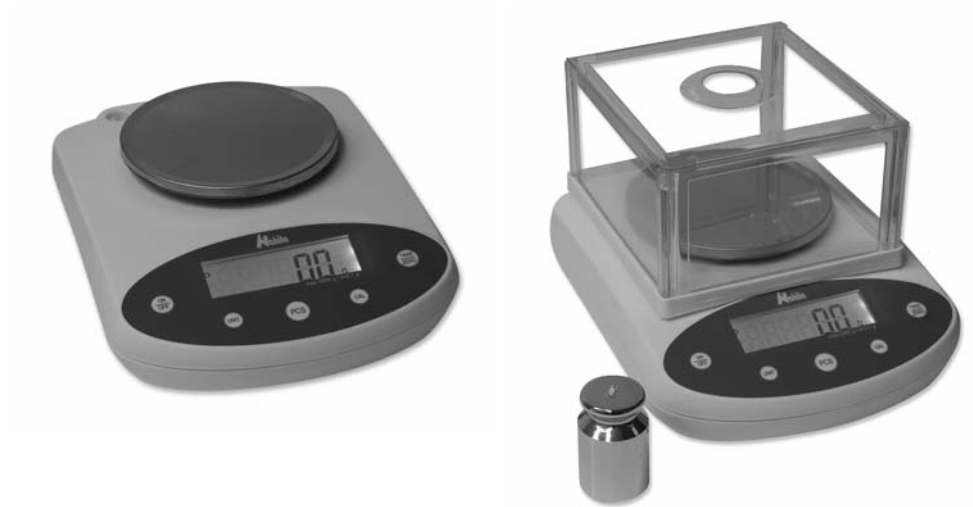SERIES | SERIES | SÈRIES 5061, 5062, 5063

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita digital balances, series 5061/5062/5063. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita digital balances 5061/5062/5063, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the balance avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the balance to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the balance.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es.
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

TROUBLE	CAUSE	SOLUTION
Value on display is unstable	Adapter is damaged The balance needs to be calibrated	Check operation with batteries, if balance works correctly then change the adapter. If the problem still persist contact NAhita Technical Service Department. Calibrate the balance
The value displayed is wrong	1 The balance needs to be calibrated 2 The balance has not been tared before weighing	1 Calibrate the balance 2 Tare the balance before weighing
Balance can not be calibrated	1 The calibration weight used is not the correct one 2 Zero is not adjusted	1 Use the proper calibration weight 2 See Tare function
The balance can not be tared	Lost of the reference zero	Calibrate the balance twice
	The weight of the object is lower than 0 g	Calibrate the balance
	The weight of the object is lower than 0 g	Calibrate the balance
The balance can not be tared	Lost of the reference zero	Calibrate the balance twice

INSTRUCTIONS ABOUT PROTECTION OF THE ENVIRONMENT

Nahita digital balances series 5061/5062/5063 can operate with batteries; once they are run down, please always deposit them in special containers used for this purpose or in a collection point. Batteries contain heavy metals (zinc, lithium, mercury, etc.) that could contaminate water, soils, etc.

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

If any liquid comes into contact with the electric parts of the balance please do immediately disconnect it from the net and send it to the technical service, as soon as possible, for its checking and adjustment.

Always use original components and supplies. Other devices can be similar but they can damage the equipment.

Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the balance and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

Note: According to the in force legislation regarding "Non-automatic weighting instruments" in which balances are included, by means of writ dating from 22nd October, 1994 (BOE 1/3rd/95), these balances must not be used for:

- Commercial transactions
- Calculating taxes, tariffs, rates, indemnities and other similar canons
- Judicial surveys
- Pharmaceutical medicine preparations, as well as analysis made in medical or pharmaceutical laboratories
- Determining the price or total amount in sale price or in pre-packaged preparations

7. TROUBLESHOOTING

TROUBLE	CAUSE	SOLUTION
No digits are displayed	1 The adapter is not connected 2 Batteries are run down	1 Connect the adapter 2 Change the batteries
Value is not displayed in a continuous way	Low battery	Change batteries

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

1. Spanish 2-11
 2. English..... 12-21
 3. French..... 22-31

INDEX OF CONTENTS


1. USES OF THE INSTRUMENT 13
 2. DESCRIPTION 14
 3. TECHNICAL SPECIFICATIONS 14
 4. INSTALLATION / SETTING UP..... 15
 5. FUNCTIONS OF THE BALANCE..... 17
 6. MAINTENANCE AND CLEANING 19
 7. TROUBLESHOOTING 20

1. USES OF THE INSTRUMENT

Nahita digital balances are a basic instrument at any laboratory to perform weighing tasks. Developed with the most advanced technology, they are reliable and durable equipments, and its solid, clean and ergonomic design is suitable also for industrial use.

2. DESCRIPTION

- 1.1 Pan
- 1.2 Digital display
- 1.3 ON/OFF
- 1.4 UNIT: selection of the weighing unit
- 1.5 PCS: piece counting function
- 1.6 CAL: calibration
- 1.7 TARE/ZERO
- 1.8 Bubble level

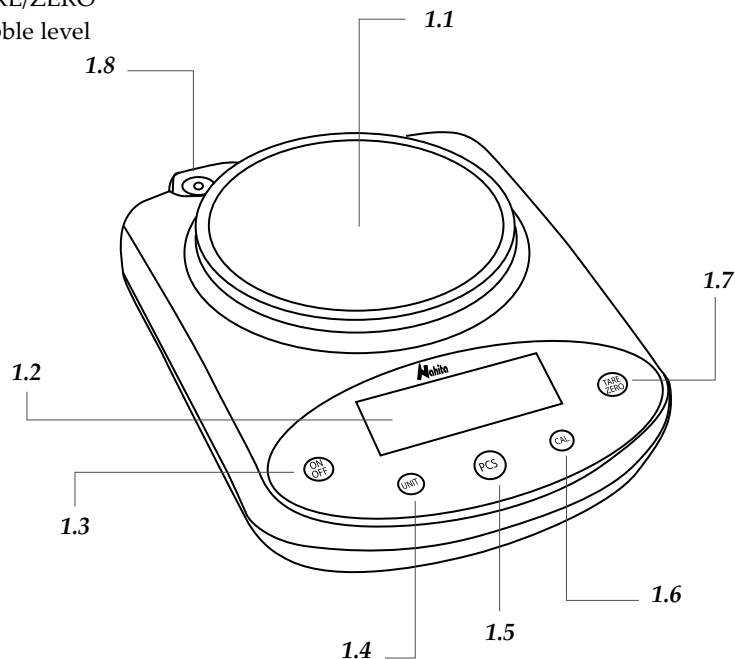3. TECHNICAL SPECIFICATIONS

Among the main features of the Nahita digital balances series 5061/5062/5063 it is worth noting:

- ◆ Quick setting up and stabilization (“o” indicator on display)
- ◆ Easy-to-use and easy-to-read display
- ◆ Weighing, tare, piece counting and calibration functions
- ◆ Weighing unit selection:: grams (g), carats (ct), ounces (oz), pounds (lb)
- ◆ Bubble level and screwed legs for a perfect leveling of the balance
- ◆ Hermetically sealed panel resistant to liquids splashes
- ◆ Operation through adapter or with battery (not included)

Security

- ◆ The balance must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the balance on top of a horizontal, plane and stable table making the enough free space around for a comfortable operation.
- ◆ Do not put the balance near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ When you are not using the balance for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece.
- ◆ Never try to repair the balance by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the equipment (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

6. MAINTENANCE AND CLEANING

To get the best results and a higher duration of the balance it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Prevent the balance from sudden movements and knocks, as well as from direct sunlight or air flows. The balance is a precision instrument, you must handle it carefully.
- ◆ The balance is supplied with an adapter. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never unplug the adapter by pulling the wire, do it from the base.
- ◆ Never use the balance in a wedged location as for example a shelf.
- ◆ Never use sharp objects as pens, etc... to press the buttons of the control panel; only use your fingers.
- ◆ Never place on the pan an object heavier than the maximum capacity of the balance, the sensor could be damaged.
- ◆ Neither submerge the balance nor spill liquids on it.
- ◆ If you use batteries, remove them when you are not using the balance for a long period of time.

1. Turn the balance on by pressing On/Off (1.3) and make sure that there is nothing on the pan (1.1).

2. Press Tare (1.7) and when "0.0 g" / "0.00 g" / "0.000 g" is displayed, press CAL (1.6) for approximately 5 seconds. -CAL- will be displayed and then the weight of the necessary calibration weight. Put the weight on the pan.

3. When the weight of the calibration weight is displayed, remove the weight; the balance will automatically register the new calibration and return to the weighing mode. "0.0 g" / "0.00 g" will be displayed.

Series 5061

Code	Calibration weight
55061250	500 g (no incluida)
55061310	500 g (no incluida)
55061320	1000 g (no incluida)
55061330	2000 g (no incluida)

Series 5062

Code	Calibration weight
55062230	200 g (included)
55062250	500 g (included)
55062310	500 g (included)
55062320	1000 g (included)

Series 5063

Code	Calibration weight
55063210	50 g (included)
55063220	100 g (included)
55063230	200 g (included)

After calibration the balance can operate as usual.

Piece counting function ("Mode PCS")

On mode PCS the balance can calculate the number of piece placed on the pan. For this function proceed as follows:

1. Press On/Off (1.3) to turn the balance on and wait until value "0.0g"/"0.00g"/"0.000g" is stable on display.

2. Press PCS (1.5); -COU- will be displayed and finally the value "10 pcs" flicks, that corresponds to the number of pieces that must be placed on the pan as a reference.

3. If any other reference value needs to be set, press CAL (1.6). The different values available (10, 20, 50, 100, 150, 200, 250, 500) will be displayed consecutively.

4. Put the selected number of pieces on the pan and press PCS (1.5) to confirm.

5. Remove the pieces (value "0" will be displayed) and put the sample on the pan to count the pieces that it contains.

6. The corresponding number of pieces will be displayed.

7. Press "PCS" to return to the weighing mode

Serie 5061 / 0.1 g

Code	55061250	55061310	55061320	55061330
Capacity	500 g	1000	2000 g	3000
Readability	0.1 g	0.1 g	0.1 g	0.1 g
Tare range	0-500 g	0-1000 g	0-2000 g	0-3000 g
Pan diameter	130 mm			
Power	Adapter / Battery (9V)			

Serie 5062 / 0.01 g

Code	55062230	55062250	55062310	55062320
Capacity	300 g	500	1000 g	2000
Readability	0.01 g	0.01 g	0.01 g	0.01 g
Tare range	0-300 g	0-500 g	0-1000 g	0-2000 g
Pan diameter	130 mm			
Power	Adapter / Battery (9V)			

Serie 5063 / 0.001 g

Referencia	55063210	55063220	55063230
Capacity	100 g	200 g	300 g
Sensibility		0.001 g	
Tare range	0-100 g	0-200 g	0-300 g
Repeatability		±0.002 g	
Linerity		±0.002 g	
Pan Diameter	90 mm		
Power	Adapter / Battery (9V)		

4. INSTALLATION / SETTING UP**Preliminary inspection**

- ◆ Unwrap the balance, take off the involving plastic and take off the polispán protection in which it comes fitted.
- ◆ Make sure that it does not present any damage because of the shipment. In case the balance presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping with all the accessories and documents supplied.

Please check that all the accessories are enclosed with the equipment:

- Windbreak removable (Series 5062)
- Calibration weight (Series 5062/5063)
- AC adapter
- User's manual

Installation

Before using the balance, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT

1. Put the balance on top of a horizontal, plane and stable table making the enough free space around for a comfortable operation.
2. Do not put the balance near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
3. Level the balance by turning the screwed legs so as the bubble level is perfectly centered (1.8)
4. Nahita digital balances series 5061/5062/5063 can operate with batteries or by connecting them to the electric current by an AC adapter.
 - a) **AC adapter installation** 
 - ◆ Connect the adapter into the plug on the back side of the balance. Then, plug the adapter into the electric current.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to people suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz \pm 10%.

b) Battery installation

- ◆ Remove the pan of the balance, and being careful not to damage the sensor, put the balance upside down and remove the lid of the battery compartment.
- ◆ Place the proper battery inside the compartment bearing in mind to put them correctly according to their polarity.

- ◆ Replace the lid of the compartment and finally, put the pan of the balance in its original position.

5. 5063 Series, before place the pan, remove the sensor's protector screw.

Setting up

- ◆ Press On/Off (1.3) to turn the balance on.
- ◆ When "0.0 g" / "0.00 g" / "0.000 g" is displayed, the balance is ready to operate.
- ◆ By pressing Unit (1.4) select the desired weighing unit.
- ◆ By pressing On/Off (1.3) again the balance will turn off.

Unit conversion

<i>Abreviature</i>	<i>Name</i>	<i>Conversion</i>
Kg	Kilogram	1000 g
g	Gram	1 g
mg	Miligram	0.001 g
ct	Carat	0.1999694 g
lb	Pound	453.592 g
oz	Ounce	28.3495231 g

5. FUNCTIONS OF THE BALANCE

Tare function

Tare function is used when a substance contained in a recipient is needed to be weight; for example liquids or solid reagents contained in a beaker to prepare a solution. For this function, proceed as follows:

- ◆ Put the empty recipient on the pan and press Tare (1.7), "0.0 g" / "0.00 g" / "0.000 g" will be displayed. When the stabilization symbol "o" is displayed the balance is tared and ready to weight the substance.

NOTE: The balance presents a tare range the same as its maximum capacity. This maximum must be taken into account since if the maximum capacity is tared nothing more could be weighted. (For example: if 450 g are tared in a balance with a maximum capacity of 500 g, only 50 g more will be weighted).

Calibration function

Before starting using the balance and whenever observing important variations in weighing or not correct weighings, balance should be calibrated. Some factors such as variation of gravity depending on latitude, change of working place as well as an inappropriate use of the equipment lead to the necessity of recalibrate the balance.

Important!! do not move the balance during calibration