4. ENTRETIEN ET NETTOYAGE

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4. MAINTENANCE AND CLEANING

Because the surface of glass fibers is coated with grease, the first time it is used, it should be heated slowly. When white smoke is seen, the power supply should be turned off till the smoke is scattered, repeat several times, until the electricity is turned on and smokeless, it can be used formally.

No flammable, explosive, corrosive liquid or gas can enter the electric heating mantle. The relative working environment humidity of the electrothermal mantle is not more than 85%.

This product must be grounded. It should be dry and not damp when stored. If there is humidity, the use should pay attention to induction electricity, do not touch with hands, after drying can restore its good insulation performance.

For proper equipment operation, it is necessary to follow some recommendations.

Note: All the usage rules mentioned above will be ineffective if continuous maintenance work is not carried out.

- » Follow the instructions and warnings provided in this manual.
- » Keep this manual on hand so that anyone can consult it.
- » Always use original components and spare parts. Other devices may appear similar, but their use can damage the equipment.
- » The heating mantle comes with a Schuko power cable; it should be connected to a grounded power outlet and kept within reach for emergency disconnection.
- » Do not attempt to repair it yourself; besides voiding the warranty, you may cause damage to the overall operation of the stirrer, as well as injuries to people (burns, wounds...), and damage to the electrical installation or nearby electrical equipment.
- » In case of a malfunction, contact your supplier for repair through the manufacturer's Technical Service.
- » For cleaning metal parts, stainless steel, aluminum, paints, etc., never use scrubbing pads or products that can scratch, as they can damage the equipment and shorten its lifespan.
- » For cleaning the equipment, we recommend using a lint-free cloth dampened with soapy water that does not contain abrasive substances.

ATTENTION! NO DEVICE FOR REPAIR WILL BE ACCEPTED IF IT IS NOT PROPERLY CLEANED AND DISINFECTED.

Attention! There is a risk of burns when the heating plate operates above 50° C. Do not touch the heating mantle, the flask, or the liquid.

Attention! For safety reasons, when heating flammable samples, the plate temperature should be lower than the sample's flash point.

NOTE: Due to the thermal transmission resistance of materials, the volume to be heated, and heat loss by radiation, the temperature in the container will always be lower than the set temperature of the equipment.

3.3. SAFETY

The multi-position heating mantle should be used by qualified personnel who are familiar with the equipment and its operation through the user manual.

- » Place the equipment on a horizontal, flat, and stable table, creating a clearance space of at least 30 cm on each side.
- » When working at temperatures above 50° C, do not touch the flask, the round-bottom flask, or the liquid, as it may cause burns.
- » Take extra precautions when working with liquids that reach temperatures above 100° C, such as oil.
- » Do not place the equipment near heat sources (lighters, torches, etc.) or expose it directly to sunlight. Avoid vibrations, dust, and very dry environments.
- » For safety reasons, when heating flammable samples, adjust the plate temperature below the sample's flash point.
- » When not using the equipment for long periods of time, make sure it is disconnected from the power supply to avoid possible accidents.
- » For any cleaning manipulation, component verification, or replacement of any component (e.g., replacing a fuse), it is essential to turn off the equipment and disconnect it from the power outlet.
- » Do not attempt to repair it yourself; besides losing the warranty, you may cause damage to the overall operation of the equipment, as well as injuries to people (burns, wounds...), and damage to the electrical installation.
- » Ensure that no water enters the control panel, even if it is properly insulated. If for any reason you suspect that water or any liquid has entered, immediately disconnect the equipment (see Maintenance).
- » Manufactured according to European directives for electrical safety, electromagnetic compatibility, and machine safety.

INDEX OF LANGUAGES

Spanish	2-6
English	7-11
French	12-16

TABLE OF CONTENTS

1. Product description	7
2. Technical data	8
3. Operation instructions.....	8
3.1. Preliminary inspection	8
3.2. Installation	9
3.3. Safety	10
4. Maintenance and cleaning.....	11

1. PRODUCT DESCRIPTION

Electric heating mantle is mainly used for liquid heating in laboratories of high schools and universities, petrochemical industry, medicine, environmental protection and other industries. It uses high temperature alkali-free glass fiber as insulating material, sealed resistance wire in the insulating layer, and then knitted into a hemispherical internal heater. This product has the advantages of large heating area, fast heating, good heat preservation effect, no open fire, and not easy to touch glassware.

2. TECHNICAL DATA

Reference	JGD010	JGD011	JGD012	JGD013
Positions	4	6	4	6
Capacity (mL)	250	250	500	500
Power (W) (for each position)	180	180	265	265
Voltage (V)	60-220V / 50-60 Hz			
Max temp	450°C (Surface temperature)			
Temp range	environment temperature ~350°C			
Work hours	continuously			

3. OPERATION INSTRUCTIONS

3.1. PRELIMINARY INSPECTION

Unpack the equipment, remove the polystyrene protection it comes packaged in, and take off the plastic wrap. Remove all protections and, without connecting the stirrer to the electrical supply, make sure it has no damage due to transport. If there is any damage, immediately notify your carrier or supplier. Keep the packaging, as returns should always be made in the original packaging with all supplied accessories.

Check the accessories that you should receive with the equipment:

- » 3x stirring rods
- » 6x double nuts
- » 2x rod holders
- » 4x clamps
- » Standard Schuko power cable
- » User manual

We only accept returns of equipment within 15 days of shipment and only if they are complete in their original packaging.

3.2. INSTALLATION

Before starting to use the instrument, it is advisable to familiarize yourself with its components and basic principles, as well as the functions of its controls.

READ THIS INSTRUCTION MANUAL CAREFULLY BEFORE OPERATING THIS EQUIPMENT IN ORDER TO ACHIEVE MAXIMUM PERFORMANCE AND LONGER DURABILITY.

Place the equipment on a horizontal, flat, and stable table. Create a clearance space of at least 30 cm on each side. Do not place the equipment near heat sources (lighters, torches, etc.) or expose it directly to sunlight, etc. During operation, no hazardous materials such as flammable liquids or pathological materials should be in this area.

Insert the AC power supply cable into the 220 V \pm 10% 50-60 Hz power socket with grounding and connect the other end to the stirrer connector.

Neither the manufacturer nor the distributor will assume any responsibility for damage to the equipment, installations, or injuries suffered by anyone due to non-compliance with the correct electrical connection procedure. The voltage should be 220 V \pm 10% 50-60 Hz.

When not using the stirrer for long periods of time, make sure it is disconnected from the electrical supply.

Startup

NOTE: Before connecting to the power supply, make sure the temperature control is turned off. To do this, the indicators should be off and the controls should be fully turned to the left.

Stirrer Connection

Connect the Schuko power cable to a grounded power outlet and keep it within reach for emergency disconnection.

Press the general/safety switch.

Sample Placement

Place a round-bottom flask with the solution to be heated in each position of the heating mantle you want to use and rotate the controller according to the required heating power.

Temperature Adjustment

To heat the solution, turn the temperature knob clockwise and the heating indicator will illuminate. Adjust the temperature intensity according to your needs. Note that the temperature knob does not have the function of setting a specific temperature; if needed, install the rod with the nut and place a thermometer (not included in the equipment) into the liquid to be stirred. This way, you can monitor the temperature.