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Tel(948) 310 513 - Fax(948) 310 500 - 312 071

e-mail: asistencia@auxilab.com <http://www.auxilab.es>

Thanks for choosing a NAHITA Balance, its high mechanical quality will allow you to make use of this equipment for a long time, nevertheless to get the best results and a higher duration of this instrument it is advisable to read carefully this manual and follow the processes of use.

In order to use this equipment, it is advisable to take into account the following points:

- ◆ This manual should be available for all the users of this equipment.
- ◆ Any doubts can be clarified getting into touch with the NAHITA
- ◆ As an accuracy instrument, it should be used very carefully avoiding the sudden movements and knocks.
- ◆ Never dismantle the different pieces of the balance to repair them, since it could produce a defective use of the whole equipment and a loss of the product warranty.

ATTENTION !!

IF AN EQUIPMENT IS NOT PROPERLY CLEAN AND DISINFECTED, IT WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

CONTENTS

1. USES OF THE INSTRUMENT	7
2. DESCRIPTION	8
3. TECHNICAL SPECIFICATIONS	8
4. INSTALLATION	9
5. OPERATING INSTRUCTIONS	9
6. MAINTENANCE	9
ANNEX I: WARRANTY INFORMATION.....	10


1. USES OF THE INSTRUMENT

NAHITA practical and comfortable balances enables a fast weighing and a direct reading. This model is of one single pan, and it is composed of jockey weights on the beams or arms, it is made of an alluminium alloy of quality. Thanks to the magnetic dumping, the time spent to weigh decreases.

It is especially designed to be used in teaching (physics and chemistry) as well as in chemical or pharmaceutical laboratories, and in those in companies.

2. DESCRIPTION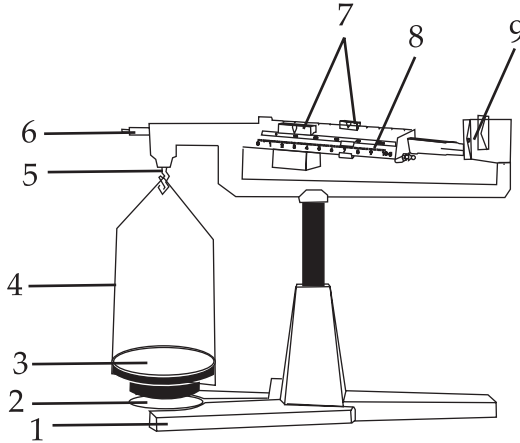

1. Base
2. Attaching for specific weight
3. Pan
4. Joke
5. Blade and bearing
6. Screw of adjustment
7. Jockey weights
8. Beams
9. Indicator

3. TECHNICAL SPECIFICATIONS

Features

<i>Capacity</i>	311 g
<i>Readability</i>	0.01g
	1st beam: 1 g. divided each 0.01 g
	2nd beam: 10 g divided each 1 g
	3th beam: 100 g divided each 10 g
	4th beam: 200 g Divided each 100 g
<i>Pan's diameter</i>	100 mm
<i>Dimensions</i>	420 x 175x 355 mm
<i>Weight</i>	1.9 Kg

Magnetic damping

Agate bearings

Steel blades

Stainless steel pan

It doesn't include additional weights

4. INSTALLATION


- ◆ Place the balance on a flat and solid surface, free of vibrations.
- ◆ Hang the jockey on the hole at the top and lay the pan on the support.
- ◆ Move the jockey weights towards the left so they are all on the holes that mark 0.
- ◆ The indicator must coincide with the mark on the right of the balance. Turn the screw of adjustment (6) in the leveling staff in the left extreme of the arm. This adjustment must be periodically done, as necessary.

5. OPERATING INSTRUCTIONS

- ◆ Weighting is done sliding the jockey weights along the beams, which are graduated and shows weighting's value.
- ◆ Put the object to be weighted the closer you can to the middle of the pan. You should first move the weight of more capacity to the right until the indicator turns down. Once the balance moves down, move the weight back one position to the left.
- ◆ Repeat this procedure with the rest of the beams as a decreasing order, what means: 1st beam: 200 g, 2nd beam: 100 g, 3th beam: 10 g and 4th beam: 1 g.
- ◆ Last beam's weight has to be slid up to a position that makes it coincide the indicator with the mark "0" on the right of the balance.
- ◆ Object's weight is the sum of jockey weight's position values.
- ◆ You should never put a weight over the maximum capacity of the balance: 311 g.

Determination of specific weight in solids

This model is provided with a platform to determine the specific weight in solids:

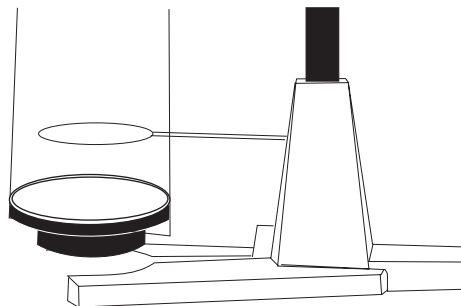Take the joke away (4) with the pan (3) of its usual position and put the specific weight platform (2) in the required height, holding it by the left extreme. After that, put the joke back (4) with the pan (3) of the sample at the top, having this way, the platform with the specific weight (2), into the joke (4) and over the pan (3).

The sample can be hung of the joke (5 pic. 1) thanks to a thin thread or a wire and a suitable recipient with water or other liquid.

Solids of a density higher than water's: the sample must be first weighted in air and then sunk into water.

$$\text{Specific weight} = \frac{\text{Weight in air}}{\text{Weight in air} - \text{Weight in water}}$$

Solids of a density lower than water's: you must do the following weightings:

- a) Body's weight in air.
- b) Total body's weight in air and container's weight in water.
- c) Total weight of both body and container in water.

$$\text{Specific weight} = \frac{a}{b - c}$$

6. MAINTENANCE

- ◆ Keep the balance clean and in good state. To have it cleaned, use a smooth gauze in a soapy solution. In the case of not accessible items, use a syringe or a dispositive blowing compressed air.
- ◆ Don't lay a weight over the maximum load.
- ◆ Never pour oil or lubricant on the blades or bearings. Try avoiding dust or any other substance.
- ◆ Don't through the package away, you can keep it away in case you move it or stop using it for a long time in future, or even to send it if it needs being repaired.
- ◆ Do not forget filling in and sending the warranty certificate in the following 15 days from purchase, to have your balance under one year warranty.
- ◆ In case the balance is damaged, contact your distributor to have it repaired by the technical service.
- ◆ We sincerely wish you enjoy your NAHITA balance. We advise you take care of it as the previous instructions. These balances are accurate instruments, so they must be properly, gently and carefully used.

WARRANTY

AUXILAB, S.L. warrants this equipment be free from defects in material and workmanship for a period of one year from date of purchase from AUXILAB S.L. or an authorised AUXILAB S.L. dealer. This warranty is limited to such purchaser and is not transferable.

During the year warranty period, AUXILAB S.L. will repair or replace, at its option, any defective equipments or pieces at no additional charge, provided that the product is returned, shipping prepaid, to AUXILAB S.L., an authorised dealer or an authorised service location. The purchaser is responsible for insuring any returned product and assumes the risk of loss during shipping. All replaced parts become the property of AUXILAB.

This limited warranty does not extend to any products which have been damaged as a result of accident, misuse, abuse or as a result of service or modification by anyone other than AUXILAB, an authorised dealer or an authorised service location.

Except as expressly set forth above, no other warranties are expressed or implied, including, but not limited to, any implied warranties of merchantability or fitness for a particular purpose. AUXILAB expressly disclaims all expressed or implied warranties not stated herein.

In order to obtain a free reparation or a free piece-replacement by our Technical Service over the above-mentioned period of time, it is absolutely necessary to submit the Warranty Certificate. This warranty doesn't apply for the maintenance of the equipment or for damaged caused by a natural, everyday use.

Some states do not allow the exclusion or limitation of incidental or consequential damages for consumer products, and some states do not allow limitations on how long an implied warranty lasts, so the above limitations or exclusions may not apply to you. This limited warranty gives you specific legal rights, and you may also have other rights which vary from state to state.

Dealer:			
Equipment:		Model:	
Serial Number:			
Date:			
Brief failure description:			

AUXILAB, S.L.

Polígono Morea Norte, 8 - 31191 BERIAIN (Navarra)

Tel(948) 310 513 - Fax(948) 310 500 - 312 071

e-mail: asistencia@auxilab.com <http://www.auxilab.es>