PRECISDIG 6001195
PRECISDIG 6001196
PRECISDIG 6001197
PRECISDIG 6001198

PRECISDIG 6001236
PRECISDIG 6001237
PRECISDIG 6001238
PRECISDIG 6001239

CON TAPA MAKROLON / WITH MAKROLON FOLDING LID:

PRECISDIG 6001186
PRECISDIG 6001188

PRECISDIG 6001187

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WATER AND OIL BATHS WITH OR WITHOUT FOLDING LID
WITH DIGITAL REGULATION AND READING OF TEMPERATURE AND TIME

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The standard equipment consist of the following components:

Pieza / Piece	Código / Code	Cant. / Qty.
Precisdig	6001195	1
Precisdig	6001196	1
Precisdig	6001197	1
Precisdig	6001198	1
Precisdig	6001236	1
Precisdig	6001237	1
Precisdig	6001238	1
Precisdig	6001239	1
Precisdig	6001186	1
Precisdig	6001187	1
Precisdig	6001188	1
Manual de instrucciones Instruction manual	80037	1

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Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Código / Code	6001195	6001196	6001197	6001198	6001236	6001237	6001238	6001239	
Temperatura máxima °C / Maximum temperature °C	100				200				
Capacidad litros / Capacity litres	5	12	20	45	5	12	20	45	
Medidas útiles cm. Tank dimensions cm	Alto Depth	15							
	Ancho Length	30	30	48	62	30	30	48	62
	Fondo Width	14	30	30	50	14	30	30	50
Medidas exteriores cm Overall dimensions cm	Alto Depth	20							
	Ancho Length	37	40	58	69	37	40	58	69
	Fondo Width	27	42	42	62	27	42	42	62
Consumo W. / Consump. W.	600	1100	1500	2400	1100	1400	2100	3800	
Peso Kg. / Weight Kg.	6	8	10	18	6	8	10	18	

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- * Remember that these are static baths.
- * Depending on the kind of fluid, the homogeneity can be altered.

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EQUIPMENT DESCRIPTION

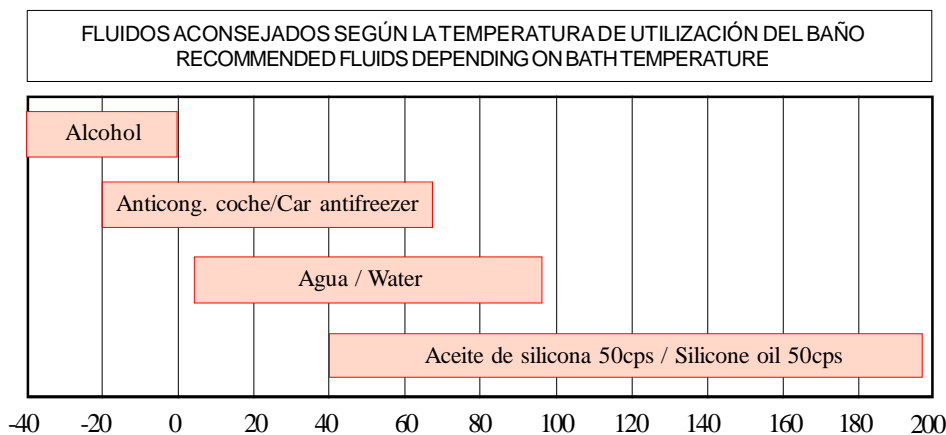
PRECISDIG baths with adjustable temperature from ambient +5°C up to 100°C or up to 200°C, depending on the model. To use the baths fill the tank with liquid until the level covers at least the heater element. The liquid or mixture to be employed must be adequated to the model of bath and the working temperature.

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The baths incorporate a drain tap for ana easy pouring.

As a guidance, here are some liquids (to use these liquids, follow the manufacturer's instructions):

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INSTALLATION

Place the bath on a flat, horizontal and level surface. Bear in mind the weight of the liquid when choosing the support. Provide for bath drainage.

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Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the equipment.

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Do not use the equipment if it is NOT earthed.

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If you have to change the plug bear in mind the following:

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- Blue cable: Neutral.*
- Brown cable: Phase.*
- Yellow/green cable: Earth.*

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It is recommended the use of decalcified water. DO NOT USE DEIONIZED OR DEMINERALIZED WATER as it has a great tendency to rust. To prolong the life of the bath and avoid rusting of the parts, dissolve 1 gr of CO_3Na_2 (Sodium Carbonate) per litre of water used in the bath.

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CONTROL PANEL:

1. *Main switch with indicator lamp.*
2. *Temperature mode indicator.*
3. *Time mode indicator.*
4. *Temperature/Time display.*

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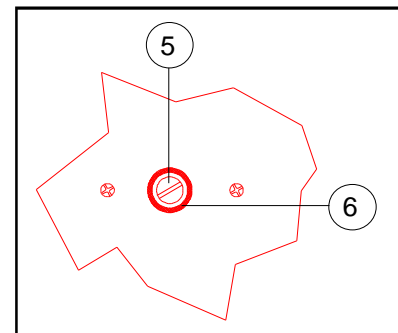
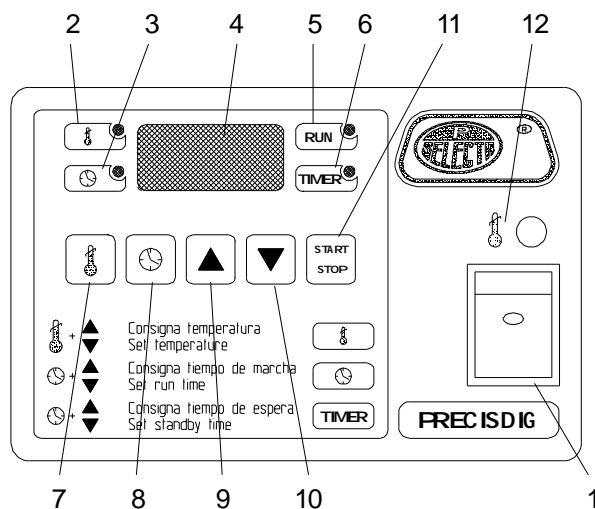
5. Starting up indicator lamp.
6. Waiting time indicator lamp.
7. Temperature selector push-button (Set of temperature from ambient +5°C up to 100°C/200°C depending on model).
8. Time selector push-button.
 - Set of starting up time: Working period between 1' and 9h 59', or up to 99.9h once the set temperature has been reached.
 - Set of waiting time: Period of time until the starting up, between 1 and 24 hours. Decimal fraction of hour (1.5 = 1hour 30 minutes).
9. Parameter increase push-button.
10. Parameter decrease push-button.
11. Start/Stop push-button.
12. Safety thermostat indicator lamp.

REAR PART

5. Safety thermostat regulation knob.
6. Safety thermostat reset knob (grey).

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The bath can work in 4 ways:

1. NORMAL MODE:

When pressing the main switch (1), placed in the lower part of the frontal, the display lights up (4). Pressing the push-button (7) and the (9) or (10) we preset the desired temperature. During this operation the indicator (2) lights up. Pressing the push-button (11) the bath will start and the indicator (5) will light up. To stop, press the push-button again (11) and the indicator will be turned off (5).

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2. TEMPORIZED MODE AT THE DISCONNECTION:

When pressing the main switch (1) the display lights up (4).

Pressing simultaneously the push-button (7) and the (9) or (10) we preset the desired temperature.

During this operation the indicator (2) lights up.

Pressing simultaneously the push-button (8) and the (9) or (10) we preset the disconnection time.

During this operation the indicator (3) lights up.

The disconnection time begins when the bath has arrived to the preset temperature.

Pressing the push-button (11) the bath will start and the indicator (5) will light up.

In this operation mode, the indicator (5) remains intermittent until the bath reaches to the preset temperature, once reached, the indicator (5) remains lighted up.

When the preset time has lapsed, the bath stops automatically and the indicator (5) will go off.

You can stop it manually pressing the push-button (11) again.

3. TEMPORIZED MODE AT THE STARTING UP:

When pressing the main switch (1) the display lights up (4).

Pressing simultaneously the push-button (7) and the (9) or (10) we preset the desired temperature.

During this operation the indicator (2) lights up.

Pressing simultaneously the push-button (8) twice and the (9) or (10) we preset the waiting time until the starting up. During this operation the indicator (6) will light up.

Pressing the push-button (11) the bath will start and the indicators (5) and (6) will light up.

In this operation mode, the indicator (6) remains lit up during the waiting time, this over, the indicator lamp (6) goes off and the bath starts automatically.

You can stop it manually by pressing the push-button (11) again.

4. TEMPORIZED WAY AT THE STARTING UP AND AT THE DISCONNECTION:

When pressing the main switch (1) the display lights up (4).

Simultaneously pressing the push-button (7) and the (9) or (10) we preset the desired temperature.

During this operation the indicator (2) lights up.

Simultaneously pressing the push-button (8) twice and the (9) or (10) we preset the waiting time until the starting up. During this operation the indicator lamp (6) will light up.

Simultaneously pressing the push-button (8) and the (9) or (10) we preset the disconnection time.

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During this operation the indicator (3) lights up. The disconnection time begins when the bath has reached to the preset temperature. Pressing the push-button (11) the bath will start and the indicator (5) will light up. In this operation mode, the indicator lamp (5) remains intermittent until the bath reaches to the preset temperature, once reached, the indicator (5) remains lighted up. Lapsed the preset time, the bath stops automatically and the indicator (5) will go off. You can stop it manually by pressing the push-button (11) again. Pressing the push-button (11) the bath will start and the indicators (5) and (6) will light up. In this operation mode, the indicator lamp (6) remains lit up during the waiting time, once over, the indicator lamp (6) goes off and the bath starts automatically.

ADJUSTMENT FOR THE SAFETY THERMOSTAT:

The low limit of working temperature for the safety thermostat is nearly 50°C. The desired temperature of the safety thermostat must be at least 5°C above the working temperature. The adjusting process of the safety thermostat is long. For that reason, the safety thermostat is useful when the users work habitually at the same temperature.

1. Turn the safety thermostat screw (5) clockwise to maximum.
2. Select with the temperature selector push-button (7) the desired temperature of the safety thermostat.
3. Turn on the bath and leave it to stabilize.
4. When the bath is stabilized, slowly turn the safety thermostat screw (5) anticlockwise until it clicks and the red indicator lamp (15) lights up.
5. Wait until the bath cools to reset the safety thermostat with (6) button.
6. Select the working temperature of the bath with temperature selector push-button (7).

MAINTENANCE

Before removing the cover disconnect the equipment from the mains.

The control elements and heating elements must only be manipulated by authorized personnel.

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The distilled or demineralized water has a great tendency to rust. To prolong the life of the bath and avoid rusting of the parts, dissolve 1 gr of CO_3Na_2 (Sodium Carbonate) per litre of water used in the bath.

CLEANING:

If normal tapwater is used calcareous deposits will appear in the bath and in the heating element. A heater covered with lime should not be used as it can lose heating power and deteriorate more quickly. Periodical cleaning of the inside of the tank and the heating element with a solution of 10% of 28% Chlorhydric acid or a solution of water and vinegar (Acetic acid) is advised. Rinse with clean water afterwards.

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with cotton duster.

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To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA, S.A.

DESCRIPCIÓN / DESCRIPTION	CÓDIGO / CODE							
Modelo / Model	6001195	6001196	6001197	6001198	6001236	6001237	6001238	6001239
1. Circuito 230V / Circuit 230V	29139				29151		29147	
2. Circuito 110V / Circuit 110V	29162		29163		29164	29165		29166
3. Resistencia / Heating element	39061	39051	39012	39123	39050	39060	39059	39124
4. Termostato / Thermostat	43031				43017			
5. Sonda Pt100 / Pt100 probe	43059							

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GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA, S.A.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, S.A. cancels the guarantee automatically.

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