DISPENSADOR+BURTEA ELECTRÓNICA DIGIPETTE SMARTFLOW
ELECTRONIC DISPENSER+BURETTE DIGIPETTE SMARTFLOW
DISPENSATEUR+BURETTE ÉLECTRONIQUE DIGIPETTE SMARTFLOW

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DIGIPETTE

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

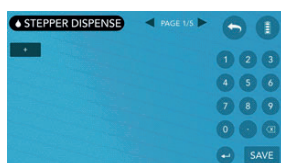
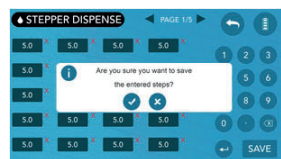
(Fig. 6)

10.4.3. PROGRAM

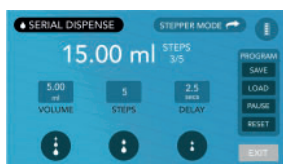
- To save programs in serial or step mode press the SAVE button in the program section (Fig. 1).
- Type in the program name and press SAVE.
- To load the stored programs (up to 10 per mode) with date and time press the LOAD function.
- To recall a stored program, press LOAD and select the program (Fig. 2).
- If the memory is full, a warning will be displayed, and it will be necessary to delete any existing program to save the new one.
- To delete programs, press the 'X' button next to the saved program.
- Use the PAUSE button to stop dispensing during operation (Fig. 3).
- Use the RESUME button to continue dispensing (Fig. 4).
- Use the RESET button to stop the program and return to the original data or exit (Fig. 5).
- Use the EXIT button to end the use; confirm rotation to recirculation mode, the piston empties the reservoir and returns to the initial position (Fig. 6). Switch off with ON/OFF button.


(Fig. 1)

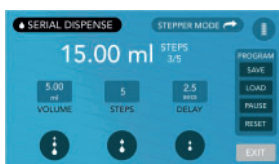
(Fig. 2)

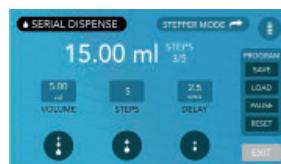
(Fig. 3)

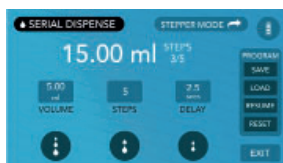(Fig. 4)

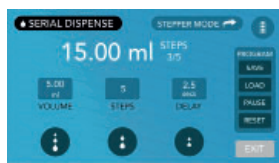(Fig. 5)

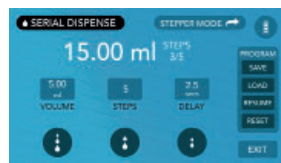(Fig. 6)

10.4.4. PIN ACTIVATION

PIN protection function to restrict user access to the configuration menu. This allows only authorised users to modify settings.

- Activate the PIN from configuration-enter password.
- Enter a 6-digit password and confirm with OK.
- After entering the password, the configuration page is accessed.
- The PIN is pre-activated, and the 6-digit password is printed on the panel box.

10.4.5. MODIFYING THE EXISTING PIN

The user can modify the existing password and set a new one by clicking on the CHANGE PASSWORD button.

- Enter the existing password in the panel box and confirm with OK.
- Enter the new password and confirm with OK.
- New password set: configuration access enabled.

10.4.6. PIN RECOVERY

In case of forgetting the PIN password, a default password is provided together with the user password, printed on the panel box. To set the password click on the FORGOT PASSWORD button.

- Enter the default password and confirm with OK.
- Enter the new password and confirm with OK.
- New password set: configuration access enabled.

10.5. SENSITIVE MEDIA

For photosensitive media (e.g. iodine solutions, potassium permanganate or silver nitrate), it is recommended to use the amber window included in the box.

- To replace the inspection window, disassemble the default window by pressing its top and removing it.
- Place the amber window in the slot at the bottom of the housing and press down until it clicks.


11. CALIBRATION

11.1. USER CALIBRATION PROCEDURE

The equipment has been laboratory calibrated to its nominal volume. However, due to possible variations in environmental conditions and the viscosity of the dispensed medium, it is recommended that a gravimetric test be performed every 3 to 12 months. The gravimetric volume test according to DIN EN ISO 8655-3 is carried out as follows:

- Enter the settings icon located in the lower right corner of the home screen.
- From the configuration menu, select Calibration.
- If the user detects a calibration problem at any dispensing speed, the user can select the specific speed to calibrate (Fig. 1).
- Fill the nominal volume with doubly distilled and deionised water by pressing the FILL button on the calibration screen of the control panel (Fig. 2).
- Dispense the loaded liquid by pressing the DISPENSE button on the calibration screen of the control panel (Fig. 3).
- Measure the dispensed liquid on a digital or electronic scale and enter the value in grams (g) using the electronic keypad available on the dispenser panel (Fig. 4).

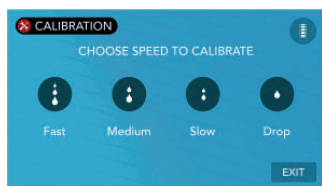
Note: This procedure is the same for calibrating the dispenser at fast, medium and slow speed for all volumes.

(Fig. 1)

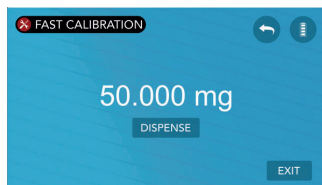(Fig. 2)

(Fig. 3)

(Fig. 4)

11.2 FACTORY RESET

- To reset the factory calibration settings, the user must access the Factory Reset option within the Factory Setup menu.
- Click on the confirmation symbol (✓) to reset the calibration to the original factory settings. The CAL label will disappear from all screens.

12. OTHER SETTINGS

12.1. CONTROL PANEL AND BRIGHTNESS ADJUSTMENT

- Click on the settings icon from the home screen.
- In the settings window, click on the BRIGHTNESS icon.
- From the BRIGHTNESS window, select the desired brightness level and click on the confirmation symbol to save it (✓).

12.2. LANGUAGE SETTING

- The user can select the preferred language from the Language setting. Select the desired language from the corresponding tab by clicking on it and confirm the selection by clicking on the confirmation symbol to save it (✓). You can switch to the following languages: English, German and Spanish.

12.3. SETTING THE DATE AND TIME

- In the configuration window, click on the Factory Settings icon.
- In the Factory Settings menu, select the date and time option.
- The user can set and modify the date and time using the available electronic keypad.

13. CLEANING AND MAINTENANCE

Whenever cleaning is required, circulate distilled water through the equipment. Completely fill the cylinder with distilled water and press the quick-dispense button to expel all the water.

Procedure for disassembling the spout:

- Unscrew the retaining nut by turning it counterclockwise and remove the spout. Clean the spout with deionised water.

Procedure for mounting the spout:

- Insert the outlet tube into the lower housing until it can no longer be moved forward.
- Screw on the retaining nut to complete the assembly, turning it clockwise.

Note: Use isopropyl alcohol applied to a cloth or cotton swab to clean the outer body.

14. TROUBLESHOOTING

Problem	Possible Cause	Solution
Difficulty moving the piston	Formation of crystals or dirt	Perform a cleaning cycle
Unable to refill	Filling valve clogged	Clean the fill valve. If the valve ball is stuck, use a 200 μ l plastic tip to loosen it.
Air bubbles in the instrument	Incomplete purge	Purge instrument again
	Loose or damaged fill tube	Firmly fix the telescopic filling tube. If necessary, cut the tube approximately 1 cm from the top or replace it.
	Filling tube is not immersed in the liquid	Fill the bottle or adjust the length of the telescopic filling tube correctly.
The delivered volume is less than indicated	The instrument is not fully primed	Re-priming the instrument
Instrument does not indicate any function	Internal error	Perform a reset
One-drop one-touch dispenser does not function	Instrument not purged	Make sure there are no air bubbles in the dispense tube or cylinder, purge correctly
	Use drop-by-drop dispensing button directly after filling instrument	After each fill, use the quick or medium button to dispense a small amount, then use drop-by-drop dispensing
Reading does not transfer from the instrument panel to the software	Driver is missing	Install the driver provided with the installation file.
	Cable is not connected or is damaged	Check the connection and, if the cable is damaged, ask your supplier for assistance