AUXILAB S.L.

Material de laboratorio  
Laboratory supplies

DECLARACIÓN DE CONFORMIDAD CE  
CE DECLARATION OF CONFORMITY  
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

Declara que el equipo | Declare that the equipment | Declare que l'appareil:

BALANZA GRANATARIO | EQUAL-ARM BALANCE | BALANCE GRANATOIRE

Código | Code | Code: 55001210, 55001220, 55001250, 55001310, 55001321, 55001322, 55001350.

Serie | Series | Série: 5001.

Cumple las siguientes Normas: | Meet the following Standards | Accomplit les normes suivantes:

EN 61010-1 | Requisitos de seguridad de equipos eléctricos de medida, control y uso en laboratorio. Parte 1: Requisitos generales.  
Safety requirements for electrical equipments for measurement, control and laboratory use. Part 1: General requirements.  
Règles de sécurité pour appareils électriques de mesurage, de régulation et de laboratoire. Partie 1: Prescriptions générales.

**BALANZAS DE 2 PLATOS  
BALANCE 2 PANS  
BALANCE 2 PLATEAUX**

**Nahita**

**Modelo / Model / Modèle 5001**

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**INDEX D'IDIOMES**

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**5. MAINTENANCE AND CLEANING**

- ◆ Preserve the balance from heavy shocks. Look for a dry place, away from dust, direct sun beams or air flows. Since it is a precision instrument, it should be handled carefully.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.
- ◆ Do not throw away the packing material. It might be useful in case you need to move it, to send it back for repairing or for long-term storage.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

*Note: According to the in force legislation regarding "Non-automatic weighing instruments" in which balances are included, by means of writ dating from 22nd October, 1994 (BOE 3/1/95), these balances must not be used for:*

- *Comercial Transactions.*
- *Calculating taxes, tariffs, rates, indemnities and othe similar canons.*
- *Juditial surveys.*
- *Pharmaceutical medicine preparations, as well as analisys made in medical or pharmaceutical laboratories.*
- *Determining the price or total amount in sale price or in prepackaged preparations.*

**INSTRUCTIONS ON ENVIRONMENT PROTECTION**

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

Thanks for choosing a NAHITA Balance, its high mechanical quality will allow you to make use of this equipment for a long time, nevertheless to get the best results and a higher duration of this instrument it is advisable to read carefully this manual and follow the processes of use.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

In order to use this equipment, it is advisable to take into account the following points:

- ◆ This manual is an integral part of the balance NAHITA and must be available to all users of this equipment.
- ◆ As it is a precision instrument, it should be handled carefully avoiding sudden movements and knocks,
- ◆ Never dismantle the different pieces of the balance to repair them yourself. It could lead to a faulty performance of the equipment.
- ◆ Any doubts can be clarified by contacting the NAHITA Technical Service by e-mail [asistencia@auxilab.es](mailto:asistencia@auxilab.es).
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

**ATTENTION!! IF EQUIPMENT IS NOT PROPERLY CLEAN AND DISINFECTED, IT WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

**INDEX OF LANGUAGES**

Spanish .....	3-6
English .....	7-10
French .....	11-14


**CONTENTS**

1. USES OF THE INSTRUMENT ..... 8  
 2. DESCRIPTION ..... 8  
 3. TECHNICAL SPECIFICATIONS ..... 8  
 4. INSTALLATION / SETTING UP ..... 9  
 5. MAINTENANCE AND CLEANING ..... 10  
 ANNEX I: CE CERTIFICATE ..... 16

**1. USES OF THE INSTRUMENT**

NAHITA balances series 5001, are very useful and handy. They are composed of two pans and a fast weighing. They are widely used in laboratories in teaching, pharmacy, industry, chemistry...

**2. DESCRIPTION**

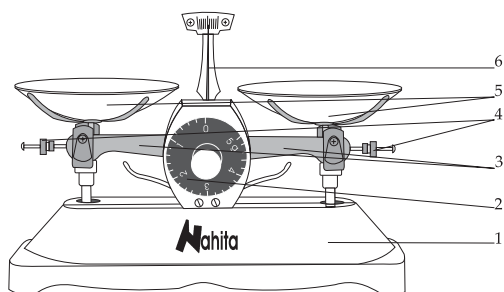1. Base
2. Pan
3. Needle
4. Screws of adjustment
5. Frontal control
6. Arm

**3. TECHNICAL SPECIFICATIONS**

<i>Code</i>	55001210	55001220	55001250	55001310	55001321	55001322	55001350
<i>Capacity</i>	100 g	200 g	500 g	1000 g	2000 g	2000 g	5000 g
<i>Sensibility</i>	0.1g	0.2g	0.5g	1.0 g	0.1 g	2.0 g	2.0 g
<i>Scale Range</i>	0 – 5 g	0 – 5 g	0 – 10 g	0 – 10 g	0 – 210 g	non-scale	0 - 50 g
<i>Scale division</i>	0.1 g	0.2 g	0.5 g	1 g	0.1 g	-	2 g
<i>Ø Pan</i>	84 mm	84 mm	110 mm	134 mm	140 mm	170 mm	210 mm
<i>Dimensions (mm)</i>	250x200x140	300x120x180	350x140x200			400x180x300	
<i>Weight</i>	1 Kg	1 Kg	1.6 Kg	2.5 Kg	6 Kg	6 Kg	12 Kg

\*Set of weights included (composition according to capacity)

**4. INSTALLATION / SETTING UP**

**Preliminary examination**

- ◆ Remove the balance from the packing material and examine it carefully to make sure no damage has occurred during shipping. If there is any damage, notify to your dealer/forwarding agency.
- ◆ Keep all packing material until you are sure that the instrument works correctly. All defective items must be returned in the original packing together with any other supplied accessories.

**We will not accept any equipment in return period unless it comes in its original wrapping.**

Verify you have received all the accessories that come with the balance:

- 2 pan
- Set of the weights
- Instructions manual

**Installation**

Before starting this instrument, carefully read this manual.

**BEFORE CONNECTING AND OPERATING THIS INSTRUMENT, READ THESE INSTRUCTIONS THOROUGHLY.**

- ◆ The instrument must be placed on a flat, steady, horizontal desk.
- ◆ Place the balance on a flat and solid surface. Put the pans (1.2) in the arm's ends (1.6).
- ◆ Get the balance adjusted. You can do it by checking the needle (1.3) points exactly the red mark on the scale (its mid point). If not, turn the screws of adjustment (1.4) located in the arm's extremes (1.6), under the pans (1.2). If you move the weight far away, you will add weight to that side, if you move it closer, it will decrease. When the needle marks the mid point of the scale, the balance is already calibrated. You should always calibrate it from the frontal control (1.5) in "0" position.
- ◆ Put the object to be weighted and the weights the closer you can to the middle of the pan.
- ◆ You should never put a weight over the maximum capacity of the balance
- ◆ These balances are supplied with a frontal control (1.5) that enables you to add up to 5 gr. in models 5001/100 and 5001/200, and 10 gr. in models 5001/500 and 5001/1000, without need to add a weight. This weight increase takes place on the pan on the right. That's why if you need to know a sample's weight, you should place the sample on the pan on the left and the weights on that on the right.
- ◆ Put the weights on the base carefully, adding weight bit by bit. As you know, the weight is between 0 and 5 gr. in models 5001/100 and 5001/200, and between 0 and 10 gr. in models 5001/500 and 5001/1000, and you will obtain the sharp weight with the frontal control (1.5).