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Referencia: QLB007

Descripción

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Code: QLB007

Description

Glass manometer tube, U-shaped, with stopcock near the end of one arm, mounted on a stand with back plate. Scale having graduations 80-0-80 with 2mm subdivisions. Size of back plate about 40×8cm (Height × Width).

Introduction

Manometer is the device used to measure the pressure. Manometer has two columns of a suitable liquid used to measure the difference in the pressure between a certain point and the atmosphere.

There are a variety of manometer design. But most commonly used design is glass tube of U shape with scale. Glass tube is filled with a liquid commonly mercury to remove the trapped air from the scale end of the tube. The glass tube is then position with the curved region at the bottom the mercury settles to the bottom.

Pressure P is defined as force per unit area

$$P = \frac{F}{A}$$

Liquid pressure is measured with respect to varying individual reference. When it is measured with respect to absolute zero or complete vacuum, it is called absolute pressure, when it is measured either above or below atmospheric pressure, it is called gauge pressure.

Manometer operate on the Hydrostatic Balance Principle

A liquid column of known height will exert a known pressure when the weight per unit volume of the liquid is known. The fundamental relationship for pressure expressed by a liquid column is

$$p = P_2 - P_1 = \rho gh$$

Where

p = differential pressure

P_1 = pressure at the low-pressure connection

P_2 = pressure at high pressure connection

ρ = Density of the liquid


g = acceleration of gravity

h = height of the liquid column

For gauge pressure P_2 is equal to zero, simplifying the equation to

$$p = \rho gh$$

If pressure is applied to the left side of the manometer the liquid recedes in the left leg of the instrument and rises in the right leg of the instrument. The liquid moves until weight of the liquid exactly balances the pressure. This is known as hydrostatic balance. The height of liquid from one surface to the other is the actual height of liquid opposing the pressure.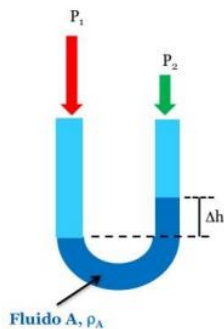

The pressure is always the height of liquid from one surface to the other surface regardless of the shape or size of the tubes.

Procedure

1. Place the manometer in the vertical position.
2. Fill the liquid in the manometer in such a way that the liquid level is same in both tubes of the manometer.
3. Close the stopcock attached at one side of the U-shaped tube.

4. Now attach the rubbing tube on the other side of the tube.
5. Now apply the air pressure either using syringe or any other method.
6. Due to the air pressure the level of the liquid changes in both tubes.
7. The difference in the height of the liquid's level in both tubes gives the height of the liquid opposing the pressure.
8. Now calculate the pressure using the equation

$$p = \rho gh$$

Advantages of manometer

- Simple in construction
- Low cost
- Very accurate and sensitive
- It can be used to measure other process variables