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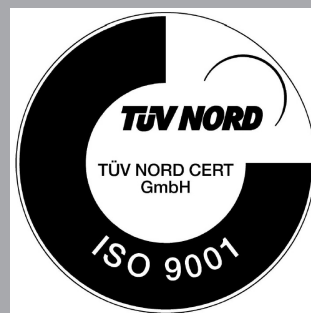
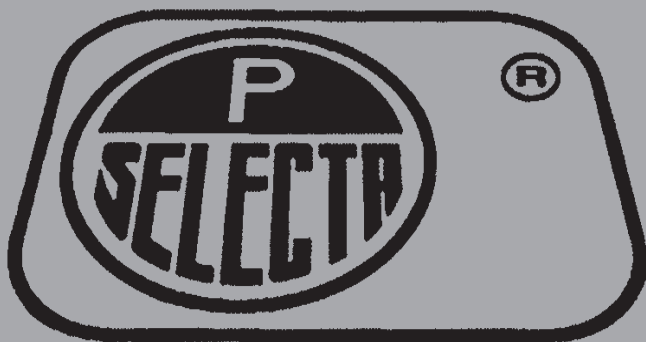
2 Installation

Unit description:

Shaker with vibrating motion action to provide homogenize effect.

Suitable for use in tubes.

Location:

- Place on flat and levelled surface. Suitable for unit weight.
- Let free 20cm, at least, around the unit.
- Near to power socket (1m).
- (The unit can cause vibrating effect on the table)

Connection to power supply:

- Check that power socket have suitable voltage, according unit plate.
- Power socket have earth connector. Unit must be grounded.
- Unit could be dangerous is not properly grounded.

3 Technical specification

Vibrating speed:	2800rpm.
Vibrating range:	6mm
Vibrating area	9mm
Power consumption:	50W
Fuse:	1A (Ø5 x 20mm)
Voltage:	230V AC ±10% (50/60Hz)
Weight:	2.5Kg
Dimensions:	170x120x170 mm

4 Operation

Press the main switch.

For shaker tube, hold tight the tube and press slightly over the vibrating surface.

PLACE DE TUBE CLOSED WITH A STOPPER iii

5 Warranty

The warranty covers the defective components or faulty assembling of the unit for 1 year since the purchasing date. Keep purchasing bill as the only valid document for the purchasing date.

The warranty does not covers the damages caused for a unit malfunction.

Return goods are not accepted without previous J.P. SELECTA,s.a. approval.

Check that the table surface is adequate to prevent displacement of the stirrer.

Hold tight the tube and press slightly over the vibrating surface

Use the suitable adapter for each sample tube.

J.P. SELECTA, s.a. 2015