**DESCRIPTION**

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*This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.*

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## DESCRIPTION

It is very useful for elementary biology class students for cutting their sections for microscopic examination of various organisms. Brass top plate with centered hole. The milled head at the base when revolved records 10 microns per click.

## INSTRUCTIONS

1. Unscrew the circular dial and set the rotating head below to the circular disc level and match the zero of circular dial with zero provided on neck pillar.
2. Now this is ready to work with organism sample. Place the organism sample merged with wax on the centre axle of hand microtome.
3. Cut out the rough organism sample from the microtome with the help of razor until it will get fully flat.
4. Now adjust the circular scale by just rotating the graduated knob to cut the organism sample as per your desired thickness.

### Least count:

The least count for the instrument is 10 microns.

### Precaution:

Keep the special attention on the time of performing the cutting of organism.

### Maintenance:

After use clean the microtome against any dust, water particles and residuals of organism sample.

### Environmental protection instructions

Do not dispose of this equipment with normal household waste at the end of its life cycle; take it to a collection point for the recycling of electrical and electronic equipment. It does not contain elements that are hazardous or toxic to humans but improper disposal will harm the environment. The materials are recyclable as indicated by the marking. By recycling materials or other forms of reuse of old equipment, you are making an important contribution to the protection of the environment. Please contact your local authority for advice on the collection points.