

VITLAB maneus®



The VITLAB maneus® Pipette Helper enables both left- and right-handers to operate all current volumetric and graduated pipettes from 0.1 to 200 ml easily and fatigue-free. Its safe and easy handling allows even inexperienced users **to adjust the meniscus precisely**.

With the design, unscrewing the adapter enables easy replacement of the hydrophobic membrane filter, which **protects the instrument against fluid penetration**.

The valve system is optimised so that liquids can be drawn up simply, without exerting pressure. The highly sensitive filling and discharge of liquids are controlled gently by the pipetting knob. Thus, the suction element provides rapid filling of the pipette (capacity: 50 ml in less than 10 seconds). The discharge bellows are used for the emptying (blow-out) of the pipette. The specially moulded intake cone ensures secure seating for all normal bulb and graduated pipettes (0.1 to 200 ml).

The VITLAB maneus® is simple to dismantle, easy to clean, and completely autoclavable at 121 °C (2 bar) according to DIN EN 285.

Included in delivery: VITLAB maneus®, replacement 3 µm membrane filter und operating manual.



Description	PU	Cat. No.
maneus®	1	1630500

Accessories for VITLAB pipeo® & maneus®

Description	PU	Cat. No.
Membrane filter, 0.2 µm, sterile, VITLAB pipeo®	1	1670647
Membrane filter, 0.2 µm, non-sterile, VITLAB pipeo®	10	1670648
Membrane filter, 3 µm, non-sterile, VITLAB pipeo®, VITLAB maneus®	10	1670650
Wall rack, VITLAB pipeo®	1	1670660

