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solution, as it may produce serious burns.

- You should act with the utmost care when working with liquids than reach temperatures higher than 100° C (for example, oil).
- During operation, do not place the magnetic stirrer near any warm supply (burners, blowlamps, etc.), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- When heating flammable substances, the hot plate temperature must be lower than the substance flashpoint.
- In case the magnetic stirrer is not to be used for a long period of time please make sure that it is unplugged from the net in order to avoid possible accidents.
- It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- Never try to repair the magnetic stirrer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the magnetic stirrer (see Maintenance).
- Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

## 5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is necessary to follow the processes of use.

**Note: All the process of use mentioned below will not have any value unless you keep a continued and careful maintenance.**

- Please follow the processes of use of this manual.
- This manual should be available for all users of these equipments.
- Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- The magnetic stirrer is supplied with a Schuko standard wire and it should be connected to a current wire provided with an earth wire, it should be handy to be disconnected in case of emergency.
- Never try to repair the magnetic stirrer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the magnetic stirrer (burns, hurts...) or damages in nearby equipments.
- In the event of breakdown please contact to your distributor to overhaul through Nahita Technical Assistance Department.

### Cleaning

- Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the magnetic stirrer and produce an early ageing of the equipment.
- Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

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Setting up

**Note: Before connecting the equipment, check that the speed and temperature knobs are off; the-  
refore, light indicators and must be switched off and the control knobs in the extreme left position.**

*1. Stirrer connection*

The magnetic stirrer is provided with a Schuko wire; it should be connected to a current wire provided with an earth wire, and it should be handy to be disconnected in case of emergency. Press the general/security switch.

*2. Placing the sample to be stirred*

In the middle of the plate place a glass with the solution to be stirred and a magnet inside, then increase gradually the stirring speed. You should do it this way to avoid the magnet to turn in an asynchronous way. Were this to happen, you should decrease the speed or -if necessary- switch the stirrer off. Then, switch it on again and increase the speed slowly.

*3. Setting the stirring speed*

Turn the speed knob clockwise and movement indicator will light up. Adjust the stirring speed according to your needs.

*4. Setting temperature*

In order to heat up the solution turn the temperature knob clockwise and the temperature indicator will light up. Then, adjust temperature's intensity according to your needs. Remember that the temperature knob does not present the function of adjusting to a specific temperature. In case you need a specific temperature you should assemble the rod and the bosshead, place a thermometer (not included with the equipment) and put it inside the solution. This way you will be able to control temperature.

Caution! There is danger of serious burns when the hot plate is operating at temperatures over 50° C. You must not touch the hot plate, nor the vessel or the solution.

Caution! When heating flammable substances the hot plate temperature must be lower than the substance flashpoint.

**Note: Due to thermal coefficient of vessel materials, solution volume and temperature loss by radiation, the solution temperature will always be lower than the temperature adjusted on the hot plate.**

Placing the thermometer

Placing the thermometer to control the temperature of the solution is an easy operation: first, you should assemble the two parts of the rod and insert it in its support at the back side of the equipment. Once the rod is placed you should put the bosshead at the desired height, hold it with the screw and put the clamp of the thermometer.

Once the stirring operation or your daily work is finished we highly recommend unplugging the equipment, as it is provided with electric components that have a little electric consumption, which involves an early ageing of the equipment, may provoke accidents due to heating production or unnecessary waste of energy.

Security

- The magnetic stirrer must be used by previously qualified staff that knows how it works thanks to the user manual.
- Put the magnetic stirrer in a horizontal, plane, stable table having a free space of at least 30 cm per side.
- When working at temperatures over 50° C you must not touch the hot plate, nor the vessel or the

**INDEX OF LANGUAGE**

Spanish .....	2-6
English .....	7-11
French .....	12-16

**INDEX OF CONTENTS**

1. USES OF THE INSTRUMENT .....	8	4. INSTALLATION / SETTING UP .....	9
2. DESCRIPTION .....	8	5. MAINTENANCE AND CLEANING.....	11
3. TECHNICAL SPECIFICATIONS.....	8		

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita magnetic stirrer with heating. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user. The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:


- This manual is inseparable from the Nahita magnetic stirrer with heating so it should be available for all the users of this equipment.
- You should carefully handle the Nahita magnetic stirrer with heating avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- Never dismantle the different pieces of the Nahita magnetic stirrer with heating to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the equipment.
- To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. This equipment is protected under the Warranties and Consumer Goods Regulation (10/2003).
- Overhaul is not covered by the Nahita magnetic stirrer with heating warranty.
- Operations made by non-qualified staff will automatically produce a loss of the Nahita magnetic stirrer with heating warranty.
- Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- Manufacturer reserves the right to modify or improve the manual or equipment.

## 1. USES OF THE INSTRUMENT

Nahita magnetic stirrers with heating series 690 are equipments perfectly suitable for daily work in laboratories that permit obtaining homogeneous solutions and improving test's speed. Thus, they are appropriate for any kind of analysis in biology, chemistry, medicine, etc.

## 2. DESCRIPTION

- |                            |                          |
|----------------------------|--------------------------|
| 1. Clamp for thermometer   | 7. Temperature knob      |
| 2. Bosshead                | 8. Movement indicator    |
| 3. Rod                     | 9. Temperature indicator |
| 4. Plate                   | 10. Speed knob           |
| 5. Control board           | 11. Power plug / socket  |
| 6. General/security switch | 12. Rod support          |

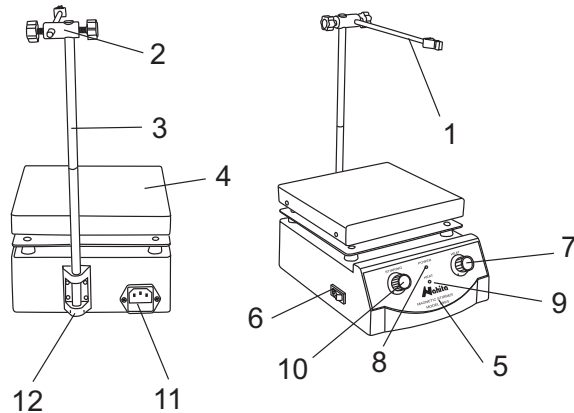## 3. TECHNICAL SPECIFICATIONS

Model	690/1
Reference	LBB001
Volume	50-2000 mL
Speed	100-2000 r.p.m
Maximum plate temperature	180° C ± 10%
Maximum sample temperature	80-85° C ± 10%
Plate	12 x 12 cm
Resistance consumption	180 W
Power consumption	200 W
Fuse	1 A

## 4. SETTING UP

### Preliminary inspection

Unwrap the magnetic stirrer, take off the polyspan protection in which it comes fitted and take off the involving plastic. Take off all the protective items and, without connecting the magnetic stirrer to the net, make sure it does not present any damage because of the shipment. In case it does, please communicate it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- Rod (31 cm y Ø 1 cm)
- Bosshead
- Thermometer clamp (20.5 cm long and 0.6 cm diameter)
- Teflonated magnet (28x9 mm)
- Schuko standard wire
- User manual
- Warranty Certificate

We will not accept any equipment for return unless it comes in its original wrapping and in the correct time limit

### Accessories

Other magnets: cylindrical with round edges, oval, octogonal. For more information please visit our website: [www.auxilab.es](http://www.auxilab.es)

### Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

**PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT**

Please put the magnetic stirrer on top of a horizontal, plane and stable table making a free space of at least 30 cm per side. Do not put the stirrer near any warm supply (burners, blowlamps...) nor expose directly to the sun. During operation, avoid inflammable fluids or pathologic material in the working area. The magnetic stirrer is supplied with a Schuko standard wire.

Please insert the wire that feeds the A/C in the base of current 220 V ± 10% 50-60 Hz provided with earth wire and to the other end to the magnetic stirrer connector.

**Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to people suffered by the improper use of the electric connection. The tension should be 220 V ± 10% 50-60 Hz.**

If you are not using the magnetic stirrer for a long period of time please make sure it is disconnected from the net.