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### 3. Autoclaving of PA, PC and PP centrifuge tube

PA begins softening at the temperature of 120°C, while PC and PP begin softening at 130°C. Generally, PA may be sterilized for 30 min at 115°C (0.7 kg/cm<sup>2</sup>), while PC and PP may be sterilized for 20 min at 121°C (1.0 kg/cm<sup>2</sup>). Too high temperature would result in deformation of centrifuge tube.

When autoclave is used, take the following steps:

- Place the centrifuge tube upright with opening facing upward. If the centrifuge tube is placed in an inclined or horizontal manner, it will deform due to the effect of gravity.
- Remove the threaded cover and inner cover to prevent deformation or crack of the centrifuge tube.
- Take the centrifuge tube only when the autoclave cools down to the room temperature.

### 4. Service life of centrifuge tube

The service life of plastic centrifuge tube depends upon the nature of sample, rotor speed and centrifugation temperature, etc. When the plastic centrifuge tube is used for centrifugation of aqueous samples (pH5-pH9), its estimated service life at the maximum rotation speed is as follows:

- High-quality centrifuge tube (PA, PC, PP): 30-50 times.
- Conventional centrifuge tube: about 10 times
- The service life of centrifuge tube is also related to the cleaning and sterilization conditions, which may reduce it.

Note: Never use any centrifuge tube with cracks.

## 11. RCF CALCULATION

Relative centrifugal force (RCF) can be calculated using the following formula:

$$RCF=1,118 \times r \times n^2 \times 10^{-5}$$

r- rotation radius, unit: cm

n- rotation speed, unit: rpm

## 12. AFTER-SALE SERVICE

To ensure safe and efficient operation of the centrifuge, periodical maintenance is required. If the centrifuge fails, don't attempt to repair it. Please contact your distributor to request technical service.

The tolerance is applied to tube pairs placed symmetrically in the rotor

**3. Precaution**

The density of sample that the centrifuge rotor can separate is less than 2.0g/ml. If the density of sample to be separated exceeds 2.0g/ml, please calculate the allowed rotation speed using the following formula:

**Allowed speed (rpm)** = Maximum speed × (2.0 (g/ml) /sample density (g/ml))½

**4. Autoclaving**

The high strength plastic rotor CANNOT be autoclaved

**10.2 Tubes**

**1. Cleaning and sterilization**

Condition		Material			
		PA	PC	PP	
Limpieza	Líquidos para limpieza	Acidic cleaning agent (pH5 or lower)	X	X	X
		Acidic cleaning agent (above pH5)	O	O	O
		Alkaline cleaning agent (above pH9)	O	X	O
		Alkaline cleaning agent (pH9 or lower)	O	O	O
		Neutral cleaning agent (pH7)	O	O	O
	70°C hot water	O	O	O	
	Ultrasónica	Neutral cleaning agent (pH7)	O	O	O
Sterilization	Autoclaving	115°C (0.7kg/cm²) 30min	O	O	O
		121°C (1.0kg/cm²) 20min	X	O	O
		126°C (1.4kg/cm²) 15min	X	X	X
	Boiling sterilization	15-30min	O	O	O
	Ultraviolet sterilization	200-300nm	X	X	X
	Gas sterilization	Ethylene oxide	O	X	O
		Formaldehyde	O	O	O

PA: Polyallomer PC: Polycarbonate PP: Polypropylene

**2. PC centrifuge tube cleaning**

PC material has relatively low chemical stability to alkaline solvent, therefore use of cleaning agent with pH value of over 9 should be avoided. Some neutral cleaning agents still have pH value of over 9 after being diluted as recommended by the vendor, therefore use of cleaning agent with pH value of 7-9 only is recommended.

**INDEX OF LANGUAGES**

Spanish ..... 2-14  
 English ..... 15-27  
 French ..... 28-40

**INDEX OF CONTENT**

Safety reminder ..... 15  
 1. Specifications ..... 17  
 2. Required operational conditions ..... 17  
     2.1 Basic operational conditions ..... 17  
     2.2 Transport and storage conditions ..... 18  
 3. Installation ..... 18  
     3.1 Location ..... 18  
     3.2 Connection of power cable and grounding ..... 18  
 4. Structure ..... 19  
 5. Control panel ..... 19  
 6. Rotor preparation ..... 20  
     6.1 Prepare the samples ..... 20  
     6.2 Place the samples into tubes ..... 20  
     6.3 Keep the tubes balanced ..... 20  
     6.4 Inspect the rotor ..... 20  
     6.5 Symmetrically load centrifuge tubes into the rotor ..... 20  
 7. Operation ..... 21  
     7.1 Normal operation ..... 21  
     7.2 Rcf operation ..... 23  
 8. Maintenance ..... 23  
 9. Troubleshooting ..... 24  
     9.1 Possible problems and solutions ..... 24  
     9.2 How to open the lid ..... 24  
 10. Rotor and centrifuge tubes ..... 25  
     10.1 Rotor ..... 25  
     10.2 Tubes ..... 26  
 11. Rcf calculation ..... 27  
 12. After-sale service ..... 27

**SAFETY REMINDER**

Common safety precautions

Carefully read the following safety precautions for a thorough understanding.

- Follow the instructions and procedures described in this manual to operate this centrifuge safely.
- Carefully read all safety messages in this manual and the safety instructions on the instrument.
- Safety messages are labeled as indicated below. They are in combination with signal words of “WARNING” and “CAUTION” with the safety alert symbol to call your attention to items or operations that could be dangerous to you or other persons using this instrument. The definitions of signal words are as follows:

**WARNING: Personal Danger**

Warning notes indicate any condition or practice, which if not strictly observed, could result in personal injury or possible death.

**CAUTION: Possible damage to instrument**

Caution notes indicate any condition or practice, which if not strictly observed or remedied, could result in damage or destruction of the instrument

**NOTE:** Notes indicate an area or subject of special merit, emphasizing either the product's capability or common errors in operation or maintenance.

- Do not operate this centrifuge in any manner not described in this User manual. When in doubt or have any troubles with this centrifuge, ASK FOR HELP.
- The precautions described in this User manual are carefully developed in an attempt to cover all the possible risks. However, it is also important that you are alert for unexpected incidents. Be carefully operating this centrifuge.

**WARNING**

- This centrifuge is not explosion-proof. Never use explosive or flammable samples.
- Do not install the centrifuge in or near places where inflammable gases are generated or chemicals are stored.
- Do not place dangerous material within 30cm around the centrifuge.
- Make sure to prepare necessary safety measures before using samples that are toxic, radioactive or contaminated with pathogenic micro-organisms at your own responsibility.
- If the instrument, rotor and/or accessories that has been contaminated by solutions with toxic, radioactive or pathogenic materials, clean it according to the decontamination procedure that you are specified.
- If you require services at site, please sterilize and decontaminate it in advance, and then notice the service center involved in the details of the particular materials.
- Do not handle the power cord or turn on or off the POWER switch with wet hands to void electrical shocks.
- For safety purposes, do not enter within 30cm around this centrifuge while it is in operation.
- While the rotor is rotating, never forcedly release the door lock.
- Unauthorized repairs, disassembly, and other services to the centrifuge except by our service center are strictly prohibited.

- This centrifuge must be located on one firm and level table.
- Make sure the centrifuge is horizontal before running.
- Make sure the angle between the door and cover is greater than 70 degrees when open the door.
- Be careful not put your fingers or hands between the door and cover when the door off.
- Do not move or relocate this centrifuge while it is running.
- If fluid spills in the rotor chamber, please promptly clean and dry with a dry cloth to avoid sample contamination.
- Ensure to remove any objects and fragments of the tubes dropped inside the rotor chamber before running this centrifuge.

2. In the event of a power cut

In case of unexpected power failure, the lid may be opened as follows:

Check whether the rotor is not moving.

Look at the rotor through the transparent lid and ensure that it is fully stopped.

Release the lock manually by the lever.

The lever is hidden under the bottom of the centrifuge.

Pull the lever to the right and release the lock. Then open the lid

**10. ROTOR AND CENTRIFUGE TUBES**

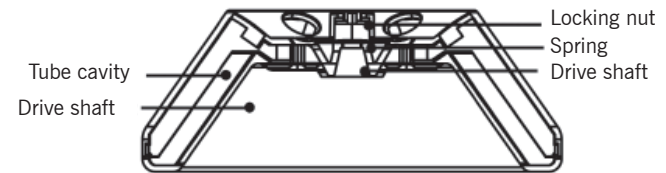
**CAUTION**

Carefully read the user manual and correctly install and use the rotor correctly. Don't exceed the maximum allowed speed of test tubes and adaptors. The maximum allowed speed of certain adaptors is lower than the maximum speed of the rotor.

**10.1 Rotor**

**1. Rotor structure**

**2. Rotor, tubes and adaptors**

Rotor	Rotor No.	Tube	Adaptor	Max speed (rpm)	Max RCF (×g)	Mass/Tube
Angular	1	Microtube 1,5/2,0 ml	White	5000	1817	2g / tube
		Ø13x75 mm	Orange	5000	2236	2g / tube
		Ø13x100 mm	Orange	5000	2236	2g / tube
		Ø16x107 mm	No	5000	2600	2g / tube
		Falcon type 15 ml	No	5000	2600	2g / tube

## Rotor

If the rotor is left unused for a prolonged period of time, remove the rotor from the centrifugal chamber store it.

Use clean water to rinse the rotor cavities if there is sample leakage from the centrifuge tubes. Dry the rotor thoroughly and apply a thin layer of silicone grease on it.

Clean the rotor regularly (every 3 months recommended). Ensure the cavities and the drive shaft hole are clean. Apply a thin layer of silicone grease.

## 9. TROUBLESHOOTING

### 9.1 Possible problems and solutions

This centrifuge is capable of onboard diagnosis. When the centrifuge fails, the time display window will indicate the failure code, leading to the immediate identification of possible failure causes.

Failure	Possible cause	Solution
No display after power ON	No power supply to the power socket.	Eliminate the failure and reconnect the power supply.
Abnormal vibration	Misfitting between the rotor and the drive shaft. The tubes are not balanced.	Mount the rotor properly again. Balance the tubes with scale and load them again.
Alarm code indicated on the time display	E-02 Lid failure	The lid is opened while running. is pressed when the lid is opened. Immediately close the cover. Close the outer lid before operation.
	E-09 Imbalanced rotor	The tubes are not balanced. Balance the tubes with scale and load them again.
	Others	Check the service manual. Contact your distributor.

Error codes E-01 / E-09 are related to erroneous operation. The centrifuge may continue running after elimination of the failure.

### 9.2 How to open the lid

- When turned ON

When the centrifuge is turned ON, the lid opens automatically.

At the end of the spin cycle, an audible signal will sound and the lid will open automatically.

If the lid is closed, press  key to open it.

### ■ Cautions on rotor

(1) Always check for damages on the rotor surface before using it. Do not use the rotor if an abnormality is found.

(2) Do not set the centrifuge speed beyond the allowable speed of the rotor kits (rotor or adapters).

(3) Do not exceed the allowable imbalance.

(4) Use the rotor and tubes within their actual capacities.

■ If any abnormal condition occurs during operation, please stop it immediately and contact your distributor to request technical service.

■ Vibrations are likely to damage the centrifuge, contact the service center if abnormality observed.

## 1. SPECIFICATIONS

Maximum rotation speed	5000rpm(300-5000rpm) step: 10rpm
Maximum relative centrifugal force (RCF)	2600×g, step: 10×g
Maximum capacity	6x15mL (Falcon type tubes)
Timing	30s-59s / 1min-99min / Continuous running (HOLD)
Drive motor	Brushless DC motor
Safety performance	Door lock, overspeed detection and automatic internal diagnosis
Power	Single phase, 100-240V, 50/60Hz
Dimensions (mm)	(Width) 300× (depth) 240× (height) 180
Weight	5.2kg
Acceleration/Deceleration	1 speed ↑ 2 speeds ↓
Noise	≤56dB
Other functions	Rotation speed/RCF switch; sound notification; 2 predefined and customizable programs

## 2. REQUIRED OPERATIONAL CONDITIONS

### 2.1 Basic operational conditions

■ Power supply: 100-240V, 50/60Hz

■ Ambient temperature: 2~40°C

■ Relative humidity: ≤80%

■ No vibration or air flow present nearby that might affect performance.

No conductive dust, explosive gas or corrosive gas exists in ambient air.

## 2.2 Transport and storage conditions

- Range of ambient temperature: -40°C-55°C
- Range of relative humidity: ≤93%
- The centrifuge must remain upright while in transit, suitably protected using wooden box.
- Lift the centrifuge by the chassis only

## 3. INSTALLATION

Users must follow the installation instructions to ensure safety and optimum performance. Before moving the centrifuge, the rotor must be removed.

### WARNING

Improper power connection might damage the centrifuge.  
Before connecting the power supply, please check the power supply for compliance with the requirements.

### 3.1 Location

- This centrifuge must be mounted on a solid, flat and tabletop with contact between the four feet of the centrifuge and the tabletop. Don't mount the centrifuge on any sliding tabletop, otherwise significant vibration might occur.
- The ideal ambient temperature is 20°C ±5°C and the ambient temperature should not be more than 30°C. Avoid direct sunlight on this centrifuge.
- Place the centrifuge properly and keep >30cm space from the rear and >10cm space on both sides to ensure effective air cooling.
- Ensure that there is no water leakage/ heat loss near the centrifuge as it may cause the rise in temperature and thereby leading to centrifuge failure.

### 3.2 Power cable connection and grounding

### WARNING

Don't touch the power cable with wet hands and avoid electric shocks.  
Ensure that the centrifuge is well grounded.

Ensure the rated current of the power socket is correct (>10A) and it complies to local electrical safety regulations before connecting the device to power supply.

## 7.2 RCF operation

Turn power switch ON and set the centrifuge to RCF (relative centrifugal force) mode.

- The relative centrifugal force set should not exceed the maximum relative centrifugal force allowed by the centrifuge tube and its adaptor.
- The relative centrifugal acceleration is calculated based on the maximum centrifugal radius and operating speed of the rotor. (See 12RCF calculation).
- Press the parameter knob to choose the rotation speed unit as xg. If the RCF value flashes it indicates that the RCF value can be set as it is in adjustment mode.
- Turn the knob to adjust the relative centrifugal acceleration, in increments of 10xg.
- Wait for 7 seconds to update the settings.
- See Section 7.1 for operation instructions.

### 7.3 Set deceleration level

Press and hold the knob for 5 seconds, after a beep sound, you can adjust the deceleration between 2 levels by rotating the knob. Level 1 is the default, and it is suitable for most applications. Set it to Level 2 for slower deceleration.

## 8. MAINTENANCE

### 8.1 Cleaning

### CAUTION

Possible damage to the centrifuge if you clean or sterilize it not following the instruction in this manual.  
Disconnect the power supply before cleaning the centrifuge.

#### Centrifuge

The color of housing might change and the label thereon might fall OFF if the centrifuge is exposed to ultraviolet for a prolonged period of time and hence cover the centrifuge with cloth to avoid exposure to light.

Clean the centrifuge using a cloth/ sponge soaked with neutral cleansing agent in case if it is dirty after use.

The centrifuge can be sterilized using cloth soaked with 70% alcohol solvent

#### Centrifugal chamber

Never pour water or other solvents directly into the centrifugal chamber as they might enter the drive unit and cause corrosion or damage to the bearings.

#### Drive shaft

It is suggested that the drive shaft be subjected to periodical maintenance by wiping it using soft cloth and applying a thin layer of silicone grease on it.

#### Outer Lid

Clean or sterilize the outer lid in the same manner as mentioned under subsection (1) centrifuge

Set the run time

- Press the parameter key  to select the time parameter and wait until its value blinks.
- Rotate the knob and set the timer to desired setting within the range of 30s-99min.
- Set the timer to HD to disable it and run the centrifuge in a mode without time limit.

Modify the predefined program

- Press the program #1 key  (or program #2 key ) to load the settings of this predefined program. Press it again and hold it for 5 seconds to edit the settings. Press it again and again to cycle between speed, RCF and timer setting. Rotate the knob to adjust the parameter of current setting. Wait for 7 seconds and the settings will be updated for the corresponding predefined program.

**3. Start run**

Press the  key to start running

- The rotor will start rotating only when the lid is locked.
- Timer starts to count down as soon as the speed reaches the set value.

Inquire and change the settings

- The setting can be modified when the centrifuge is running at the set speed.
- Press the knob to go back to settings screen. Follow the instructions in 7.1.2 to review and modify the settings. Wait for 7 seconds, the settings will be updated, the display will go to the running screen again and the centrifuge will run on the new settings.
- The timer will continue to count down without reset

Error message

- The centrifuge will automatically stop if any failure occurs while running, with the error code shown in the timer section on the LCD display. Please look up the cause of failure and take appropriate action in Table 9.1.

**4. Stop running**

When the time is up or you press the Start/Stop key  while running, the centrifuge will stop running.

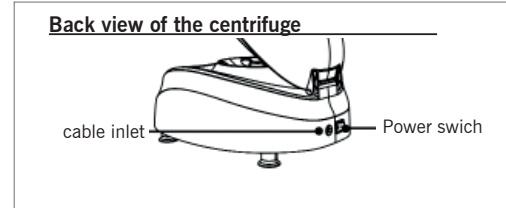
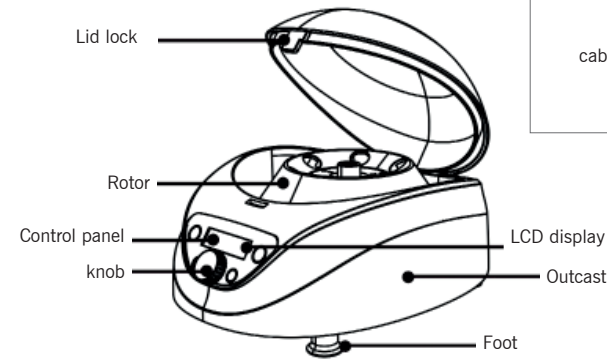
- The centrifuge will beep to notify the user when the rotor fully stops.

The lid lock will be released when the rotor fully stops.


- Press the  key to release the lock if the lid is still locked.
- The current settings will be stored. It will be loaded automatically as soon as the centrifuge is switched ON next time.

Retrieve the samples.


**4. STRUCTURE**

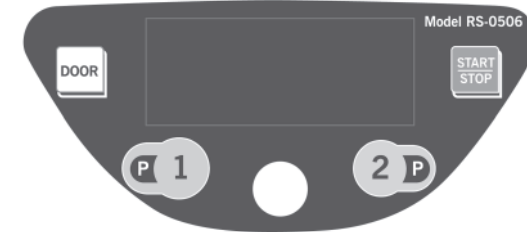**5. CONTROL PANEL**

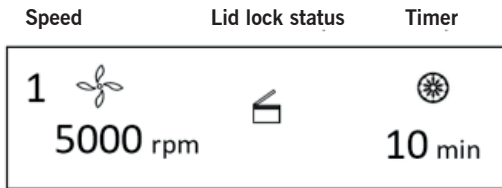Button	Name	Function
	Door lock open key	When the rotor is fully stopped, press this key to release the door lock. When the rotor is still running, the door will remain locked.
	Program #1 key	Press this key load/edit the predefined program #1.
	Parameter knob	Rotate this knob to adjust the parameter setting of speed, RCF, timer or deceleration.
	Program #2 key	Press this key load/edit the predefined program #2.
	Start/Stop key	When the centrifuge is ready to run, press this key to start running; While the centrifuge is running, press this key to stop.

The main interface of LCD display is shown in Figure 5.2. In this example, the speed is set at 5000rpm, the set running time is 10 minutes and the lid is open.

When the speed icon rotates, it indicates that the machine is in run mode. The icon rotates quicker at higher centrifuge speed.

The timer section shows the ratio of elapsed time to the total time.

## 6. ROTOR PREPARATION

### 6.1 Prepare the samples

### 6.2 Place the samples into tubes

The amount of sample should not exceed the working volume capacity of the centrifuge tube.

#### CAUTION

Adding excessive samples into the centrifuge tube will result in leakage, therefore don't add excessive samples.

### 6.3 Keep the tubes balanced

Although this centrifuge may be used with visually confirmed balance, it is suggested that samples be weighed using a balance to ensure balanced centrifuge tube in order to prolong the service life of centrifuge.

Although the imbalance to some extent is allowed (See 10.1.2 Rotor and adaptor), don't run this centrifuge under poor balance conditions.

### 6.4 Inspect the rotor

Check the rotor for any damage before use.

Avoid using the rotor with scratches.

Never use the rotor of other brands on this centrifuge.

Do not expose the rotor to direct sunlight/ultraviolet.

### 6.5 Symmetrically load centrifuge tubes into the rotor

Ensure to tighten the rotor to the main shaft firmly and the lid is secured properly on the rotor. Otherwise, the rotor might fall OFF while the centrifuge is in operation, resulting in centrifuge or rotor damage.

## 7. OPERATION

### 7.1 Normal operation

Turn ON the power switch, the most recent settings will be loaded and shown on the LCD display as in Fig. 7.1.

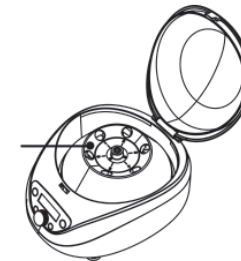
The speed set to 5000rpm and the timer set to 10min.

The lid lock is released.

### 1. Rotor installation and replacement

Locking nut

When mounting the rotor, ensure the full contact between the rotor and the motor shaft.

After placing the rotor in place, rotate the rotor gently with hands to check for normal operation of the rotor. If necessary, adjust the rotor position once again.

Hold the rotor with one hand, and rotate the locking nut clockwise using the rotor wrench with the other hand. Tighten the rotor and main shaft firmly.

Load the tubes, close the lid and run the centrifuge.

Dismantle the rotor in a similar manner with opposite rotating direction (counterclockwise).

Ensure the rotor is properly mounted on the centrifuge.

### 2. Set operating parameters

The parameter knob is used to adjust the parameter settings. Press the knob gently to select the setting, make it blinking on the display and ready to be adjusted. Rotate the knob clockwise to increase and counterclockwise to decrease the parameter. Rotate the knob quicker to adjust the parameter more dramatically. The minimum rotation speed increment is 10rpm, the minimum RCF increment is 10g and the minimum timer increment is 1second.

#### Set the rotation speed

Press the parameter knob to choose the rotation speed parameter value in rpm. The speed parameter should blink on the display.

Rotate the knob to adjust the rotation speed value to desired setting.

The minimum speed is 300rpm, and the minimum increment is 10rpm.

The parameter increases or decreases in a cyclic way. If you continue rotating the knob when the parameter hits the upper limit, it will start from the lower limit again.