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English version

1 Safety

The Equipment comes complete with safety features.

This manual indicates areas of possible risk.

1.1 Safety Icons

Identifies risk situations and safety measures that should be adhered to. The icons make reference within the paragraph, marked as a lined grey.

Risk danger

Risk danger

Beware, follow instructions as indicated.

Electrical risk

Risk of electrical shock if accessing zones are shown with this icon.

Beware, follow instructions as indicated.

Important information

- Important information in obtaining the best results or optimum performance of the equipment.
- Important information to extend the life of the equipment and maintain its optimum performance.

1.1 Risks which the operator is expoused:

- Risk of infection by handling of samples
- Risk of parts ejected
- Electric risks

1.2 Personnel training

This kind of equipment only must be used by personnel conveniently trained who read the instruction manual.

The biological samples handling requires special skills.

1.3 Use of hazardous products

When is known that the products to centrifuge are toxic, radioactive or when them be contaminated by pathogenic microorganismit, personnel must observe the following safety measures:

- Containers which ensure the sealing of these products and specified by the manufacturers must be used.
- When Risk Group II products be handled, as the identification of the «Biological safety for laboratories manual» edited by the World Health Organization, follow its instructions.
- In case of work with products belonging to higher groups it is necessary a protection level higher.
- It is forbided to use the following products wits this centrifuges:
 - Flammable or explosive products.
 - Products which may produce violent chemical reactions.

2 General information

- 2.1 Handle the equipment with care. Unpack and check that all items coincide with the delivery note. If you see any discrepancy or damage of any kind, notify your distributor as soon as possible.
- 2.2 Read this manual before using the equipment.
- 2.3 These instructions are an important part of the equipment and should be made available to all users.
- 2.4 If in any doubt or clarity on the use of this instrument, contact your local distributor or the service department of J.P. Selecta, s.a.
- 2.5 **ATTENTION, EQUIPMENT NOT CLEAN AND DISINFECTED SHALL NOT BE REPAIRED.**
- 2.6 Any unauthorised modification, lack of maintenance or deliberate damage to the machine, transgresses the directive 89/655/CEE where the manufacturer will not be responsible for damages that could occur due to undue care and attention.
- 2.7 Do not use the equipment for fluids that can produce vapours that can form inflammable or explosive mixtures.

3 Technical specifications.

Voltage supply 115-230V 50/60Hz according to the characteristics plate indications.

- Code:	7002356
- Dimensiones (cm):	
High:	20
Width:	28
Depth:	34
- Power (W):	170
- Speed max. (rpm):	5000
- RCF max.:	2515
- Weight (kg):	6

4 Contents list

4.1 Standard contents

The standard equipment consists of the following items:

- Centrifuge:	7002356
- Angle head for 8 tubes 15ml	30098
- 8 Aluminium buckets of 15ml	7001500
- 8 Glass tubes of 15ml	7000772
- Power supply cord	7140
- Instruction manual	80224

5 Introduction

The centrifuge Centro-8-BL is a tabletop centrifuge with a free maintenance induction motor and electronic digital control of speed and working time, controlled acceleration and deceleration, temporized mode from 1 minute to 60 minutes and safety electromechanic lock.

6 Equipment description


6.1 Equipment

The Centro-8- BL centrifuges have the following features:

- Electronic digital control circuit board
- Induction motor (brushless)
- Angle rotor 8 tubes of 15 ml
- Electromagnetic safety lock


6.2 Features

- Regulable speed until 5000 rpm
- Max RCF 2515 g
- Speed reading on the display (resolución 10 rpm)
- Working time reading on display (resolución 1 min)
- Instant start/stop key to the set speed
- Electronic lid opening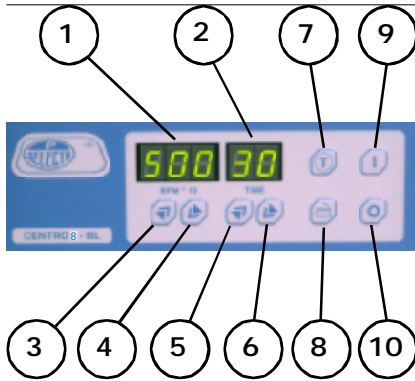

6.3 Control panel

1. Speed display (rpm)
2. Time display (min)
3. Increase speed value key
4. Decrease speed value key
5. Increase time value key
6. Decrease time value key
7. Instant start/stop key
8. Start key
9. Stop key

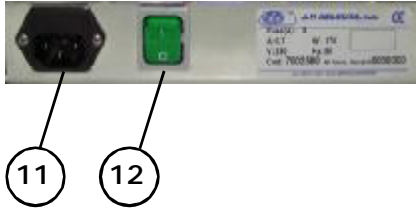6.4 Rear part

10. Power supply socket
11. Main switch

FREE SPACE AROUND THE CENTRIFUGE

Safety zone:

Leave a free space of 30 cm around of the centrifuge.

When the centrifuge is working:

- Do not place dangerous products in the safety zone.
- The personnel must be out of safety zone.

7 Installation

7.1 Placement

- Place the piece of equipment near of a socket keeping in mind the its amperage on a leveled surface enough resistant to support its weight.
- The surface of placement must allow the sticking of the legs of the centrifuge.
- Left a free spece of 30 cm around the piece. (See danger indication)
- This kind of equipment may produce vibrations which could affect nearby equipment.

DO NOT INSTALL THE CENTRIFUGE IN AREAS WHERE INFLAMMABLE LIQUIDS ARE STORED OR SPECIAL PROTECTION AREAS.

7.2 POWER SUPPLY CONEXION

Make sure that the voltage of given by the socket is the indicated on the feature label attached on the rear part.

Do not use the equipment unless it is correctly earthed.


If you replace the plug keep in mind the following:

Blue wire:	Neutral
Brown wire:	Phase
Yellow/green wire:	Ground

8 Operation

STARTING UP

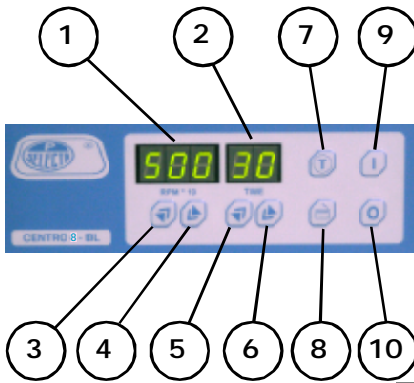- Connect the equipment to the power supply
- Turn on the main switch (12). The displays (1) and (2) will light.
- Press key (8) to open the lid
- Insert the tube in the rotor in balanced compensate way
If only use 2 tube insert them in opposite sides of the rotor(not contiguous)

Do not use the centrifuge:

If the speed showed on the screen does not stabilize never.

(In this case the real speed could be different than the showed by the screen)

Contact with the technical service of J.P. Selecta, S.A.

Never place one or three tube. Always two or four.

- Close the lid pressing on it until heard the sound of the lock.
- Pressing the keys (3) and (4) increases or decreases the speed set which is shown on the display (1).
- Pressing the keys (5) and (6) increases or decreases the set time which is shown on the display (2).

WORKING FOR A NOT DEFINED TIME

When the display (2) shows «00» by pressing the key (6) the display shows: «--». If the centrifuge starts with this set time it only will stop manually.

- Start pressing the key (7). The centrifuge will stop once finished the set time.
During the working the display (1) shows the speed read by the tachometer.
The work can be stop pressing the key (10).

INSTANT START/STOP

- Do the previous operation as is explained on the section «Starting up». Once the speed and time of the working are set and different of 0, by pressing the key (7) the rotor starts to turn at the speed set. By stop pressing the centrifuge will stop.

EMERGENCY OPENING

In case of a cut of the power supply and it would be necessary opening the door, do the following:

- Remove the plastic stopper on the right lateral of the apparatus.
- Pull carefully from this stopper and the lid will open.

Do not use this emergency opening device when the machine is working.

ERRORS LIST:

Error 1: Eprom failure
 Error 3: Tachometer speed failure
 Error 4: Overcurrent
 Error 5: Low voltaje
 Error 6: High motor feedback voltaje.

9 Maintenance

Before of remove the cover to handle inside, unplug the the equipment from the socket.

The handle of the electronic circuit boards by unauthorized personnel may cause damages. Make sure to give the equipment to an authorised technical assistance.

9.1 Cleaning and decontamination

External cleaning

We recomend to use a cotton cloth wet with a water-alcohol disolution.

Inner cleaning and decontamination of sampler or products run out into the centrifuge vessel

The user has the responsibility to do the decontamination according to each dangerous product that could be run out inside the centrifuge vessel.

In case of samples splashed in the vessel clean it with a biological disinfectant.

9.2 Maintenance operation

The following parts make the user protection up againts impacts of unexpected parts thrown so a quarterly inspection of them are required:

- Lid, hinges, locks and enclosures: Inspect visually possible craking formations, deformations, rust stains, etc which might sign a possible dropping material resistance.
- Rubber legs: Check that the rubber legs keep the flexibility and grip on the working surface.

9.3 SPARE PARTS

The spare parts more commons:


Spare description	7002356
Electromagnetic lock	10040
Main switch	20017
Motor	24226
Tachometer sensor circuit board	29230
Control panel circuit board	29364
Power circuit board	29366
Mains filter	45011
Transformer 115-115/18V 1.2A	45048

10 Guarantee

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA,S.A.

Any manipulation of the apparatus by personnel not authorized by J.P. SELECTA,S.A. automatically cancels the guarantee.

11 Electric diagram

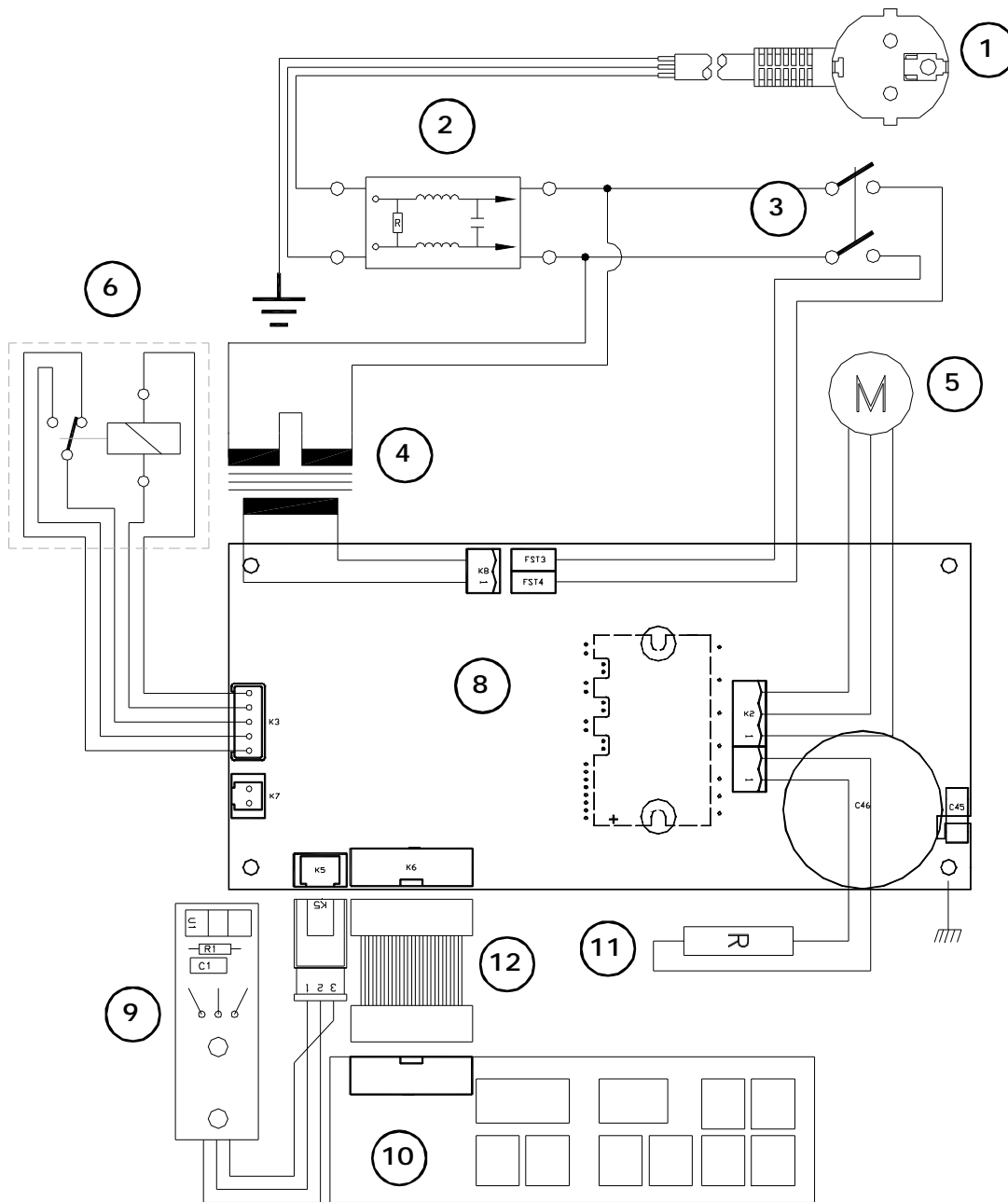Pos	Description	Code
1	Power supply cord	7140
2	Mains filter	45011
3	Main switch	29017
4	Transformer 115-115/18V 1.2 A	45048
5	Motor	24226
6	Electromagnetic lock	10040

Pos	Description	Code
8	Power circuit board	29366
9	Tachometer sensor circuit board	29230
10	Control panel circuit board	29364
11	Break resistance	
12	Board connecting wire	

DECLARATION DE CONFORMITE "CE" "EC" CONFORMITY DECLARATION

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PROGRAMME DE FABRICATION / MANUFACTURING PROGRAMME

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