pH-metro de bolsillo

REF - KDB003

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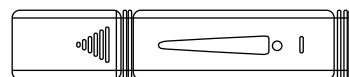
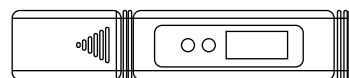
CODE- KDB003

1. SPECIFICATIONS

Range: 0.00 – 14.00 pH
Resolution: 0.01 pH
Accuracy: ± 0.05 pH
Without temperature compensation
Calibration in acid or basic pH range
Operating temperature: 0-60°C
Batteries: 2 x 1.5V (LR44 button cell)
Dimensions: 155 x 31 x 18 mm
Weight: 50 g

Note: Soak the electrode in distilled water for about 10 minutes before using it for the first time.

2. CALIBRATION

1. Remove the protective cap.
2. Rinse the electrode with distilled water and wipe it with filter paper.
3. Turn the pH meter on by pressing the ON/OFF key.
4. Immerse the electrode in pH 6.86 standard buffer solution. Wait for the reading to stabilize.
5. Press the CAL button for 5 sec. approx. (an increment of values will be displayed on the screen) and release the button.
6. The value of 6.86 will be fixed and then will flash.
7. Wait until it stops flashing.
8. Rinse the electrode with distilled water and wipe it with filter paper.
9. Put the electrode in the pH 4.00 solution.
10. Press the CAL button for 5 sec. approx. (an increment of values will be displayed on the screen) and release the button. The value of 6.86 will be reflected. Press the CAL button once to get the value of 4.00 (or 9.18 if you press the CAL button again to calibrate with it if necessary).
11. Wait until it stops flashing.
12. Turn off the instrument by pressing the ON/OFF key.
13. Clean the electrode with distilled water and wipe it with filter paper.
14. Place the protective cap.

Note: If your test solution is known to be above pH 7.00, calibrate using 6.86 and 9.18 standard buffer solutions.

3. OPERATION

1. Remove the protective cap.
2. Rinse the electrode with distilled water and wipe it with filter paper.
3. Turn the pH meter on by pressing the ON/OFF key.
4. Immerse the electrode in the sample to be tested. Wait for the reading to stabilize.
5. Turn off the instrument by pressing the ON/OFF key.
6. Clean the electrode with distilled water and wipe it with filter paper.
7. Place the protective cap.