AGITADOR DIGITAL 9-15 POSICIONES MULTIPOSITION MAGNETIC STIRRER, 9-15 POSITIONS AGITATEUR MAGNÉTIQUE MULTIPOSITIONS, 9-15 POSITIONS

Ref. | Code | Réf.

LBJ003 - Mod. 610/9

LBJ004 - Mod. 610/15

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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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TABLE OF CONTENTS

1. INTRODUCTION	4
2. TECHNICAL SPECIFICATIONS	4
3. OPERATION	5

1. INTRODUCTION

Nahita-Blue multi-position magnetic stirrers are characterized by a compact and robust structure, combined with a refined design and high versatility, making them reliable and easy-to-use instruments, ideal for scientific research environments.

This equipment is especially suitable for liquid mixing applications in the laboratory, allowing the simultaneous stirring of low-viscosity liquids or solid-liquid mixtures. It can be configured to operate either synchronously or asynchronously, depending on the user's needs.

2. TECHNICAL SPECIFICATIONS

Code	LBJ003	LBJ004
Model	610/9	610/15
Number of stations	9	15
Working plate material	PET anticorrosion plate	
Max stirring volume	9x500 mL	15x250 mL
Motor	Brushless motor	
Mixing direction	Counterclockwise direction	
RPM	50-1500 rpm	
Stirring bar size	30 mm	
Display	LCD Display	
Timing	1 min-23h59min	
Voltage	100-240V 50/60 Hz	
Permissible temperature and humidity	5-40°C / 80% HR	
Dimensions	435x330x85 mm	576x330x63 mm

3. OPERATION

■ Using steps:

- Connect the device to the power supply and press the power switch.
- Turn the left knob to select the station, then briefly press the right knob to adjust the desired parameters.
- Press and hold the right knob for 2 seconds to start/stop operation.
- Press and hold the left knob for 2 seconds to activate synchronization.

■ Instrument use environment:

- Indoor use only
- Altitude: ≤ 2000 m
- Working temperature range: $+5$ °C a $+40$ °C
- Humidity range: $\leq 80\%$
- No vibrations or airflow that could affect performance.
- No conductive dust, explosive gases, or corrosive gases in the surrounding air.
- Maintain a safety distance of 10 cm around the instrument while it is in operation.

■ Safety precautions:

- Do not unplug or plug in the power connector or toggle the power button with wet hands.
- Do not unplug the power plug while the instrument is powered on.
- Do not perform maintenance or cleaning while the instrument is powered on.
- Do not install the instrument on an uneven, swaying, or vibrating work surface.

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