**ESPECTROFOTOMETRO**

**ESPECTROFOTOMETER**

**REF: HJD007**

**MODELO: 4260/50**

# Español

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# English

<b>English</b> .....	<b>1</b>
<b>Safety</b> .....	<b>3</b>
<b>Package Contents</b> .....	<b>3</b>
<b>Unpacking</b> .....	<b>3</b>
<b>Installation</b> .....	<b>4</b>
3. <i>Environment Required</i> .....	4
4. <i>Install Spectrophotometer</i> .....	4
<b>Overview</b> .....	<b>4</b>
<b>Symbols</b> .....	<b>5</b>
<b>Main Specifications</b> .....	<b>5</b>
<b>Description of Appearance and Keys</b> .....	<b>6</b>
4. <i>Appearance</i> .....	6
5. <i>Keypad</i> .....	7
6. <i>Description of Keys</i> .....	8
<b>Functions</b> .....	<b>9</b>
<b>Getting Started</b> .....	<b>9</b>
<b>Important Guidelines</b> .....	<b>10</b>
<b>General Operating</b> .....	<b>11</b>
<b>Measuring</b> .....	<b>12</b>
8. <i>Photometry</i> .....	12
9. <i>Quantitation</i> .....	14
10. <i>Wavelength Scan</i> .....	17
11. <i>Kinetics</i> .....	18
12. <i>DNA/Protein Mode</i> .....	20
13. <i>Multi Wavelength Mode</i> .....	21

14. System Utility.....	22
Wavelength Reset .....	22
Printer Setup.....	23
Lamp Service.....	23
Clock Setup.....	25
Refresh Dark Current.....	25
Connect to PC.....	26
Beeper on/off.....	26
Refresh System Baseline.....	26
Deleted Entire Saved Files.....	26
Restore Default Settings.....	26
<b>Troubleshooting .....</b>	<b>27</b>
<b>Repair and Maintenance .....</b>	<b>27</b>
3. Daily Maintain.....	27
4. Spare Parts Replacement.....	28
<b>Warranty .....</b>	<b>32</b>
<b>Equipment Disposal.....</b>	<b>33</b>

# Safety

Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

We recommend against the use of Spectrophotometer.

- Do not open the device.
- Disconnect the device from the mains supply before carrying out maintenance work or changing the fuses.
- The inside of the device is a high-voltage area Danger!
- Do not use the device if it is damaged, especially if the main power cable is in any way damaged or defective.
- Repairs may only be carried out by the service technicians from us and authorized contractual partners.
- The device must be connected to a power outlet that has a protective ground connection.
- If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

- Do not allow any liquid to enter into the device.
- Do not operate the device in a hazardous location or potentially explosive environment.

# Package Contents

Description	Quantity
Spectrophotometer	1PC
10mm Glass Cuvette	4PCS
10mm Quartz Cuvette	2PCS
Power Cord	1PC
User's Manual	1PC
Dust Cover	1PC

# Unpacking

Open the package, according to carefully check the packaging packing list items, if found inside the packaging are missing or damaged items

please contact us and authorized contractual partners.

## Installation

### 3. Environment Required

To ensure the best performance, the following conditions are required:

- The best working temperature range is 16—35°C and the humidity is 45—80%.
- Keep it as far as possible away from the strong magnetic or electrical fields or any electrical device that may generate high-frequency fields.
- Set the unit up in an area that is free of dust, corrosive gases and strong vibrations.
- Remove any obstructions or materials that could hinder the flow of air under and around the instrument.
- The power requirement is  $110\pm 11\text{V}/60\pm 1\text{Hz}$  or  $220\pm 22\text{V}/50\pm 1\text{Hz}$ .
- Use the appropriate power cord and plug into a grounded outlet.
- If the local voltage is not stable, a voltage regulator is required.
- Be away from direct sunlight.

### 4. Install Spectrophotometer

#### Placement

Place the instrument on the stable table carefully.

#### Install Printer (Printer is Optional Accessories)

Check to confirm instrument power switch is turned off, connect the printer's data cable to the instrument's parallel port.

#### Link the Power Cord

Check to confirm instrument power switch is turned off, the power cord plug into two separate power interface and power supply socket apparatus.

## Overview

Spectrophotometer is an electrical measure instrument which is widely used in the laboratories.

- Use Frequency: Intermittence
- Excessive Voltage(Current): No
- Pollution Class: Class 1

# Symbols

The following chart is an illustrated glossary of the symbols that are used in this manual.

Caution, Danger!

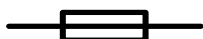
Caution, High Voltage!

Caution, Hot!

Ground

Fuse

Recycle, this instrument will be called back by the appointed Electrical Treatment Department or by the original Manufacturer when wasted.

# Main Specifications

· <b>Optical System</b>	Double beam
· <b>Wavelength Range</b>	190—1100nm
· <b>Wavelength Accuracy</b>	±0.3nm
· <b>Wavelength Repeatability</b>	0.2nm
· <b>Photometric Range</b>	-0.3—3A, 0—200%T
· <b>Photometric Accuracy</b>	±0.5%T
· <b>Photometric Repeatability</b>	0.3%T
· <b>Spectral Bandwidth</b>	0.5/1/2/4/5nm
· <b>Stray Light</b>	0.05%T@220nm&360nm
· <b>Stability</b>	±0.001A/h@500nm
· <b>Work Mode</b>	Photometry, Quantitation, Kinetics,

- |                            |   |
|----------------------------|---|
|                            | Wavelength Scan, DNA/Protein,<br>Multi Wavelength |
| · <b>Interface</b>         | USB, Parallel (printer)                           |
| · <b>Power Requirement</b> | AC 110/220V, 50/60Hz                              |
| · <b>Dimensions</b>        | 600x450x260                                       |
| · <b>Weight</b>            | 20kg  |
| · <b>Work Environment</b>  | 15—35°C, 15—70% relative humidity                 |
| · <b>Store Environment</b> | -10 — 50 °C , 15 — 70% relative<br>humidity       |

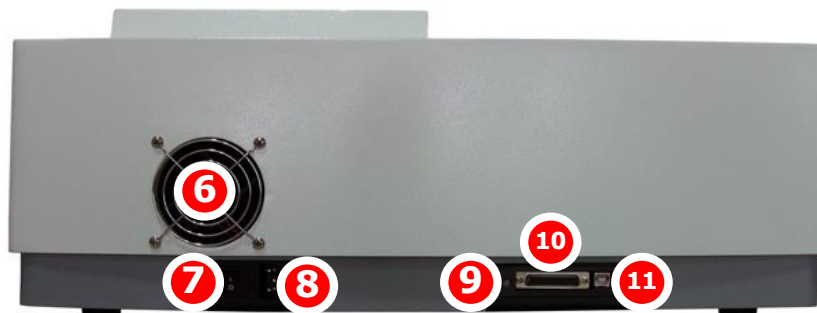
## Description of Appearance and Keys

### 4. Appearance

Front View


Back View


- |                                     |                 |
|-------------------------------------|-----------------|
| 1 LCD Display                       | 2 Keypad        |
| 3 Lid of Sample Room                | 4 Sample Room   |
| 5 USB Port (Type A, USB Memory)     | 6 Cover of Fan  |
| 7 Power Switch                      | 8 Power Socket  |
| 9 LCD Contrast Adjust               | 10 Printer port |
| 11 USB port (Type B, Communication) |                 |


## 5. Keypad
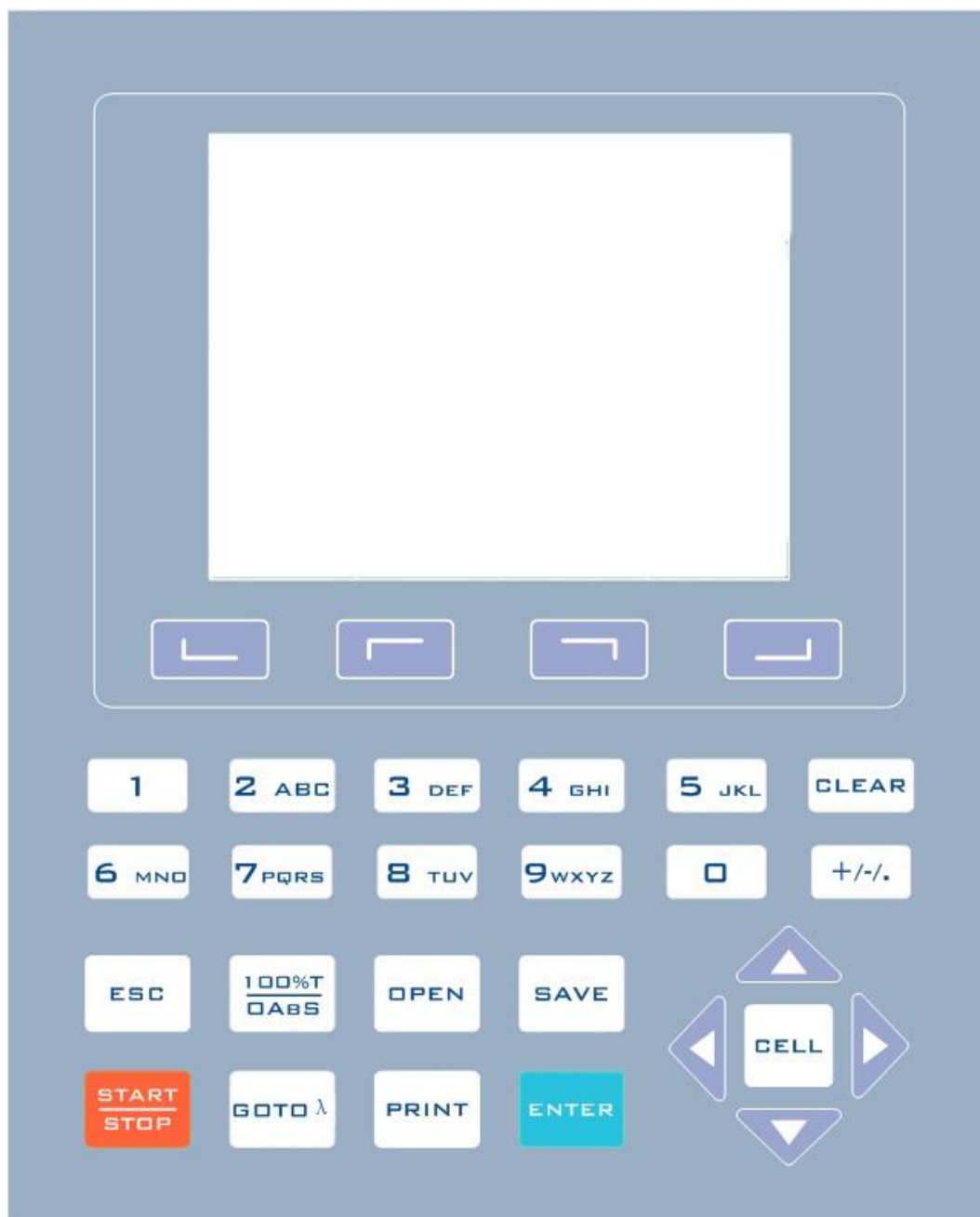
## 6. Description of Keys

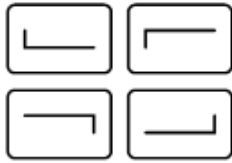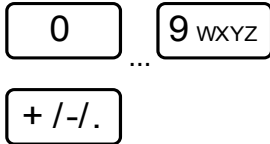**Function Keys:** Functions on-screen prompts


**Numeric Keys:** Enter numbers and letters

**CLEAR Key:** Delete the input value or stored data

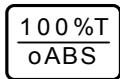**ESC Key:** Return to previous Interface

**100%T/0Abs Key:** Blank

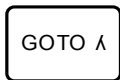
**OPEN Key:** Open files stored in internal memory

**SAVE Key:** Save files to internal memory

**START/STOP Key:** Start/Stop testing

**GOTO λ Key:** Set wavelength

**PRINT Key:** Print measuring result

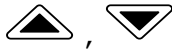
**ENTER Key:** Confirm operation

**CELL Key:** Select/Deselect Auto-cell Holder

**RIGHT, LEFT Keys:** Search peak/valley and set X scale

**UP, DOWN Keys:** Scroll menu/data and set Y scale

## Functions

### Photometry

Display results as Abs, %T or Concentration.

### Quantitation

We provide 2 methods to establish a Standard Curve:

- 2 methods to set up the regression curve.  
Up to 10 standard samples to calibrate the regression curve.
- 3 methods for curve fit.  
Linear fit, Quadratic fit and Cubic fit.

### Wavelength Scan

- Allow user to set scan step (0.1, 0.2, 0.5, 1.0 and 5.0nm).
- Spectrum display mode can be changed (Wavelength-%Transmittance or Wavelength-Absorbance).

### Kinetics

- Allow user to set scan Interval (0.5, 1.0, 2.0, 5.0, 10, 30 and 60s).
- Spectrum display mode can be changed (Time-%Transmittance or Time-Absorbance).

### DNA/Protein Measurement

- Wavelength points and ratios can be set up.
- Results will be grouped into a table format automatically.

### Multi-wavelength



- Up to 20 wavelength points can be set up.
- Results will be grouped into a table format automatically.

## Getting Started


The following chart describes the basic operation of the instrument.
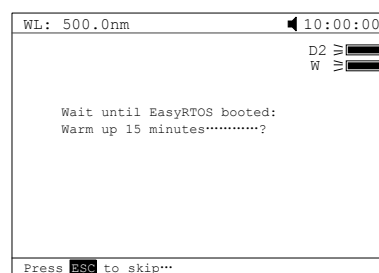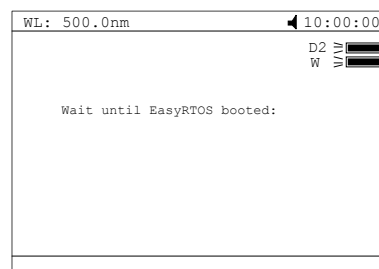
## Turn On and Self-check


Switch on the power. Then the instrument begins to self-check and 15 minutes' warm up. Self-check includes the following steps: Turn on lamps → Check Sensor → Initialize AD → System position → Get Dark Current → Warm up.

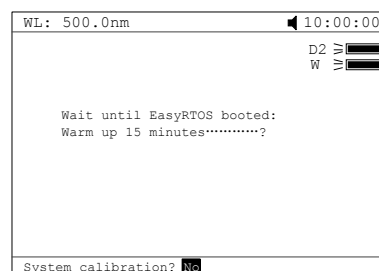Finished self-check, "System calibration" interface displays, press the keys ,  to select to do calibration or not.

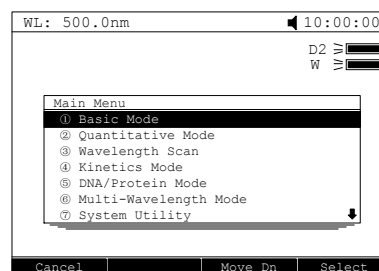After system do or skip calibration, instrument displays Main Interface.

## Important Guidelines

- Reagents and dilution buffers can cause cauterization and other damage to health.
- Samples (nucleic acids, proteins, bacteria cultures) can be infectious and cause serious damage to health.
- During sample preparation, measuring procedures and maintenance and cleaning work, observe all local laboratory safety precautions (e.g.

wear protective clothing and gloves, use of disinfectant) regarding the handling of sample material.

- Dispose of measuring solutions and cleaning and disinfectant materials in accordance with the relevant local laboratory regulations.

## General Operating

### Select Application

Main menu, press numeric keys or use the key ▲, ▼ to choose corresponding menu, then press  to enter into.

### Set Wavelength

Press  to set wavelength, use numeric keys to input the values, press  to confirm and go to the point you set, then do blank automatically.

### Set Parameters

Press function key to set parameters, press ▲, ▼ to choose or input the values by numeric keys, press  to enter into, press  to return.

### Set Auto-cell Holder

Press  to active the auto cell holder and press the numeric key (1-8) to make corresponding cell position at the light path. Press  again to inactive the auto cell holder.

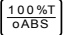
### Delete the Input Value

Press  to delete a character, press  to delete all the characters.

### Delete the Test Results and Stored data

Press  to delete the test result or stored data.

### Blank

Put the Reference in the light path, press  to do blank.

### Measure Samples

Put the samples in the light path, press  to measure.

### Print the Test Results

Press  to print the test results.

### Store the Test Results

Press  to store the test results, input the file name by numeric keys and press  to save. If the USB Disk was installed, the data will save in the USB Disk, or the data will save in the internal memory.

### Load the Stored File

In the test interface, press  to go into file selecting interface, press ,  to choose the file you want, press  to open. If the USB Disk was installed, the data will open from the USB Disk, or the data will open from the internal memory.

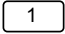
### Select Bandwidth


Main menu, press , press ,  to choose the Bandwidth, press .

## Measuring

### 8. Photometry

#### Step 1. Start Photometry

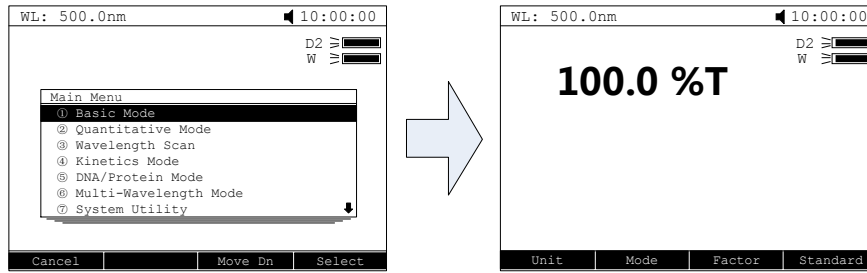Main menu, press numeric key  or ,  to choose "Basic Mode", then press .

## Step 2. Set Photometric Mode

Press  to set photometric mode. Press ,  to choose "Abs.", "T%" or "Conc./Factor" and press  to confirm. If users choose "Abs." or "T%", please go to step 5 directly.

## Step 3. Set Concentration Unit

Press  to set concentration unit. Press ,  to choose unit followed with  pressed to confirm. You can also choose "other" to input the unit.

## Step 4. Set "Factor" or "Standard"

Two methods are under your choice:

### Method 1: Input Factor F

Press  to set F. Input the value of F by numeric keypad, press  to confirm. Then the F value would display on the screen.

### Method 2: Use Standard Sample

Put the reference sample in the light path and calibrate 100%T/0Abs;

Put the standard sample in the light path, press  to start the mark. Input the concentration value of the standard and press  to confirm, then it displays on the screen.

## Step 5. Set Wavelength

Press  to set wavelength, input the value by the numeric keypad followed with  pressed to confirm.

## Step 6. Measure samples

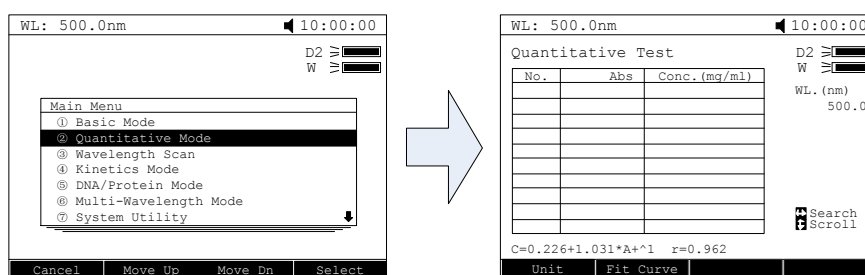
Put the sample to be measured in the main light path and put Reference in the reference light path, then the result displays on the screen automatically.

# 9. Quantitation

## Step 1. Start Quantitation

Main menu, press  or press ,  to choose "Quantitative Mode"

followed with  pressed to confirm.


## Step 2. Set Unit

Press  to set concentration unit, press ,  to choose and press  to confirm.

## Step 3. Establish Standard Curve or load the stored curve

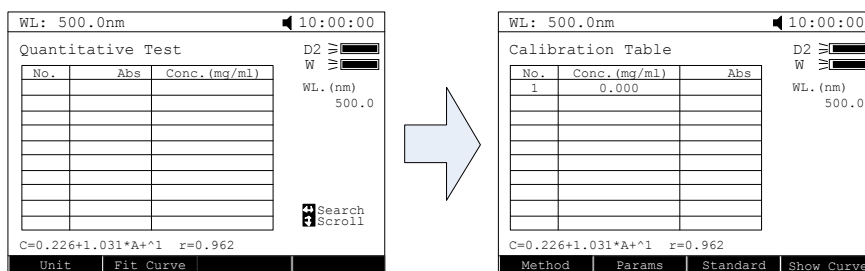Press  to go into set up interface, 2 methods are under your choice.

### Establish Standard Curve:

#### Method 1: Input Regression Equation

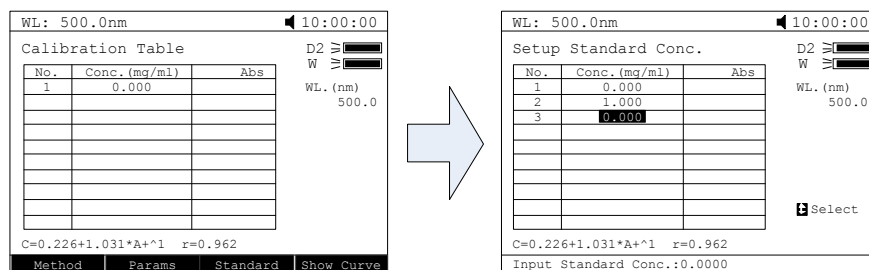
4) **Set Fit Curve Method.** Press  to set Fit method, use ,

to choose the method and press  to confirm.

- 5) **Set Wavelength.** Press  to set wavelength. Use ,  to choose measure method, then press  to confirm. Input the wavelength value you need and press  to confirm.
- 6) **Input the Factor of the Regression Equation.** Press  and input the factors, press  to confirm.


### Method 2: Use Standard Samples

- 5) **Set Fit Method.** Press  to set fit method, press ,  to choose fit method, then press  to confirm.
- 6) **Set Wavelength.** Press  to go into wavelength setting interface, press ,  to select measure method and press  to confirm. Input the value of the wavelength and press  to confirm.
- 7) **Setup Standard Samples.** Press  to setup standard, input the concentrations of corresponding standard samples according the indication and press  to confirm. Users can use ,  to choose the value you just input and press  to delete, then input a new value, press  to confirm. Press  to cancel after all the input.


- 8) **Calibrate Standard Samples.** Put the corresponding standard samples in the main light path and put Reference in the reference light path as the screen indicates and press  to

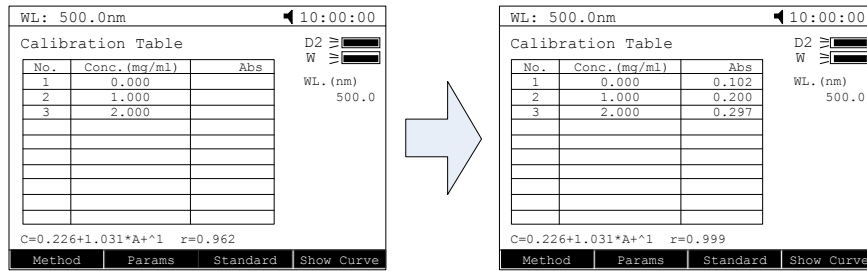measure. Then the Abs. value would appear in the corresponding table.


### Load the Stored Curves

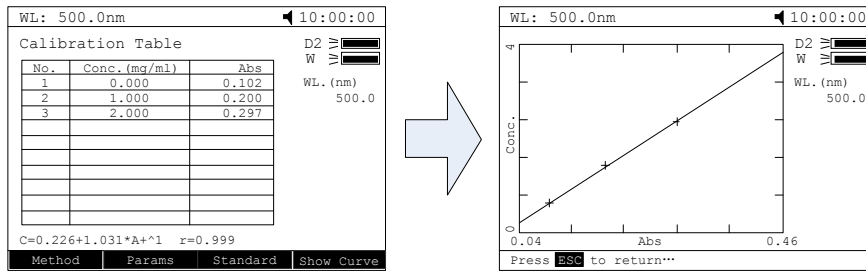In the "Calibration Table" interface, press **OPEN** to go into files select interface. Use **▲**, **▼** to select the curve you need and press **ENTER** to load. Users can press **□** to view the curve.

### Step 4. Return the sample measurement interface

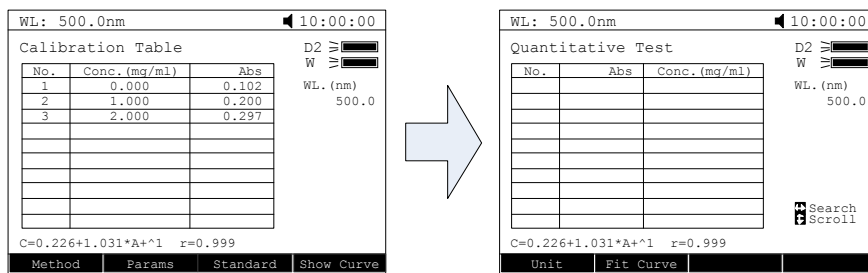In the "Calibration Table" interface, press **ESC** to return the Quantitative Test interface.

### Step 5. Measure Samples

Place the sample to be tested in the light path, press **START STOP** to measure. Then the test result will display in the data sheet. Repeat this step to finish measuring all the samples.

WL: 500.0nm 10:00:00

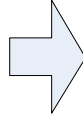
Calibration Table

No.	Conc. (mg/ml)	Abs
1	0.000	0.102
2	1.000	0.200
3	2.000	0.297

D2 W W  
W W  
WL. (nm)  
500.0

$C=0.226+1.031*A+1$   $r=0.999$

Method Params Standard Show Curve

WL: 500.0nm 10:00:00

Quantitative Test

No.	Abs	Conc. (mg/ml)
1	0.158	1.601
2	0.155	1.598
* 3	0.160	1.614

D2 W W  
W W  
WL. (nm)  
500.0

Search  
Scroll

$C=0.226+1.031*A+1$   $r=0.999$

Unit Fit Curve

## 10. Wavelength Scan

### Step 1. Enter into Wavelength Scan

Main menu, press numeric key **3 DEF** or , to choose "Wavelength Scan" and press **ENTER** to enter.

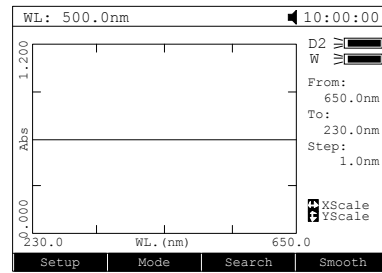
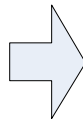
WL: 500.0nm 10:00:00

D2 W W  
W W

Main Menu

- ① Basic Mode
- ② Quantitative Mode
- ③ **Wavelength Scan**
- ④ Kinetics Mode
- ⑤ DNA/Protein Mode
- ⑥ Multi-Wavelength Mode
- ⑦ System Utility

Cancel Move Up Move Dn Select

### Step 2. Parameters Setup

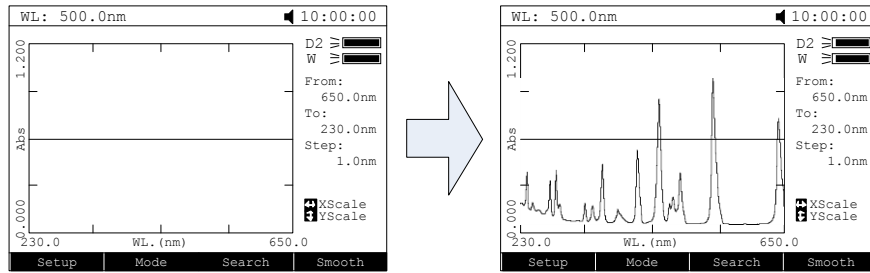
Press to set parameters, set "Scan From" , "scan to", "scan step" and "scan speed", press **ENTER** to confirm.

### Step 3. Set Photometric Mode



Press to set photometric mode, choose "T%", "Abs." or "E" and press **ENTER** to confirm.


### Step 4. Scan Samples

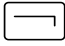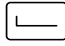Put the sample to be measured in the main light path and put Reference in the reference light path, press **START STOP** to scan the sample, press **ESC** to cancel.


### Step 5. Search Peaks

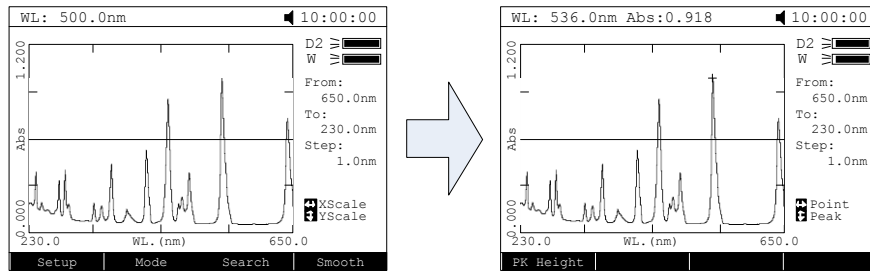After scanned, press  to go into peak search mode. Press  to set peak height, input the peak height and press  to confirm. Press ,  to display the value of every wavelength point. Press ,  to display the value of every peak.


### Step 6. Smooth the Curve

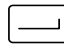After scanned, if there are many burrs on it, press  to smooth the curve.


## 11. Kinetics
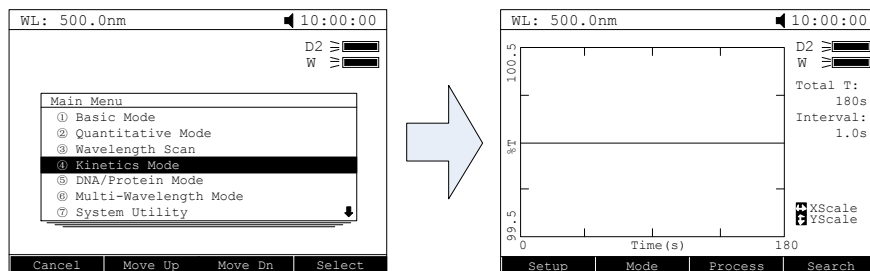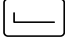
### Step 1. Enter into Kinetics


Main menu, press  or ,  to select "Kinetics Mode" and press  to confirm.

### Step 2. Setup Parameters

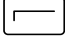Press  to set parameters, input the corresponding values of "Total Time", "Delay Time" and "Time Intervals" according the screen indicates. Press  to confirm.

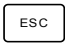
### Step 3. Set Photometric Mode


Press  to set photometric mode, choose "T%" or "Abs." and press  to confirm.

### Step 4. Set Wavelength

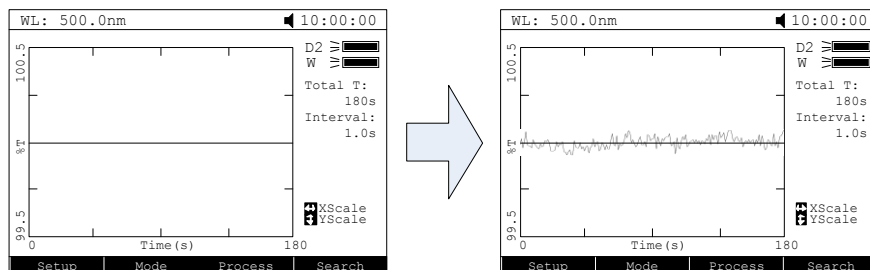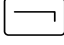
Press  to set wavelength, input the value of the wavelength by numeric keypad and press  to confirm.


### Step 5. Measure Samples

Put the sample to be measured in the main light path and put Reference in the reference light path, press  to begin the test, press  to cancel.



### Step 6. Calculate Response Rate

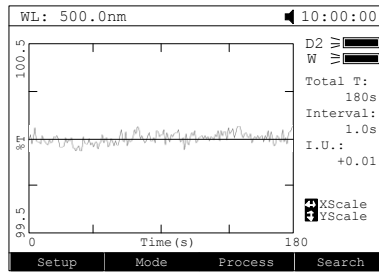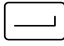After time scanned, if users want to calculate the response rate of some period, you can press  to go into "Process" interface. Input the values of "Begin Time", "End Time" and "Factor" separately and press  to confirm. Then the value of "I.U." would display on the screen.


### Step 7. Search Peaks

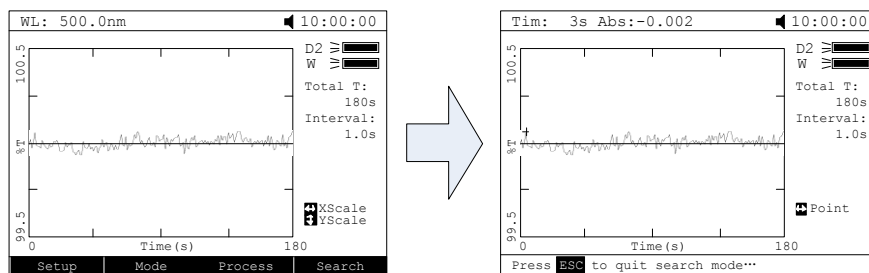
After scan finished, press  to go into search mode. Press ,  to search the value of every point.


## 12. DNA/Protein Mode
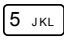
### Step 1. Enter into DNA/Protein Mode

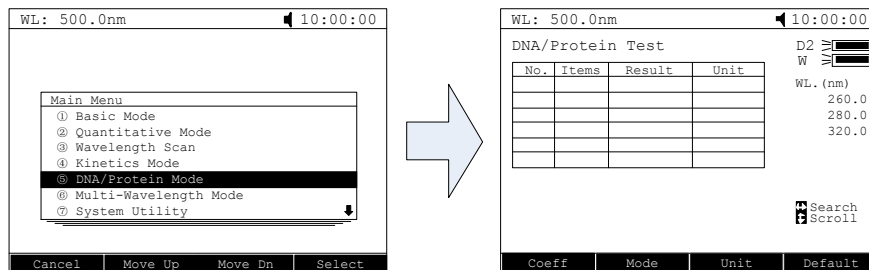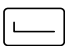Main menu, press  or ,  to choose "DNA/Protein Mode" and press  to confirm.


### Step 2. Setup Parameters

Press  to set coefficients, input all the values of f1 to f4 by numeric keypad according the indication and press  to confirm.

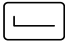
### Step 3. Choose Measure Method

## Step 2. Setup Wavelength

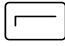
Press  to go into wavelength setting interface, input all the wavelength value one by one by the numeric keypad. Press  to confirm, press  to return.

## Step 3. Set Photometric Mode

Press  to set the photometric mode, use ,  to select "Abs." or "T%" mode, press  to confirm.

## Step 4. Measure Samples

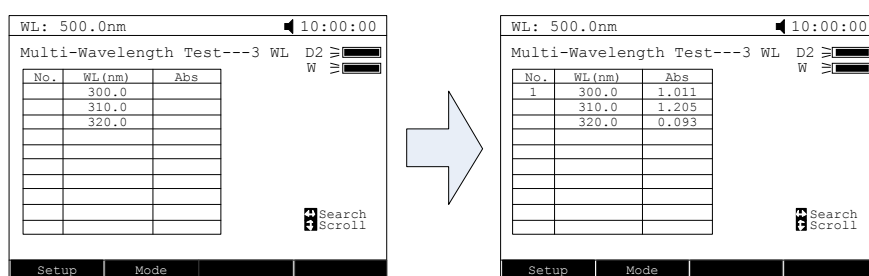Put the sample to be measured in the main light path and put Reference in the reference light path, press  to measure, the test result will display in the data table.


The diagram shows two screenshots of the 'Multi-Wavelength Test' interface. The left screenshot shows a table with headers 'No.', 'WL (nm)', and 'Abs' and three rows of empty data. The right screenshot shows the same table with the following data:

No.	WL (nm)	Abs
1	300.0	1.011
	310.0	1.205
	320.0	0.093

# 14. System Utility

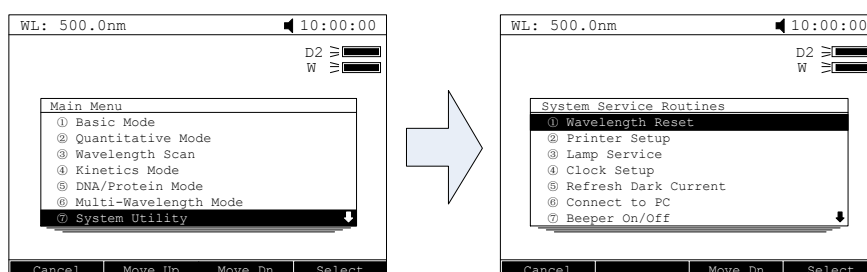Main menu, press  or use ,  to select "System Utility" and press  to confirm.


The diagram shows two screenshots of the system utility menu. The left screenshot shows the 'Main Menu' with 'System Utility' selected. The right screenshot shows the 'System Service Routines' menu with 'Wavelength Reset' selected.

Main Menu	
①	Basic Mode
②	Quantitative Mode
③	Wavelength Scan
④	Kinetics Mode
⑤	DNA/Protein Mode
⑥	Multi-Wavelength Mode
⑦	System Utility

System Service Routines	
①	Wavelength Reset
②	Printer Setup
③	Lamp Service
④	Clock Setup
⑤	Refresh Dark Current
⑥	Connect to PC
⑦	Beeper On/Off


## Wavelength Reset

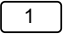Press  or use ,  to choose "Wavelength Reset" then press  to begin the calibration. During the course, opening the lid of the

compartment is prohibited.

### Printer Setup

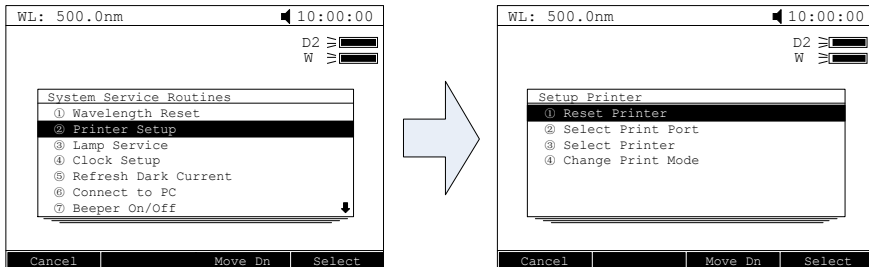Press **2 ABC** or **▲**, **▼** to select "Printer Setup" then press **ENTER** to confirm.

### Reset Printer

Press **1** or use **▲**, **▼** to choose "Reset Printer" and press **ENTER** to confirm. Then the printer will resume the initial condition.

### Select print port

Press **2 ABC** or use **▲**, **▼** to choose "Select print port" and press **ENTER** to confirm. Use **▲**, **▼** to choose "LPT" or "Comm." and then press **ENTER** to confirm.

### Select Printer

Press **3 DEF** or use **▲**, **▼** to choose "Select printer" and press **ENTER** to confirm. Use **▲**, **▼** to select the printer's model and then press **ENTER** to confirm.

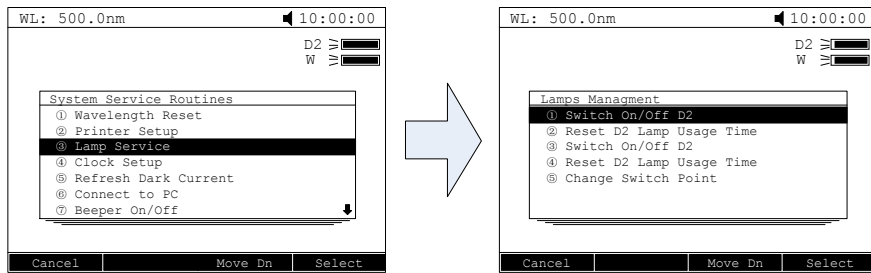
### Change Print Mode

Press **4 GHI** or use **▲**, **▼** to choose "Change print mode" and press **ENTER** to confirm. Two modes are under your choice, one is "Print Data Sheet", the other is "print the display interface".

### Lamp Service

Press **3 DEF** or use **▲**, **▼** to choose "Lamp Service", press **ENTER** to go

into "lamps management" interface.

### Switch On/Off D2 lamp

Press  or ,  to choose "Switch On/Off D2 lamp" then press  to switch on or switch off the D2 lamp.

### Reset the usage time of D2 lamp

Press  or use ,  to choose "Reset D2 lamp usage time" and press . Then users can find the D2 lamp's usage time and the system will ask if you are sure to reset the usage time. Press ,  to choose "Yes" and then press  to confirm. The system will record the usage time from zero.

### Switch On/Off W lamp

Press  or use ,  to choose "Reset W lamp usage time" and press . Then users can find the W lamp's usage time and the system will ask you if are sure to reset the usage time.

### Reset the W lamp usage time

Press  or use ,  to choose "Reset W lamp usage time" and press . Then users can find the W lamp's usage time and the system will ask if you are sure to reset the usage time. Press ,  to choose "Yes" and then press  to confirm. The system will record the usage time from zero.

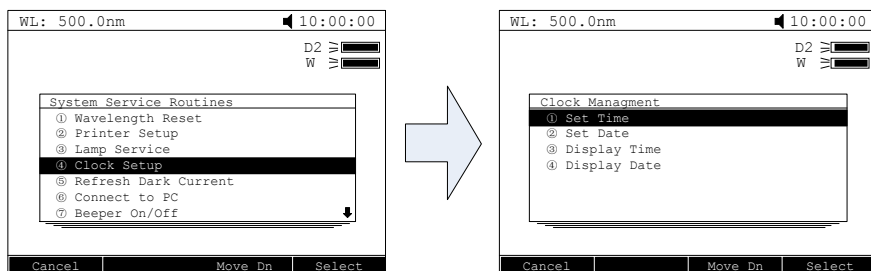
### Change Switch point

Press  or use ,  to choose "Change Switch point" and

press  to confirm. Input the wavelength point value(325–375nm)and press  to confirm.

### Clock Setup

Press  or use ,  to choose "Clock Setup", then press  to go into "Clock Management" Interface.

### Set Time

Press  or ,  to choose "Set Time" and press  to confirm. Input the time (Hour, Minute, Second) by numeric keypad, press  to confirm and return automatically.

### Set Date

Press  or use ,  to choose "Set Date" and press  to confirm. Input the Date(Year, Month, Day) by numeric keypad, press  to confirm and return automatically.

### Display Time

Press  or use ,  to choose "Display Time" and press  to confirm. Then the time will display on the Right Top corner.

### Display Date

Press  or use ,  to choose "Display Date" and press  to confirm. Then the Date will display on the Right Top corner.

### Refresh Dark Current

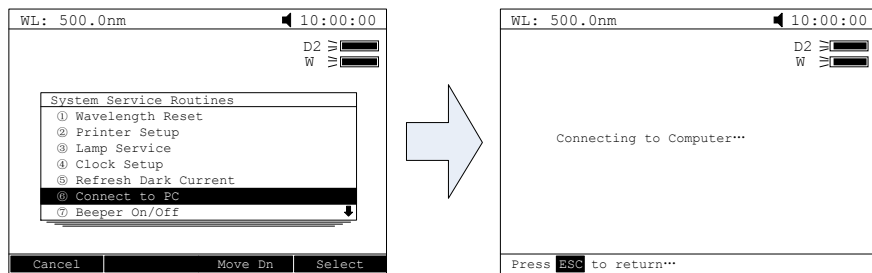
Press  or use ,  to choose "Refresh Dark Current" and press

**ENTER** to confirm. Then the system begins to refresh Dark Current.

**Note: During the course, open the lid of the compartment is prohibited.**

### Connect to PC

Press **6 MNO** or use **▲**, **▼** to choose "Connect to PC", and press **ENTER** to go into waiting interface. When the instrument was connected to PC, it displays "Controlled by PC".

### Beeper on/off

Press **7 PQRS** or use **▲**, **▼** to choose "Beeper on/off" and press **ENTER** to switch on or switch off the beeper.

### Refresh System Baseline

Press **8 TUV** or use **▲**, **▼** to choose "Refresh System Baseline" and press **ENTER** to confirm. Then the system will scan the baseline. During this course, opening the lid of the compartment is prohibited.

### Deleted Entire Saved Files

Use **▲**, **▼** to choose "Deleted Entire Saved Files" and press **ENTER** to confirm. Then the system ask you "Delete entire files, are you sure?", use **▲**, **▼** to choose "Yes" , then all the files in the RAM will be deleted.

### Restore Default Settings.

Use **▲**, **▼** to choose "Restore Default Settings" and press **ENTER** to Confirm, then the system will restore the initialization.

# Troubleshooting

Review the information in the table below to troubleshoot operating problems.

Problem	Cause	Solution
Power on, no response	Power cord connection is not reliable	Improve connectivity
	Fuse burning	Replace fuse
Measurement uncertainty	Warm up is not enough	Warm up more time
	Glass cuvettes used in UV Range	Use quartz cuvettes
	Sample is not Stable	Improve the sample
	The concentration of sample is too high	Diluted sample
	Power Supply Voltage Low or not Stable	Improve the Power Supply
	Lamp damage or lamp life maturity	Replace lamp
Dark Current Error when self-check	The lid of the compartment is open during self-check	Close the lid, restart
System Calibrate Failed	Something block the Light path	Remove it, calibrate again
Power on, back light is OK, but nothing display on the screen or display is not clear	Display Contrast problem	Adjust the contrast potentiometer
Measurements inaccurate	Cuvettes were contaminated	Clean cuvettes
	Samples were contaminated	Improve samples
	Worse matching of the cuvettes	Improve the matching of the cuvettes
	Dark current error	Resample dark current

## Repair and Maintenance

### 3. Daily Maintain

#### Check the Compartment

After measurement, the cuvettes with sample solutions should be taken

out of the compartment in time. Or the volatilization of the solution would make the mirror go moldy. Users must pay more attention to the corrosive sample and liquid easy to volatilize. Any solution remains in the compartment should be wiped off immediately.

#### **Surface Clean**

The cover of the instrument is with paint. Please use wet towel to wipe off the drips on the surface immediately. Organic solution is forbidden to be used to clean the cover. Please wipe off the dirt on the cover timely.

#### **Clean the Cuvettes**

After every test or after a solution change, the cuvettes should be cleaned carefully, or the remains on the surface would cause measuring error.

## **4. Spare Parts Replacement**

### **Replace Fuse**

**Danger! Be sure to switch off the power and unplug the socket before replacement!**

#### **Step 1. Tools preparation**

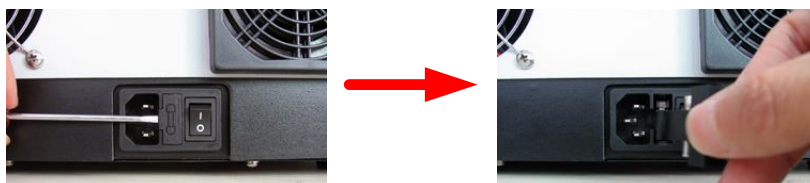
Prepare a 3×75 Flat Blade screwdriver.

#### **Step 2. Switch Off the power supply**

Switch off the power supply, and unplug the socket.


#### **Step 3. Take out the Fuse Seat**

Take out the fuse seat by the screwdriver.

#### **Step 4. Replace a new fuse**

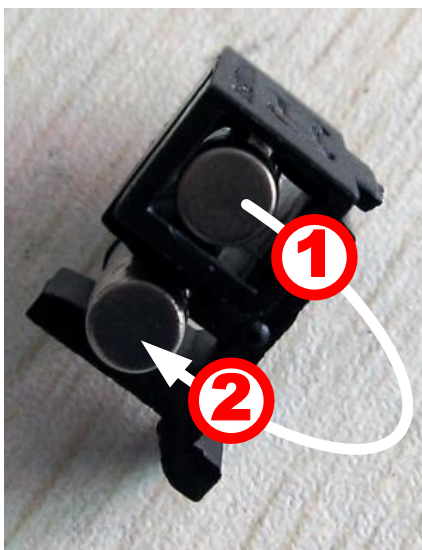Pick out the spare fuse (3.15A/250V) and replace it to the working position.

**Step 5. Reset the fuse seat**

Replace the fuse seat in the power socket.

**Step 6. Switch on the power**

Plug the socket and switch on the power.

**Replace Lamps**

**Hot! Wait 20 minutes before open the lamp chamber after power off to avoid scald!**

**Step 1. Tools preparation**

Prepare a 6×150mm Cross Blade screwdriver and a pair of glove.

**Step 2. Power Off**

Switch off the power supply and unplug the socket.

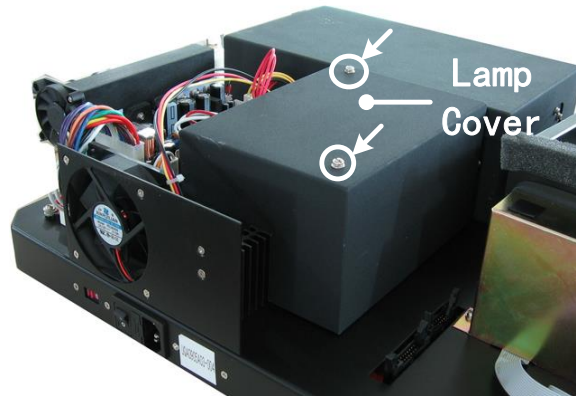
**Step 3. Open the cover**

Unscrew the 4 screws indicated(Each side with 2 screws) and remove the cover.


**Step 4. Open the cover of the light chamber**

Unscrew the 2 screws on the light chamber cover and remove it.


**Step 5. Replace the D2 lamp**

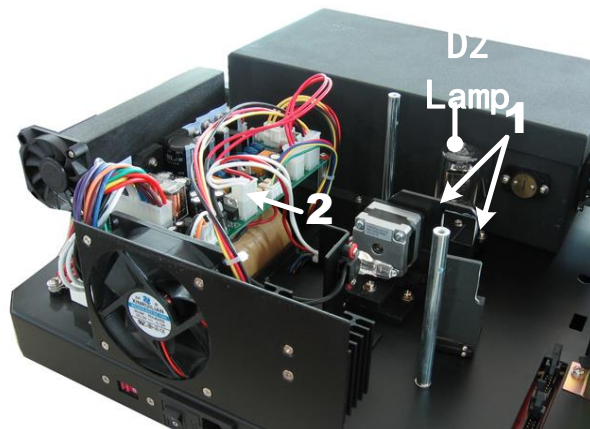Unscrew the 2 screws on the D2 Flange (No.1), unplug the connector in the Power Board (No. 2) and remove the D2 lamp. Draw on the cotton glove and replace a new lamp. Fix the 2 screws and plug the connector again.


**Step 6. Replace W lamp**

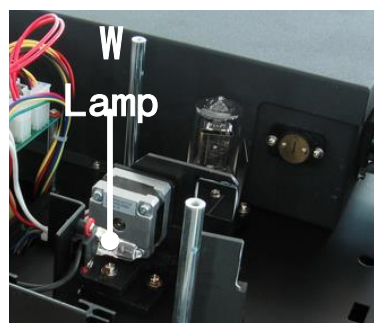Pull out the defected W lamp and draw on the cotton glove. Insert the new W lamp as deep as possible on the lamp seat.

**Step 7. Adjust the position of the W lamp**

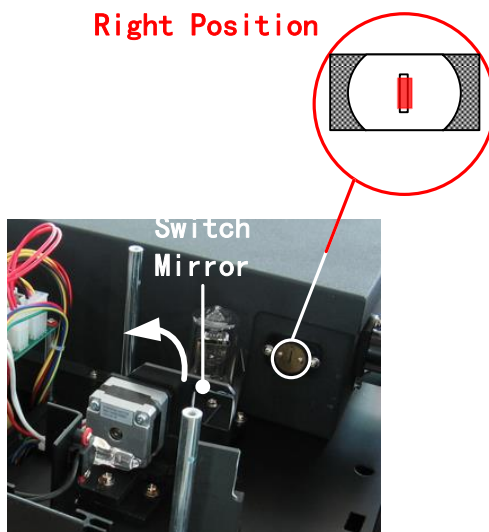Switch on the power(the Switch Mirror should be placed to the position as indicates). Observe the entrance facular, adjust the W Lamp until the light is in the center of the entrance hole.

**Step 8. Finish**

Reset the cover of the light chamber and fix the screws. Reset the cover of the instrument and fix the screws. Recover the Pole in the compartment, then the course finished.

**Replace the Battery**

**Danger! Be sure to switch off the power and unplug the socket before replacement!**

**Step 1. Prepare the tools**

Prepare a 6×150mm Cross Blade Screwdriver.

**Step 2. Switch off the power supply**

Switch off the power supply and unplug the socket.

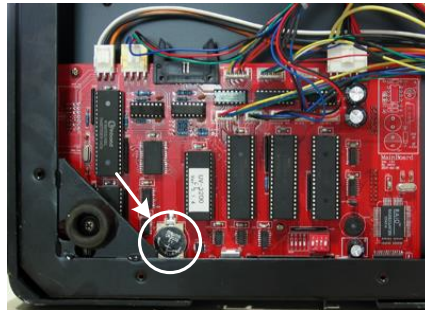
**Step 3. Open the Bottom cover plate**

Unscrew all the screws indicated then remove the bottom plate.

#### **Step 4. Replace the Battery**

Pick out the old battery and replace a new one.

#### **Step 5. Finish**

Recover the bottom plate and fix all the screws, then the course finishes.

## **Warranty**

**We** warrant that this product will be free from defects in material and workmanship for a period of one (1) year from date of purchase. If a defect is present, we will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

# Equipment Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you!