# ESPECTROFOTÓMETROS SPECTROPHOTOMETERS SPECTROPHOTOMÈTRE

## Modelos / Models / MODÈLES 4211/20 Y 4211/50

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Zuzi spectrophotometers models 4211/20 and 4211/50. We highly recommend looking after this equipment according to what is stated in this manual.

Zuzi develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the Zuzi spectrophotometer model 4211/20 and 4211/50, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the spectrophotometer avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the spectrophotometer to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the spectrophotometer.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Zuzi Technical Assistance Department by email to [asistencia@auxilab.es](mailto:asistencia@auxilab.es).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Zuzi Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

### INDEX OF LANGUAGES

Spanish .....	2-16
English.....	17-30
French.....	31-45

### INDEX OF CONTAINS

1. USES OF THE INSTRUMENT .....	18
2. DESCRIPTION .....	18
3. TECHNICAL SPECIFICATIONS .....	19
4. INSTALLATION / SETTING UP .....	19
5. MAINTENANCE AND CLEANING .....	29
6. TROUBLESHOOTING .....	30
ANNEX I: CE CERTIFICATE .....	46

## 1. USES OF THE INSTRUMENT

Zuzi spectrophotometers are versatile instruments that perform precise quantitative and qualitative analysis in the visible and ultraviolet spectral range (depending on model). They are very useful in conventional laboratories and in a wide range of fields such as petrochemical and mining industry, agricultural engineering, environmental preservation, biochemistry, hospitals and research laboratories.

## 2. DESCRIPTION

Figure 1

- 1.1 LCD display and control panel
- 1.2 Sample compartment
- 1.3 Cell changer manual knob
- 1.4 Power ON/OFF
- 1.5 Fuse holder
- 1.6 Socket
- 1.7 USB and parallel ports
- 1.8 Manual power switch 110/220V (see detail)


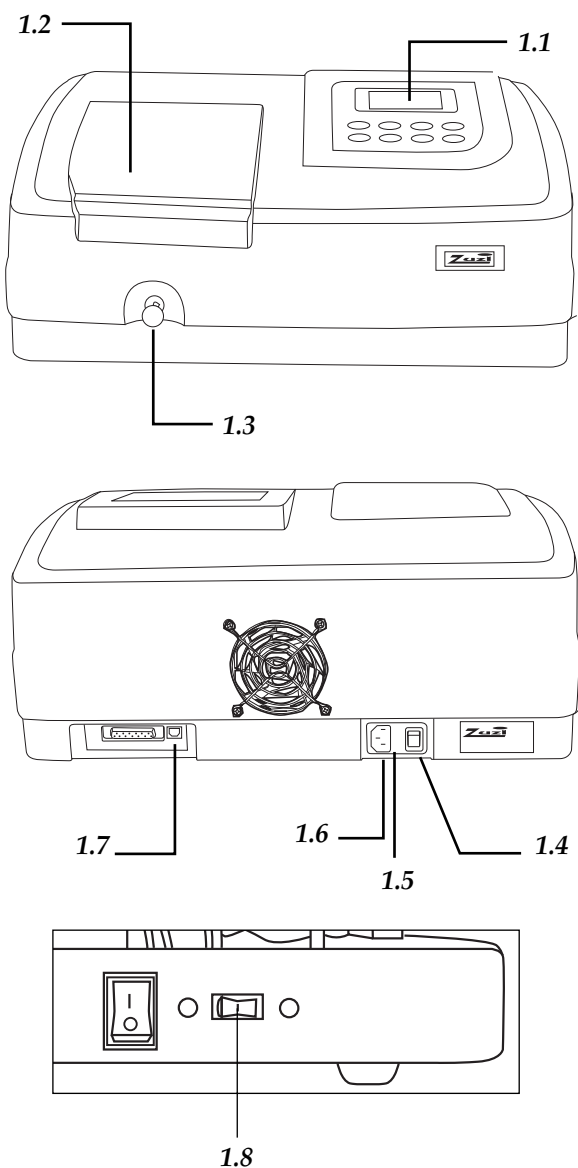
Figure 2

- 2.1 GOTO $\lambda$ : selection of wavelength
- 2.2 ZERO: blank
- 2.3 SET: selection of the different functions
- 2.4 PRINT: Printing of results
- 2.5 (-) left (F1): confirmation of the left command on display
- 2.6 (-) right (F2): confirmation of the right command on display
- 2.7 ▼: parameter decrease
- 2.8 ▲: parameter increase

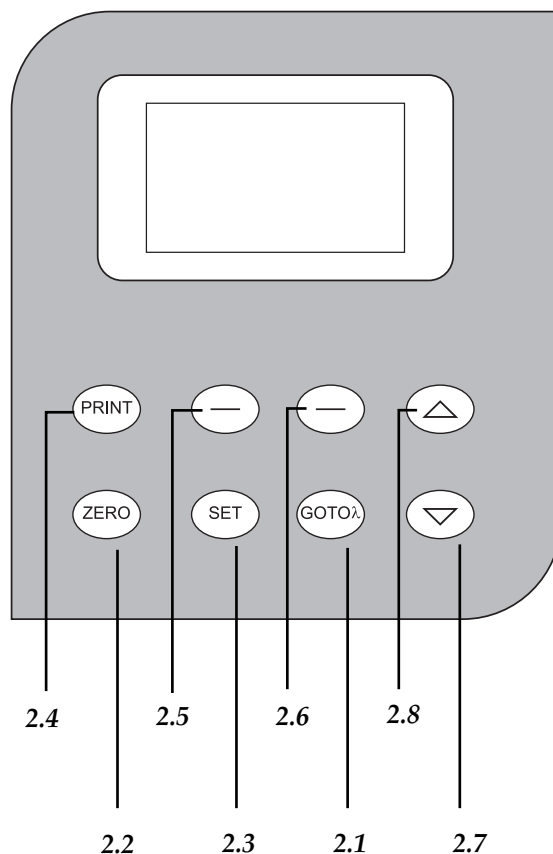### 3. TECHNICAL SPECIFICATIONS

Code	54211020	54211050
Wavelength range	325-1000 nm	200-1000 nm
Spectral bandwidth	4 nm	
Optical system	Simple beam, grating 1200 lines/mm	
Wavelength accuracy	± 1nm	
Wavelength repeatability	0.5 nm	
Wavelength resolution	± 1 nm	
Photometrical range	-0.097/2.5 A, 0/125% T	
Photometrical accuracy	± 0.5 %T	
Photometrical repeatability	± 0.3 %T	
Stray light	0.3% T	
Stability	± 0.002 A/h at 500 nm	
Detector	Silicon photodiode	
Sample compartment	4 cells of 10 mm	
Lamps	Tungsten	Tungsten and deuterium
Output	USB port, parallel port (printer)	
Power	AC 220V/50Hz or AC110V/50HZ	
Dimensions	480x360x160 mm	

### 4. INSTALLTION / SETTING UP

Unwrap the spectrophotometer, take off the involving plastic and take off the protection in which it comes fitted. Take off all the protective items and, without connecting the spectrophotometer to the net, make sure that it does not present any damage because of the shipment. In case the spectrophotometer presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- 4 glass cells
- 2 quartz cells (only model Zuzi 4211/50)
- Software MWave Basic (Compatible with MWave Professional software (Ref. 94251000) not included)
- USB wire
- Schuko wire
- User's manual
- Warranty

**We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.**

#### Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

**PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT.**

Please put the spectrophotometer on top of a horizontal, plane and stable table making a free space of at least 30 cm per side. Do not put the spectrophotometer near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid inflammable or toxic substances in the working area.

The spectrophotometer is supplied with a Schuko standard wire. Please insert the wire that feeds the AC electric current in the base of current 220 V, 50 Hz ± 10% provided with earth wire and to the other end to the spectrophotometer connector.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz  $\pm$  10%.

Setting up

- ◆ Press On/Off (1.4).
- ◆ The equipment initializes and makes an autotest. This process can spend a few minutes; during this time *Zuzi self test...* will be displayed. After this time, *Warm up 20 minutes any key to skip* will be displayed; wait for 20 minutes before operating with the equipment.


*NOTE: Do not open the sample compartment during autotest.*

**BASIC FUNCTIONS**

1. Wavelength adjustment (Fig. 3)

- ◆ Press GOTOλ (2.1) to enter the wavelength adjustment display. Select the desired value of wavelength by pressing ▼ ▲ (2.7) (2.8).
- ◆ Select OK with F1 (2.5) to save the wavelength value.
- ◆ The equipment automatically returns to the main menu.
- ◆ If there is no need of changing the wavelength value, select Return with F2 (2.6) to go to the main menu.


Figure 3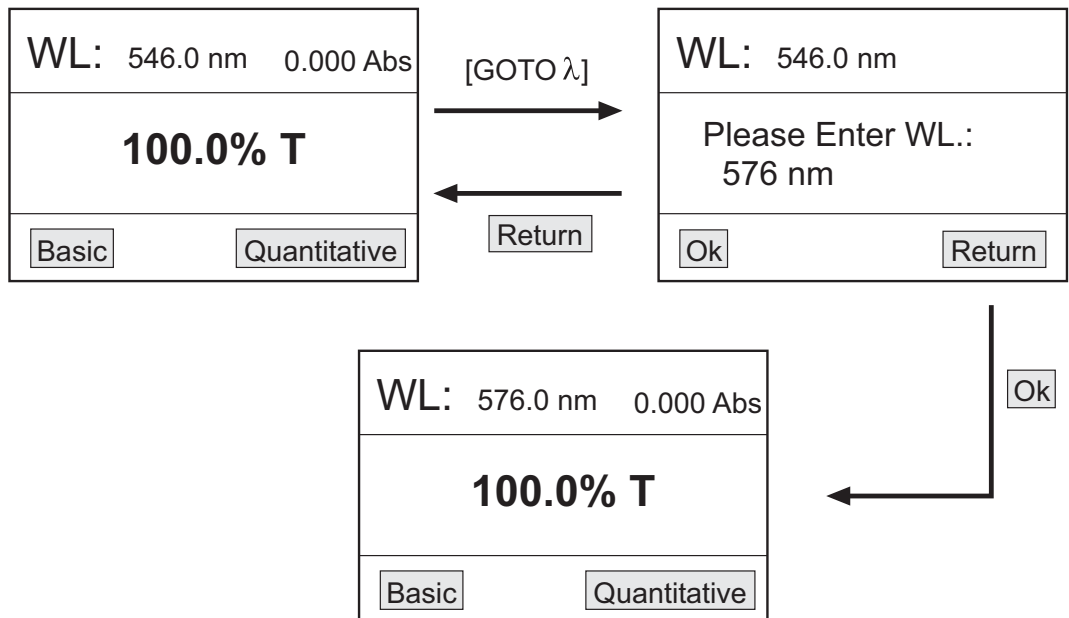

2. Blank (Fig. 4)

- ◆ Put the cell with the reference sample into the cell holder compartment.
- ◆ Press ZERO (2.2); the equipment will make the blank automatically.


Figure 4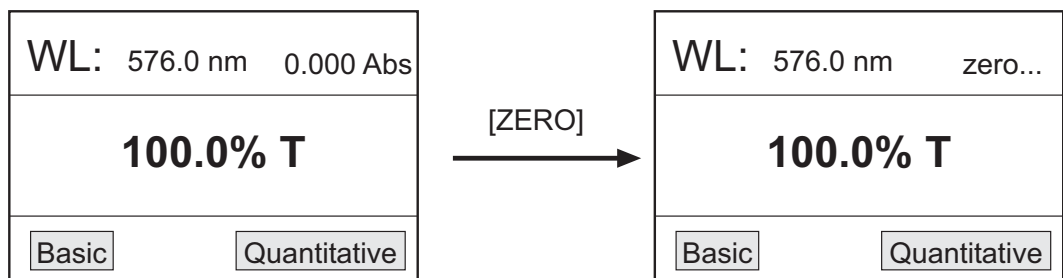

### 3. Utilities (Fig. 5)

- ◆ Press SET (2.3) to enter the function selection display.
- ◆ Use ▼ ▲ (2.7) (2.8) to move along the different functions and select OK with F1 (2.5) to enter the selected function.
- ◆ If no function needs to be selected, select Cancel with F2 (2.6) to return to the main menu.

Figure 5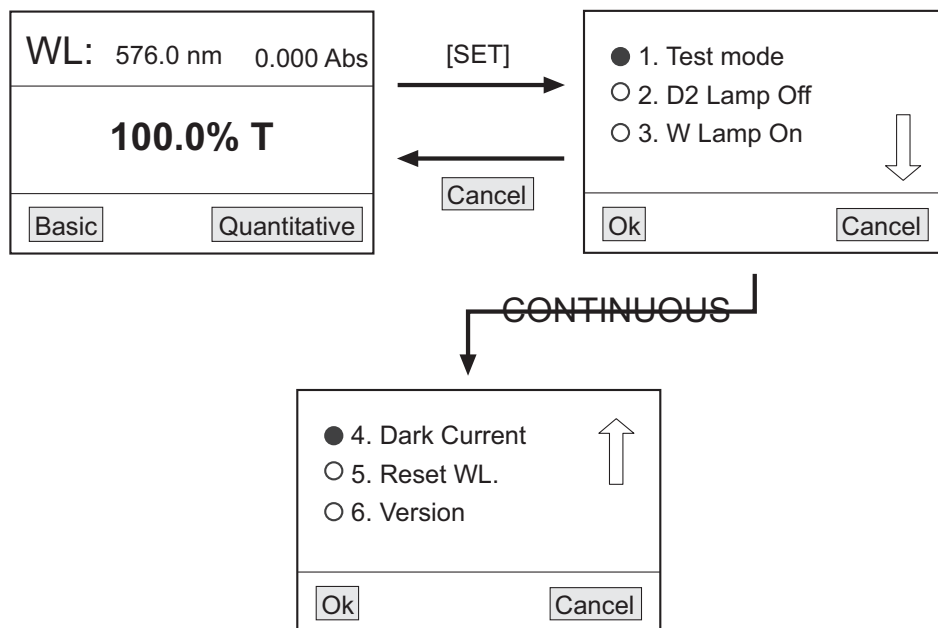


#### 3.1 Test Mode (Fig. 6)

To select the measuring mode, absorbance or transmittance, proceed as follows:

- ◆ At the function selection display, move the cursor with ▼ ▲ (2.7) (2.8) up to the Test Mode function.
- ◆ Confirm OK with F1 (2.5) to enter the test mode selection display.
- ◆ Select the desired measuring mode with ▼ ▲ (2.7) (2.8) : Abs (absorbance), T% (transmittance) or Energy. Confirm OK with F1 (2.5).
- ◆ The symbol ✓ will be displayed beside the selected measuring mode.
- ◆ Confirm Cancel with F2 (2.6) to return to the function selection display.

Figure 6

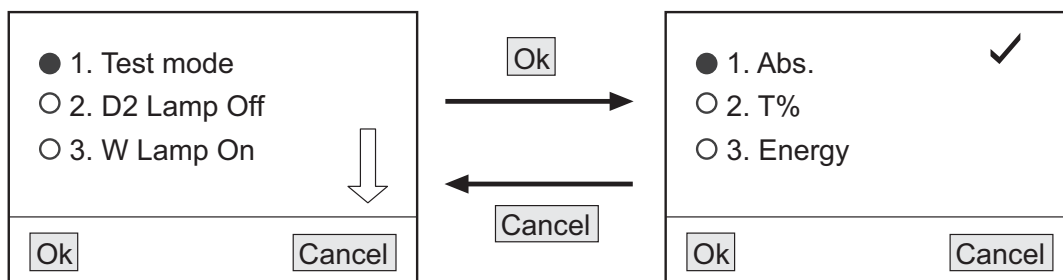#### 3.2 Deuterium lamp on/off (D2 Lamp On/Off) (Fig. 7). Only for model 4211/50


**Note:** For model 4211/50, Zuzi recommends to have the deuterium lamp switched off when working out of the ultraviolet range.

- ◆ At the function selection display, move the cursor with ▼ ▲ (2.7) (2.8) up to the D2 Lamp function.
- ◆ Confirm OK with F1 (2.5) to enter the deuterium lamp state selection display.
- ◆ With ▼ ▲ (2.7) (2.8) select On or Off and confirm OK with F1 (2.5). Beside the selected state the symbol ✓ will be displayed.
- ◆ Confirm Cancel with F2 (2.6) to return to the function selection display.

**Note:** When the deuterium lamp is switched on, it will take a few minutes to warm up.

Figure 7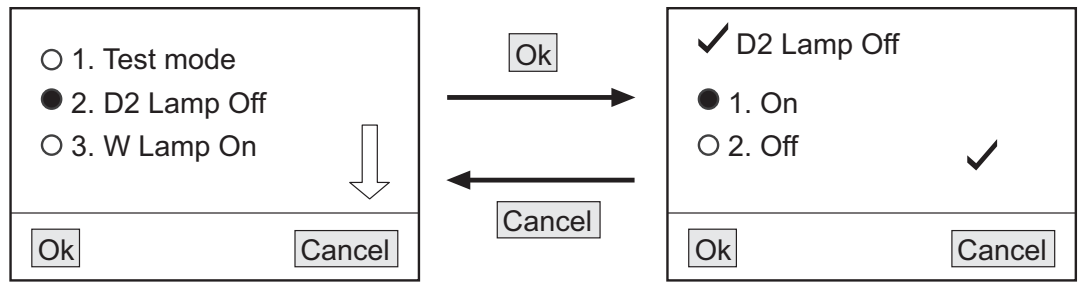


3.3 Tungsten lamp on/off (W Lamp On/Off) (Fig. 8)

Note: For 4211/50, Zuzi recommends to have the tungsten lamp switched off when working out of the visible range.

- ◆ At the function selection display, move the cursor with ▼ ▲ (2.7) (2.8) up to the W Lamp function.
- ◆ Confirm OK with F1 (2.5) to enter the tungsten lamp state selection display.
- ◆ With ▼ ▲ (2.7) (2.8) select On or Off and confirm OK with F1 (2.5). Beside the selected state the symbol ✓ will be displayed.
- ◆ Confirm Cancel with F2 (2.6) to return to the function selection display.

Figure 8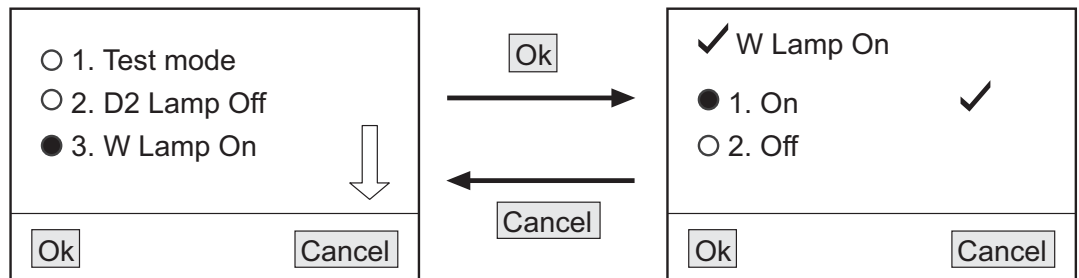


3.4 Dark Current (0%T) (Fig. 9)

Zuzi recommends to carry out this function whenever changing the environmental conditions (changes in temperature or equipment location, peaks in electrical current, etc.). For this purpose:

- ◆ At the function selection display, move the cursor with ▼ ▲ (2.7) (2.8) up to the Dark current function.
- ◆ Confirm OK with F1 (2.5); Get dark current will be displayed.
- ◆ After a few seconds, the equipment will automatically return to the function selection display and will be ready to operate.


Figure 9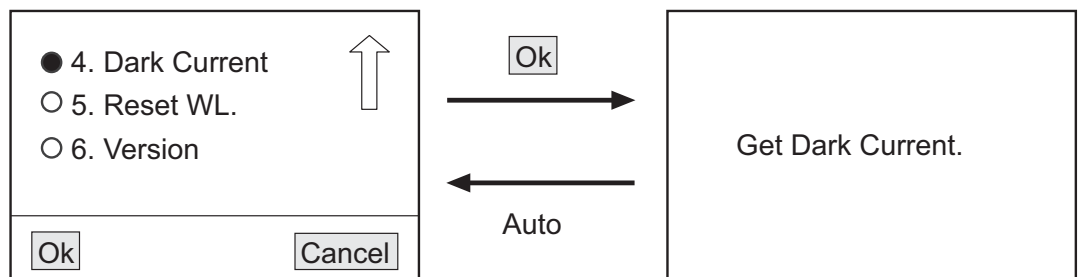

3.5 Wavelength calibration (Reset WL) (Fig. 10)

- ◆ At the function selection display, move the cursor with ▼ ▲ (2.7) (2.8) up to the Reset WL function.
- ◆ Confirm OK with F1 (2.5); Calibrating WL... will be displayed.
- ◆ After a few seconds, the equipment will automatically return to the function selection display and will be ready to operate.

Figure 10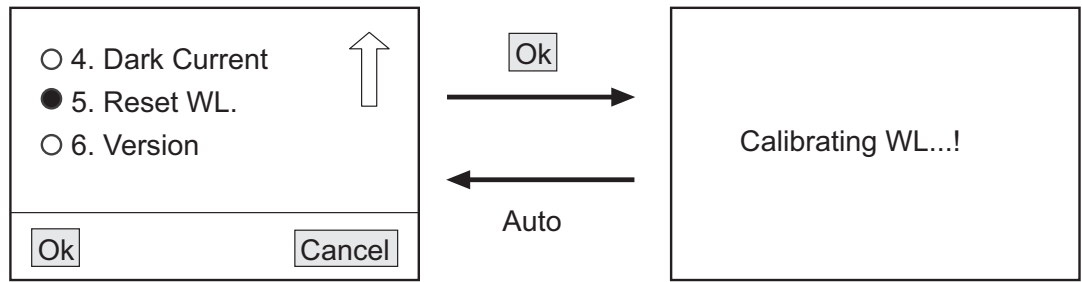

**4. Sample analysis**


Before proceeding to analysis clean the cell with distilled water. Wipe the cell with a cellulose paper to remove drops of water and fingerprints.

**4.1 Basic mode**

To determine the absorbance or transmittance of different samples at a certain wavelength:

- ◆ Put the cell with the blank reference into the cell holder and place it into the light path.
- ◆ Select Basic with F1 (2.5). The equipment will make the blank automatically and then will enter the measuring mode display (Fig. 11).

Figure 11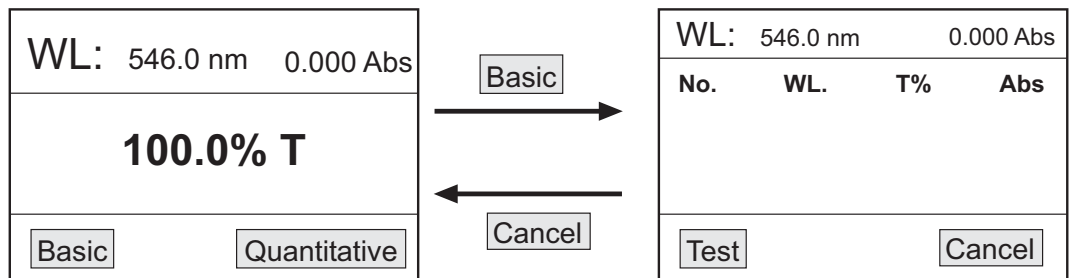


- ◆ On the upper part, the selected wavelength (WL) and 100% T or 0.0000 Abs depending on the working measuring mode will be displayed. Immediately below a chart headline will appear:

No.	WL	T%	Abs
Number assigned in order to each result	Wavelength	Transmittance value	Absorbance value

The results of the successive analysis will be shown on this chart.

- ◆ Put the cell with solution to be analyzed into the light path. Select Test with F1 (2.5); the result of the measurement will be displayed

Figure 12

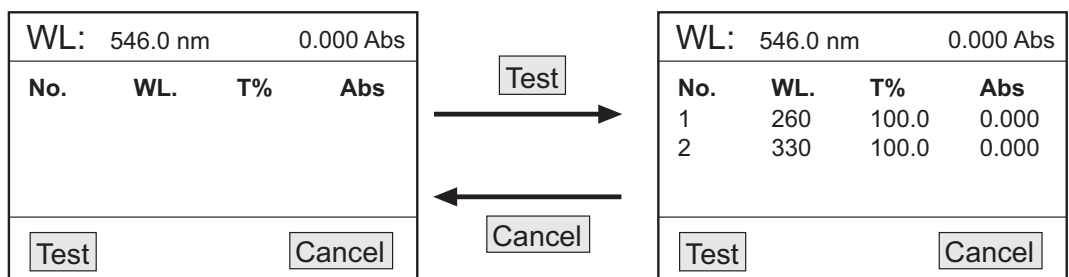- ◆ Repeat this process with the different samples. The spectrophotometer can save up to 200 data.
- ◆ To return to the main menu select Cancel with F2 (2.6).

*Note: At the basic mode, the wavelength can be adjusted and the blank can be made as many times as needed by pressing GOTO λ (2.1) and ZERO (2.2) respectively.*

#### 4.2 Quantitative mode

To determine the concentration of different samples from a curve equation in which the absorbance and concentration are related:

$$C=K \cdot A+B$$

*C: concentration*


*A: absorbance*

*K and B: curve coefficients*

◆ Select Quantitative with F2 (2.6) to enter the quantitative mode main menu; the following options will be displayed (Fig. 13):

1. Create Curve
2. Load Curve: to load a curve previously created and saved.
3. Remove Curve: to delete a function previously saved.

Figure 13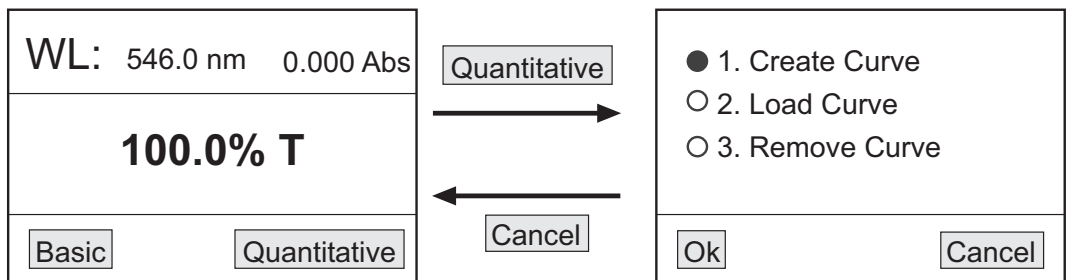


#### Create a curve

◆ Select the option Create curve with ▼ ▲ (2.7) (2.8) and then confirm OK with F1 (2.5). The two methods that can be followed will be displayed (Fig. 14):

- Coefficient: this method is used when the equation of the curve is known. First of all, the equation coefficients must be introduced and then, proceed to the analysis of the samples of unknown concentration.
- Standard curve: this method is used when the equation of the curve is unknown. First of all, the curve must be created from standard samples of known concentration and then, proceed to the analysis of the samples of unknown concentration.


Figure 14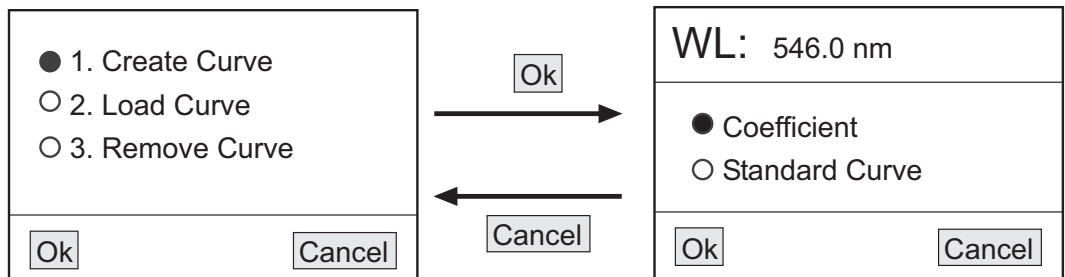

#### 1. Coefficient method


◆ Select the option Coefficient with ▼ ▲ (2.7) (2.8) and then confirm OK with F1 (2.5) (Fig. 15).

Figure 15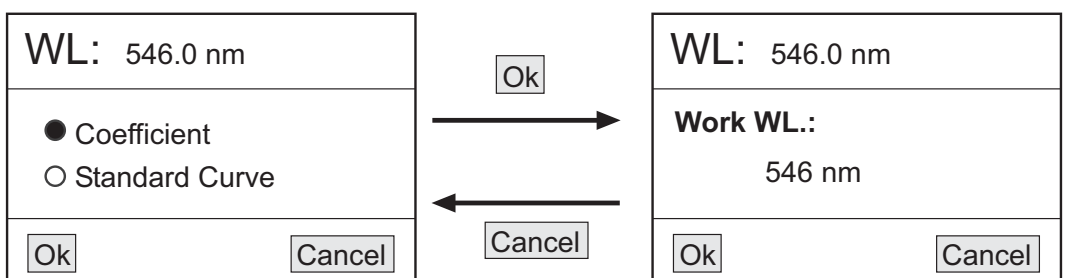


◆ The wavelength selected value will be displayed. To modify this value press ▼ ▲ (2.7) (2.8) and then confirm OK with F1 (2.5) (Fig. 16).

Figure 16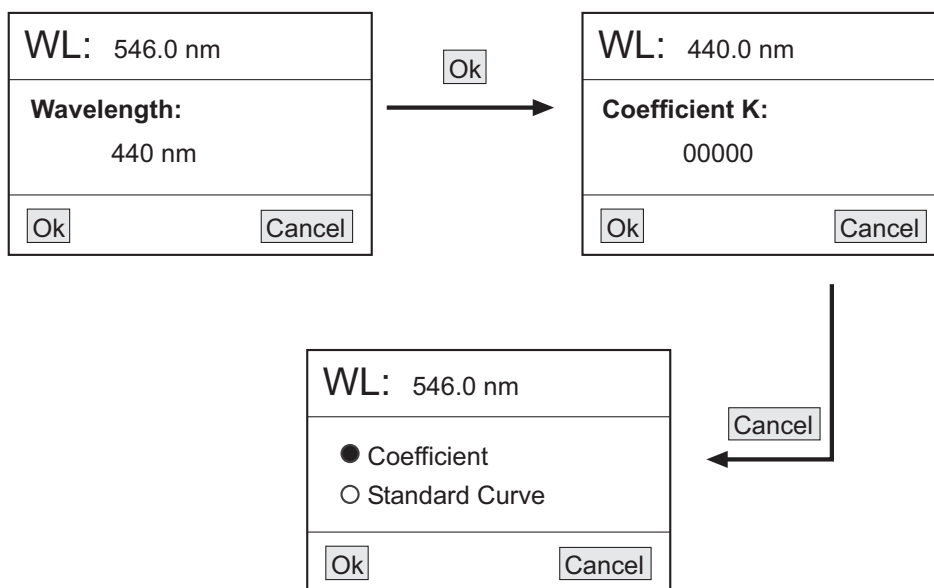

◆ Enter the value of the K coefficient on the new display. Adjust the value digit by digit with ▼ ▲(2.7) (2.8); to go from one digit to the next one select OK with F1 (2.5). When the value of the last digit is adjusted select OK with F1 (2.5) (Fig. 17).

Figure 17

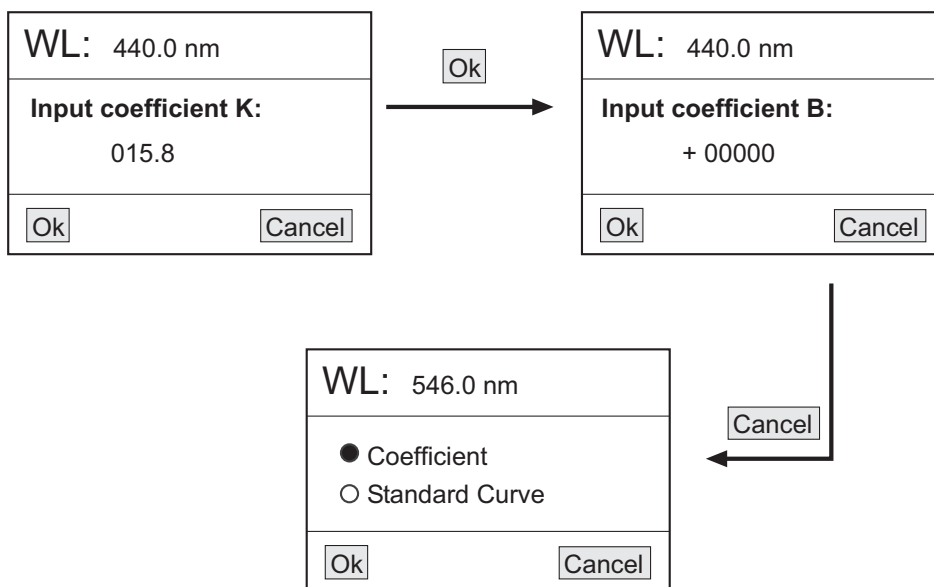◆ The equipment will enter the B coefficient value adjustment display. Adjust the value digit by digit with ▼ ▲(2.7) (2.8); to go from one digit to the next one select OK with F1 (2.5). When the value of the last digit is adjusted select OK with F1 (2.5) (Fig. 18).

◆ The new display will show the curve and its equation.  
 ◆ To determine the concentration of the different samples to be analyzed, put each sample into the light path and select Test with F1 (2.5). The results will be shown in order in a chart (Fig. 19).

2. Standard curve method

◆ Blank  
 - Select the option Standard curve with ▼ ▲(2.7) (2.8) and then confirm OK with F1 (2.5). Please Insert Blank will be displayed (Fig. 20).

Figure 18

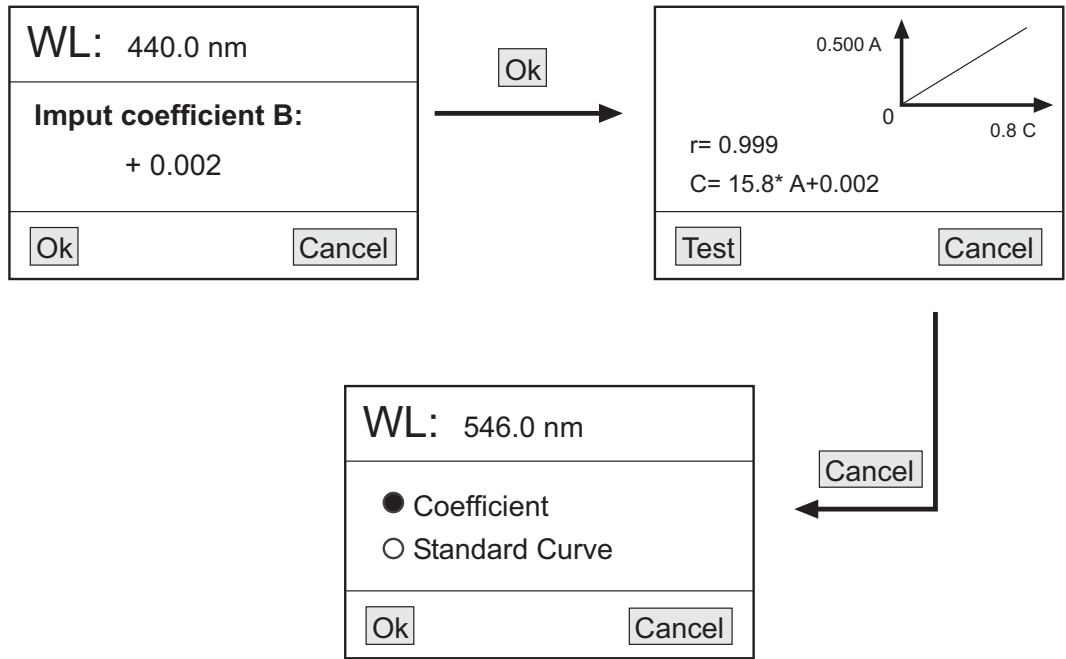
Figure 19


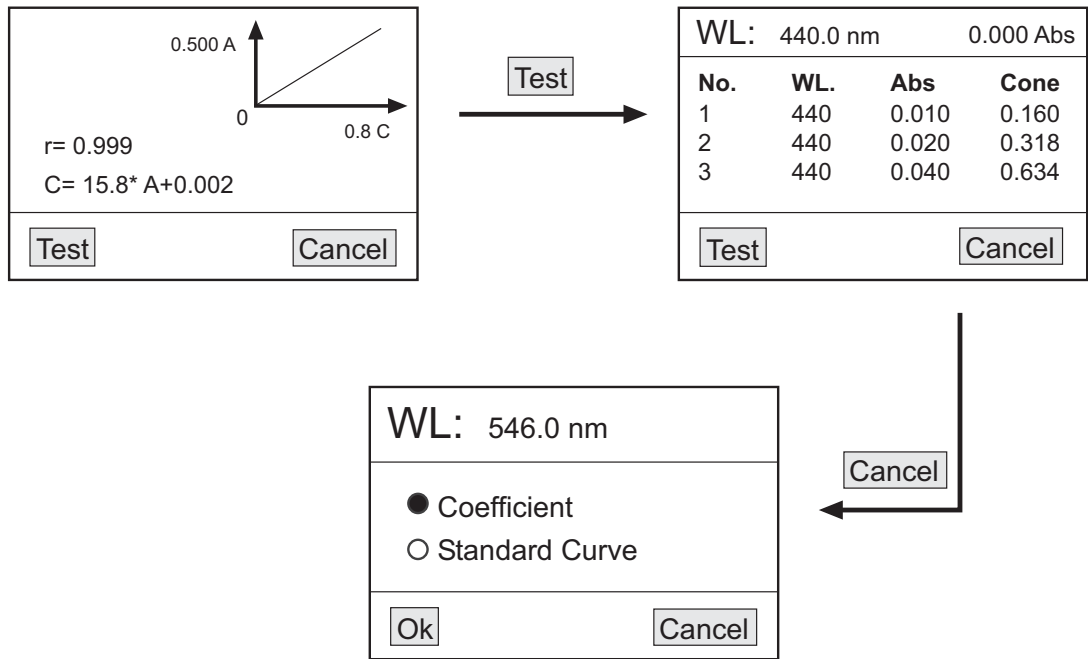
Figure 20

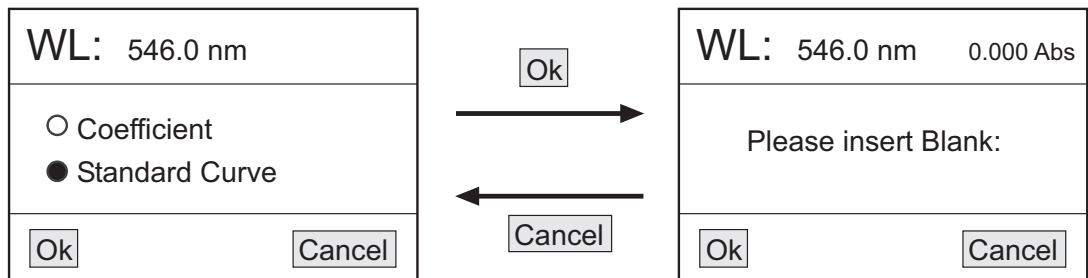- Put the cell with the blank sample into the cell holder and place it into the light path.
- Press GOTO  $\lambda$  (2.1) and adjust the working wavelength value.
- Select OK with F1 (2.5) to make the blank.

- ◆ Input the number of samples
  - After blanking, the system will ask for the number of standard samples that are going to be used to create the curve.

**NOTE: The maximum number of standard samples is 9.**

- Input the number of samples with  $\blacktriangledown \blacktriangle$  (2.7) (2.8) and then select OK with F1 (2.5) (Fig. 21).

Figure 21

WL: 546.0 nm	0.000 Abs
<b>Number: 3</b>	
Ok	Cancel

- ◆ Input the concentration of the standard samples
  - Once selected the number of Standard samples, their concentration must also be introduced.

- Put the first sample into the cell holder and place it into the light path.

- Input the concentration value of the sample. Adjust the value digit by digit with  $\blacktriangledown \blacktriangle$  (2.7) (2.8); to go from one digit to the next one select OK with F1 (2.5). When the value of the last digit is adjusted select OK with F1 (2.5)

- Proceed in the same way for the rest of the standard samples.

- ◆ Curve visualization

- Once the concentration of all the standard samples has been input, the curve and its equation will be displayed (Fig. 23).

**Note: The curve will be automatically saved in the equipment memory. The equipment can save up to 200 different curves.**

- ◆ Measurement of the concentration of the samples to be analyzed

- Put each sample into the light path and select Test with F1 (2.5). The results will be shown in order in a chart (Fig. 24).

Load a curve

- ◆ Select the option Load curve with  $\blacktriangledown \blacktriangle$  (2.7) (2.8) and then press OK with F1 (2.5). All the saved curves will be displayed (Fig. 25).

**Note: The equipment can save up to 200 curves in its memory.**

- ◆ Select the desired curve with  $\blacktriangledown \blacktriangle$  (2.7) (2.8) and then confirm OK with F1 (2.5). The graph and curve data will be displayed (Fig. 26).

Figure 23

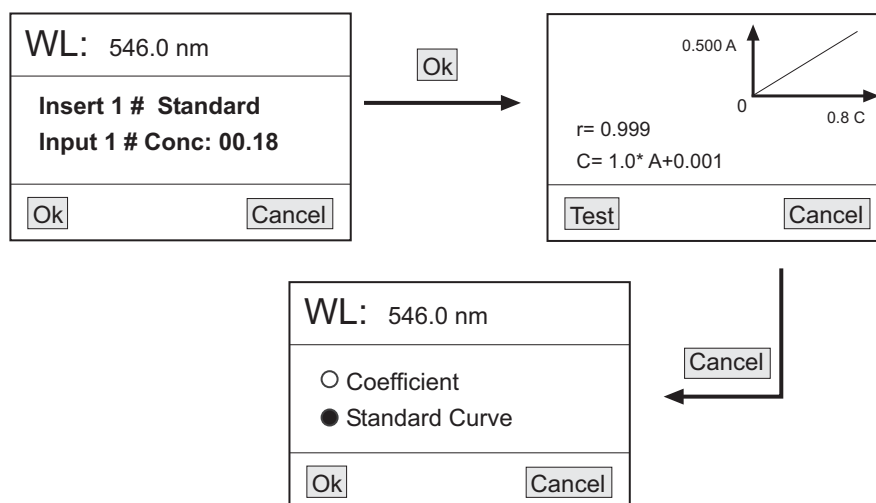
Figure 24

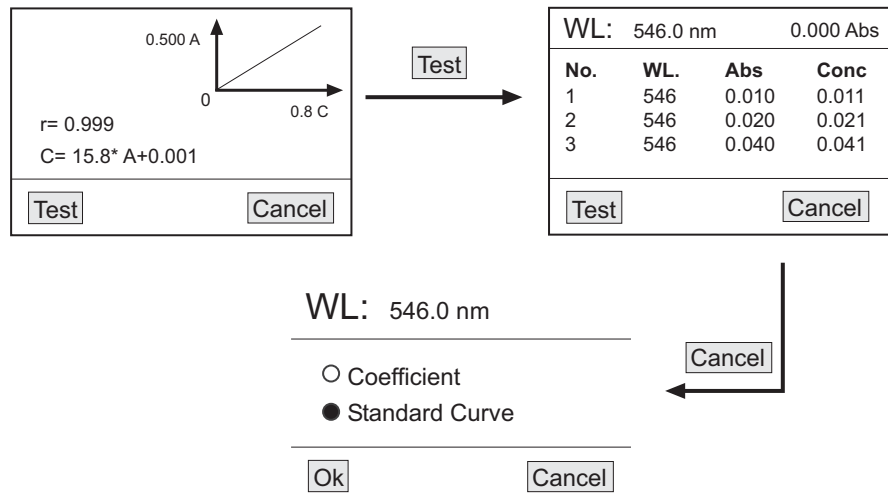
Figure 25


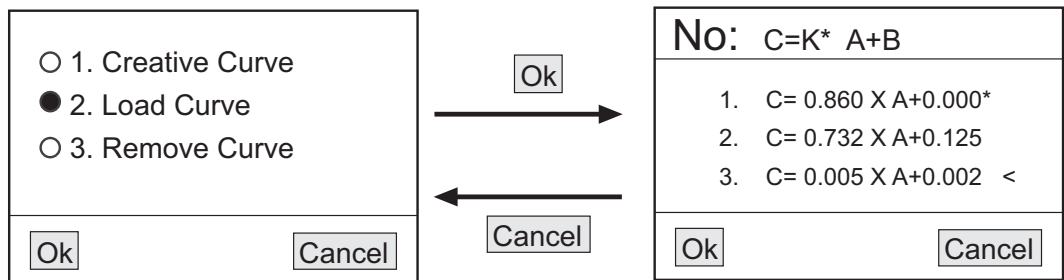
Figure 26

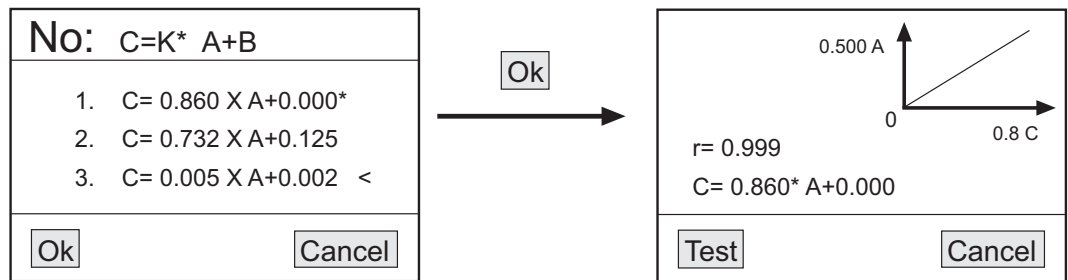◆ To make new measurements using that curve, put the sample into the light path and select Test with F1 (2.5). The results will be shown in order in a chart.

Delete a curve


◆ Select the option Remove curve with ▼ ▲ (2.7) (2.8) and then confirm OK with F1 (2.5). All the saved curves will be displayed.

◆ Select with ▼ ▲ (2.7) (2.8) the curve to be deleted and confirm OK with F1 (2.5). The curve will be automatically deleted from the screen and from the memory of the spectrophotometer.

4.3 Deleting and printing results

◆ At the result chart of any measuring mode, press PRINT (2.4) and the following options will be displayed (Fig. 27).

Figure 27

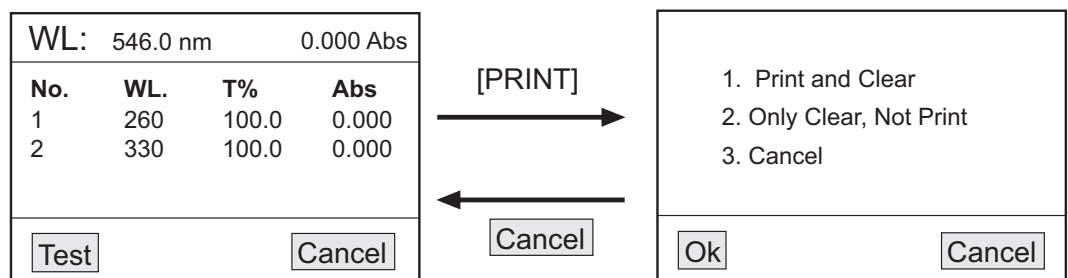1. Print and clear: for printing and then, deleting data
2. Only clear, not print: for deleting data but not printing
3. Cancel

Select the desired option with ▼ ▲ (2.7) (2.8) and then confirm OK with F1 (2.5).

If you are not using the spectrophotometer for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

### Security

- ◆ The spectrophotometer must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ You should put the spectrophotometer on top of a horizontal, plane, stable table, having a safety area of at least 30 cm per side.
- ◆ Do not place the spectrophotometer near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- ◆ During operation dangerous materials such as flammable or pathological substances must be out of the safety area.
- ◆ When you are not using the spectrophotometer for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the spectrophotometer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the spectrophotometer (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

## 5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

*Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.*

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The spectrophotometer is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never try to repair the spectrophotometer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the spectrophotometer (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Zuzi Technical Assistance Department.

### Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the spectrophotometer and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

## 6. TROUBLESHOOTING

TROUBLE	CAUSE	SOLUTION
The equipment does not turn on	<ul style="list-style-type: none"> <li>- The wire is not connected</li> <li>- The power source is broken</li> <li>- Voltage error</li> <li>- Fuse blown or defect in an electric component</li> </ul>	<ul style="list-style-type: none"> <li>- Connect the wire</li> <li>- Change the power source</li> <li>- Check the voltage of the electric installation</li> <li>- Contact the Zuzi Technical Assistance Department</li> </ul>
100% T (0.000A) can not be selected	<ul style="list-style-type: none"> <li>- The sample is not well placed in the cell holder</li> <li>- The lamp is worn out or defective</li> <li>- The lamp is turned off</li> <li>- Defect in an electric component</li> </ul>	<ul style="list-style-type: none"> <li>- Put correctly the sample into the cell holder</li> <li>- Change the lamp</li> <li>- Look up the user's manual, lamp management section</li> <li>- Contact the Zuzi Technical Assistance Department</li> </ul>
The equipment strays and there is noise	<ul style="list-style-type: none"> <li>- Not enough period of warm up</li> <li>- Use of glass cell in the UV range</li> <li>- The lamp is worn out or defective</li> <li>- Instable power source</li> <li>- Significant change of temperature</li> <li>- The lamp is not adjusted</li> <li>- Defect or dirtiness in detector or defect in an electric component</li> </ul>	<ul style="list-style-type: none"> <li>- Leave 20 minutes to warm up</li> <li>- Use a quartz cell</li> <li>- Change the lamp</li> <li>- Stabilize the voltage</li> <li>- Adjust correctly the lamp</li> <li>- Contact the Zuzi Technical Assistance Department</li> <li>- Contact the Zuzi Technical Assistance Department</li> </ul>
Incorrect readings	<ul style="list-style-type: none"> <li>- Not enough volume of sample</li> <li>- Equipment out of calibration</li> <li>- Bubbles or particles in sample</li> <li>- Lost of vapours of sample</li> <li>- Wrong blanking</li> </ul>	<ul style="list-style-type: none"> <li>- Fill the cell with sample</li> <li>- Contact the Zuzi Technical Assistance Department</li> <li>- Check sample preparation, analytical procedure</li> <li>- Adapt the sample to the equipment</li> <li>- Use the adequate ventilation</li> <li>- Check the wavelength and the analytical procedure</li> <li>- Make Dark current and calibrate the equipment</li> </ul>

**At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.**

**The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.**

**Please inquire at the community administration for the authorized disposal location.**

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