

INSTRUCTION MANUAL

Digital Stopwatch

Code KKB001



READING FUNCTION

Normal display: Hour, minutes, seconds, and day of week

Press A: For alarm time

Press B: Mode

Press C: For month and day of week


STOP FUNCTION

Press B for stop function

Press C for count & stop

Press A for lap & reset

ALARM & CHIME ON/OFF

Chime ( /.....) on and off.

In normal display hold A and then press B.

Chime signal appears, chime on. Chime signal disappears, chime off.

Alarm () / ) on and off.

In normal display hold A and then press C.

Alarm signal appears, alarm on. Alarm signal disappears, alarm off.

ALARM TIME SETTING

Hold B about 3 seconds or press B repeatedly (day of week signal above Monday flashing).

Press C or A to change hour / minutes function, then advance figure by pressing A or C.

Press B to normal display.

NORMAL TIME SETTING

Hold B about 3 seconds and then press B once, or push B repeatedly (day of week signal above Tuesday flashing).

Select flashing digit to be set by press C or A.

Advance figure by pressing A or C.

Select 12/24 hour cycle option in normal time.

Hold A then press B or when time (hour) setting, 12/24 hour cycle option will appear alternatively on every 24 hour cycle during hour advance.

Month and date change by holding C and then pressing A or cannot be changed.