DECLARACIÓN DE CONFORMIDAD CE
CE DECLARATION OF CONFORMITY
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

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8. RECOMMENDED ADAPTABLE ACCESSORIES

CODE	DESCRIPTION
90404351	Sodium vapour lamp
90404910	Test tube of 100 mm
90404920	Test tube of 200 mm
90404901	Glass joints for test tubes, 15 mm Ø
90404691	Circular rubber washer for test tube
90404692	Test tube stopper
90400001	Quartz piece for polarimeter calibration, value +5°
90400002	Quartz piece for polarimeter calibration, value -5°
90400003	Quartz piece for polarimeter calibration, value +17°
90400004	Quartz piece for polarimeter calibration, value -17°
90400005	Quartz piece for polarimeter calibration, value +35°
90400006	Quartz piece for polarimeter calibration, value -35°

**INSTRUCTIONS ON ENVIRONMENT PROTECTION**

At the end of its life cycle, please, do not dispose of this polarimeter by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Zuzi polarimeter model 430. We highly recommend looking after this equipment according to what is stated in this manual.

Zuzi develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the Zuzi polarimeter model 430, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the polarimeter avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the polarimeter to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the polarimeter.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by email to info@auxilab.com or by telephone: +34 807 117 040 (1.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Zuzi Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

- ◆ Remove the lid with the grilles (5.1) placed at the left side of the polarimeter; this way, the sodium lamp (5.2) can be seen easily.
 - ◆ With the help of a screwdriver loosen the screw (5.3) that holds the base of the lamp.
 - ◆ Holding it by the base and with the help of both hands, take the lamp out.
 - ◆ Put the new lamp and tighten the holding screw again.
- Replacement of the fuse
- ◆ Fuses (1.6) are located at the back side of the polarimeter.
 - ◆ With the help of a screw driver take the fuse compartment out and replace the fuse.

Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the polarimeter and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

7. TROUBLESHOOTING

TROUBLE	CAUSE	SOLUTION
Problem with light source	Light source is dirty Light source is damaged	Open the lid and clean the lamp Replace the lamp
After pressing the on/off switch, the equipment does not turn on	Energy does not reach the equipment The fuse is blown	Check the cable and the socket Check the 3 A fuse and replace if necessary
The DC switch is on but the sodium lamp is off	The lamp is not warmed The lamp is damaged or run down	Let the lamp warm for 5 min more Change the lamp
Instable readings	The lamp is not working on direct current (DC) The sodium vapour of the lamp is in bad conditions	Put the AC/DC switch in DC position Change the lamp
Too high noise	Mechanical friction	Gears are not well lubricated

- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the polarimeter (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

6. MAINTENANCE AND CLEANING


To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The polarimeter is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never try to repair the polarimeter by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the polarimeter (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Zuzi Technical Assistance Department.

IMPORTANT: before changing the lamp or the fuse, make sure that the polarimeter is disconnected from the net. Check you are using the proper lamps, as other types can provoke malfunction.

Replacement of the lamp (Figur 5)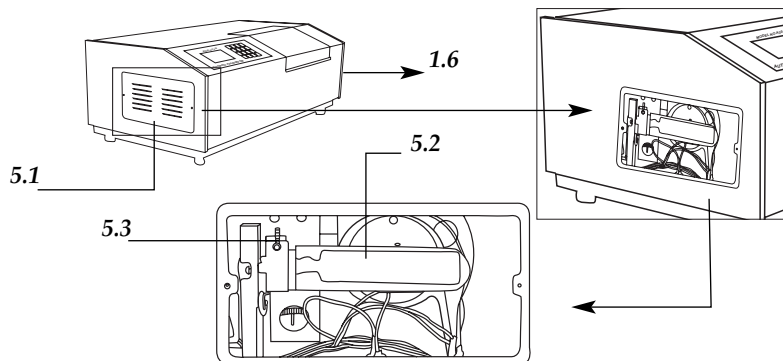

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

Ingles	14-24
Castellano	3 -13
Francés	25-35

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT	15
2. DESCRIPTION	16
3. TECHNICAL SPECIFICATIONS	17
4. THEORETICAL FOUNDATIONS	17
5. INSTALLATION / SETTING UP	19
6. MAINTENANCE AND CLEANING	22
7. TROUBLESHOOTING	23
8. RECOMMENDED ADAPTABLE ACCESSORIES	24
ANNEX I: CE CERTIFICATE	

1. USES OF THE INSTRUMENT


The Zuzi polarimeter model 430 is a complete equipment that allows determination of the optical rotation, specific rotation, concentration and international sugar degrees of optically active substances. Thus, it is very useful in several fields of industrial and quality control processes as well as in research. Among the main application of the polarimeter there are:

- Research: isolation of crystallized substances, evaluation and characterization of optically active compounds, kinetics and changes in concentration.
- Pharmaceutical industry: determination of the optical rotation of medicines (antibiotics, codeine, analgesics) and drugs.
- Food industry: determination of the sugar and condiment content.
- Chemical industry: study of biopolymers, synthetic and natural polymers, etc.


2. DESCRIPTION

- 1.1 Control panel.
- 1.2 Sample compartment.
- 1.3 Power on/off.
- 1.4 AC/DC switch.
- 1.5 Socket.
- 1.6 Fuses.
- 1.7 RS232 output.
- 1.8 Grille.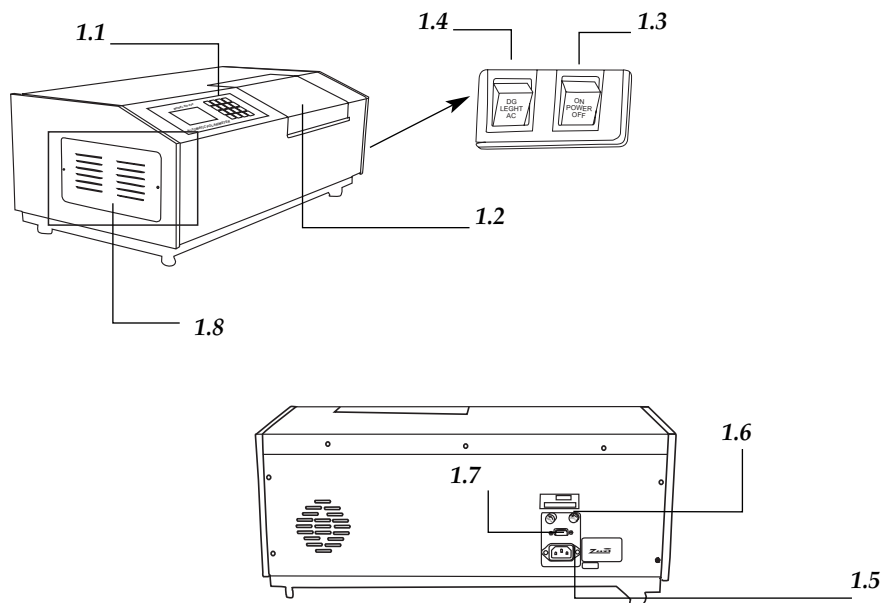

- 2.1 Display.
- 2.2 Zero.
- 2.3 →
- 2.4 Measure.
- 2.5 ↵ Enter for parameter confirmation.
- 2.6 [0]-[9]: numeric keypad.
- 2.7 Dot.
- 2.8 Reset: repetition of the measurement.

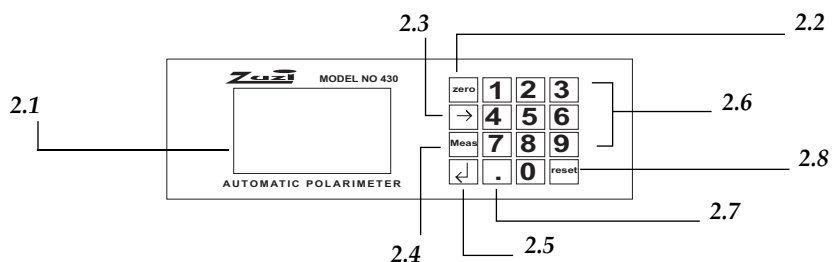- ◆ By using the numeric keypad (2.6) select the desired measurement mode and then:
 - If there is no need to change any parameter press Measure (2.4) and then ↵ (2.5); the equipment will automatically enter the measurement mode.
 - If it is necessary to change any parameter press ↵ (2.5) to go from one parameter to the next; adjust the parameters by using the numeric keypad (2.6). Once the last parameter has been adjusted press ↵ (2.5) and the equipment will enter the measurement mode.
- ◆ “0.000” will be displayed.
- ◆ Put the tube with distilled water or the blank solution into the sample compartment (1.2), let the displayed reading stabilize and press Zero (2.2).
- ◆ Remove the blank solution or distilled water from the sample compartment and put the sample to be measured; the equipment will automatically measure and display the result.

Note: the equipment will perform the number of measurements (n) previously selected and will finally display the average value of all measurements.

Note: if it is necessary to repeat the measurement, press Reset (2.8).

- ◆ Before measuring a new sample, it is necessary to make the zero. For this:
 - Remove the sample with the already analyzed sample from the sample compartment.
 - Put the blank sample or distilled water, let the displayed reading stabilize and press Zero (2.2).
 - Put the new sample and the equipment will automatically measure and display the result.
- ◆ To change the measurement mode press ↵ (2.5) and proceed as previously described.

Note: before turning the equipment off, remove the last analyzed sample from the sample compartment, let the displayed reading stabilize and press Zero (2.2); then turn the equipment off.

Security

- ◆ The polarimeter must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the polarimeter on top of a horizontal, plane, stable table, having a safety area of at least 30 cm per side.
- ◆ Do not place the polarimeter near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- ◆ During operation dangerous materials such as flammable or pathological substances must be out of the working area.
- ◆ When you are not using the polarimeter for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of the lamp or a fuse).
- ◆ Never try to repair the polarimeter by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz \pm 10%.

Operation mode

Turning the equipment on

- ◆ Make sure that the AC/DC switch (1.4) is in AC position and press the power on/off (1.3); check by looking through the grille (1.8) that the sodium lamp is turned on.
- ◆ Let the lamp warm up for 5 min. Important: the AC/DC switch (1.4) must be in AC position to let the vapour of the lamp warm.
- ◆ After pre-warming time, put the AC/DC switch (1.4) in DC position to start measuring samples. **Important:** if the switch is in AC position during measurement, the shown reading will fluctuate notably.

Note: if when putting the switch in DC position the lamp lights off, press several times the AC/DC switch (1.4) until the lamp lights on again.

Measurements

The equipment presents different measuring modes in each one of which it is necessary to establish some parameters. The following chart shows the different measuring modes and parameters required in each one of them:

Reading		Parameters required
Mode 1	Rotation angle (α)	n
Mode 2	Specific rotation ($[\alpha]$)	L, C, n
Mode 3	Concentration (C)	L, α , n
Mode 4	Sugar degrees ($^{\circ}$ Z)	n

Being:

n: number of measurements to perform up to a maximum of 6.

L: sample length (dm.)

C: concentration of the optically active substance (g/100 mL).

Note: when the number of measurements (n) is higher than 1, the reading shown on display will be the average of all the measurements performed and will be displayed as α_{AV} , $[\alpha]_{AV}$, C_{AV} or Z_{AV} depending on the selected measuring mode.

To perform measurements proceed as follows:


- ◆ Press \downarrow (2.5); at the upper side the measurement mode and the required parameter will be displayed.

3. TECHNICAL SPECIFICATIONS

Code	50410030
Model	430
Range	45 $^{\circ}$; \pm 130 $^{\circ}$ Z \pm
Min. displayed value	0.001 $^{\circ}$
Accuracy	\pm (0.001 $^{\circ}$ + measured value \times 0.05%); \pm (0.003 $^{\circ}$ Z + measured value \times 0.05%)
Reproducibility	\leq 0.003 $^{\circ}$; \leq 0.003 $^{\circ}$ Z
Min. Sample transmittance	10%
Monochromatic light source	Sodium lamp and filter (λ : 589.44 nm)
Tubes	100 mm and 200 mm
Output	RS232
Power	220 V, 50 Hz \pm 10%
Dimensions	60x32x20 cm
Weight	28 Kg

4. THEORETICAL FOUNDATIONS

Figure 3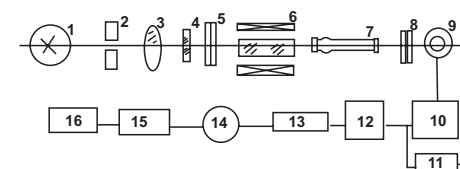

3.1 Light source.

3.2 Aperture.

3.3 Lens.

3.4 Coloured filter.

3.5 Polarizer A.

3.6 Magnetic rotation coil.

3.7 Sample.

3.8 Polarizer B.

3.9 Photo-multiplier.

3.10 Pre-amplifier.

3.11 Automatic high voltage.

3.12 Frequency amplifier.

3.13 Amplifier power.

3.14 Servo motor.

3.15 Worm-gear.

3.16 Counter

Zuzi polarimeter 430 is based on the optical null principle to obtain accurate readings.

For this purpose, it uses a 20 W sodium vapour lamp as the light source. Firstly, the aperture (3.2) and the lens (3.3) form a parallel light beam, that, after passing through the polarizer A (3.4) becomes a polarized light beam whose vibrating plane is $00'$ (Figure 4a). When the polarized light beam passes through the Faraday coil (3.5), its vibrating plane will swing in a $\pm\beta$ angle at a frequency of 50 Hz. After passing through the polarizer B (3.8) the light beam is projected into the photo-multiplier (3.9) so as an alternating current electric signal is generated.

According to the optical null principle, the equipment takes as the optical null position ($\alpha = 0^\circ$) the point in which the axes of the polarizers A and B are orthogonally arranged ($00' \perp PP'$) and therefore the light beam is completely blocked and does not reach the photo-multiplier.

When an optically active sample is placed, the vibrating plane will rotate into an α angle (Figure 4b) (optical rotation of the sample) at a certain frequency F that finally will become an electric signal of the same frequency F. This signal makes the servo motor spin (3.14), which at the same time makes the polarizer A turn in an α angle until the optical null position is restored and the frequency F signal disappears (Figure 4c).

The α angle that the polarizer A has rotated will be displayed as the reading of the optical rotation of the analyzed sample.

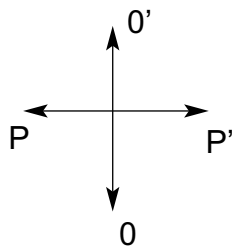
Figure 4a

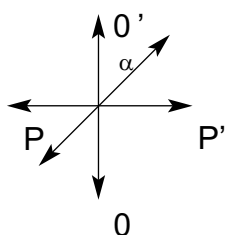
Figure 4b

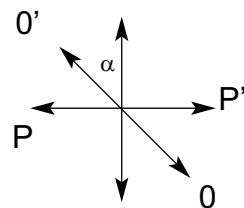
Figure 4c

NOTE: $00'$ is the polarization axe of polarizer A and PP' is the polarization axe of polarizer B

Figure 4a: The polarized light generated by the polarizer A vibrates in the plane $00'$.

Figure 4b: When light passes through an optically active sample its vibrating plane swings in a α angle.

Figure 4c: The polarizer A rotates in a α angle to compensate the optical rotation of the sample and to restore the optical null position; the α value will be displayed with the correct sign.

5. INSTALLATION / SETTING UP

Preliminary inspection

◆ Unwrap the polarimeter, take off the involving plastic and take off the protection in which it comes fitted. Take off all the protective items and, without connecting the polarimeter to the net, make sure that it does not present any damage because of the shipment. In case the polarimeter presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

◆ Please check that all the accessories are enclosed with the equipment:

- 100 mm-long tube and 200 mm-long tube
- 8 Rubber washers
- 8 Glass joints
- Fuses
- Screw driver
- Schuko wire
- User's manual
- Warranty certificate

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT.

◆ Please put the polarimeter on top of a horizontal, plane and stable table making a free space of at least 30 cm per side.

◆ Do not put the polarimeter near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations in the working area.

◆ Avoid the presence of inflammable or toxic substances in the working area.

◆ The polarimeter is supplied with a Schuko standard wire. Please insert the wire that feeds the AC electric current in the base of current 220 V, 50 Hz \pm 10% provided with earth wire and to the other end to the polarimeter connector (1.5).