INDEX OF LANGUAGE-ÍNDICE DE IDIOMAS

ESPAÑOL.....	2
ENGLISH.....	5

CALORÍMETRO DE JOULE

JOULE'S HEATING UNIT

REF. QLJ003

CALORÍMETRO DE JOULE

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Theory

When electric current is made to flow through a conductor, rise in temperature of conductor can be observed. This is due to the heat generation because of resistance offered by the conductor to the flow of electric current and is analogous to the heat produced on account of motion of a body against friction.

The laws governing production of heat on passage of electric current (i.e., heating effect of electric current) were given by Joule. On passing a known amount of current for a known time through a conducting coil of wire of

known resistance, immersed in a calorimeter containing water, whose initial temperature and thermal capacity were known, he noted the rise in temperature. From the observations, he calculated heat produced. From the results obtained through a series of experiments by varying the different parameters, he concluded that.

- a) Amount of heat produced gets doubled on doubling the resistance of coil, keeping current and time constant.
- b) Amount of heat produced becomes four times, if current is doubled, provided resistance of coil and time of flow of current remained constant.
- c) If time of flow of current through the coil is doubled, the amount of heat generated also becomes twice for same coil resistance and current.

Thus, if "H" is the heat produced on account of flow of current "I" through a conductor of resistance "R" for a time period "t", then mathematically

$$H = I^2 R t$$

Same result can also be derived from Ohm's law. Joule was not aware of Ohm's law, when he experimentally deduced his law for production of heat due to flow of electric current.

By definition, the potential difference between two points of a conductor is one volt if one joule of work is done in carrying a unit charge (of a coulomb) from one point to the other. If "I" is the magnitude of current that flows through a conductor of time "t", the quantity of charge "Q" that has flown through the conductor is given as

$$Q = It$$

The work done in carrying charge “ Q ” against a potential difference of one volt is = It joules.

Therefore, work done against a potential difference of “v” volts (in Joules) is

$$W = VIt$$

.....(1)

Equation (1) gives the energy spent in overcoming the resistance offered by the conductor. This complete energy is converted into heat.

If “J” is the mechanical equivalent of heat, then by definition,

$$\text{Heat produced} = H = \frac{W}{J} \text{Calories} = \frac{VIt}{4.186} \text{Calories}$$

Here, “J” is the Joule’s constant = 4186 J/Cal

Also, from Ohm’s law, $V = IR$

$$\therefore H = \frac{I^2 R t}{4.186} \text{Calories}$$

$$\Rightarrow H \propto I^2 R t$$

For better results, it is more desirable to use equation (2) than equation (3), since, resistance of conductor “R” in equation (3) varies significantly with the change in temperature.

Observations and calculations

1. Water equivalent of calorimeter =W=
2. Mass of empty calorimeter (in grams)=M=
3. Mass of Calorimeter filled with water (in grams) =M’=
4. Mass of calorimeter filled inside calorimeter (in grams)=m=M’ -M=
5. Ammeter reading (in Amperes)= I =
6. Voltmeter reading (in Volts) = V=
7. Initial temperature of water (in °C) =1=
8. Final temperature of water (in °C)=2=
9. Time for which current is passed (in seconds)=t=
10. Decreased temperature when it is allowed to fall for the same time duration “t” (in °C) =3=
11. Corrected final temperature after applying radiation correction (in °C) =4=

$$= \theta_2 + \frac{\sigma_3}{0} t$$
12. Rise of temperature = =4-1=
13. Mechanical equivalent of heat = $\frac{VIt}{(m + W)\theta}$?

Precautions

1. The turns of heating coil should not touch each other.
2. Constantly stir the water with stirrer when electric current is flowing through the coil.
3. In order to obtain better results, some other liquid of known specific heat such as oil can be used. In case of water, certain amount of electrolysis takes place which is quite insignificant when the voltage applied does not exceed 8-10 volts and resistance of coil is less than 2-3 ohms. Other advantages of using oil are (i) losses due to evaporation are small due to their low vapor pressure; and (ii) rise in temperature for a given amount of heat is more on account of low specific heat.

Description of apparatus

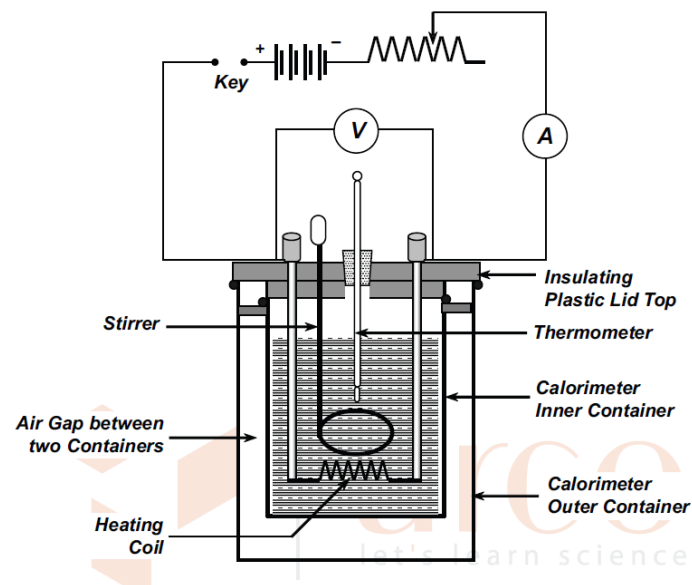
Diagram 1: Experimental setup with Joule's Calorimeter

Joule's apparatus is used for determination of Mechanical equivalent of Heat "J" (also called Joule's constant) by electrical method using the principle of calorimetry. The apparatus consists of two cylindrical shaped metal containers, having sea miles construction and rolled rims, both of different diameter and size. With the smaller container located inside larger one, the air gap between the walls of smaller and larger container provide sufficient insulation, which significantly reduces the heat loss. A high resistance coil is mounted to the underside of plastic top lid of containers through two binding posts, which ends in two color-coded 4mm socket terminals on upper side of lid and are used for connecting electrical

power source to the coil. Plastic lid also has holes for stirrer and mounting of thermometer.

If “M” is the mass of calorimeter (with stirrer) , W is the water equivalent of calorimeter “m” is the mass of water contained in calorimeter and, 1 and 2 are initial and final temperatures of water then heat taken by calorimeter and its contents “H” is given as

$$H = (m + U)(\theta_2 - \theta_1) \quad \dots\dots\dots(4)$$

Hence, from equation (2)

$$J = \frac{VI t}{H} = \frac{VI t}{(m + U)(\theta_2 - \theta_1)} \quad \dots\dots\dots(5)$$

If “S” is the specific heat of calorimeter, then MSW=

For determination of water equivalent or specific heat of calorimeter, refer to operating instructions of calorimeter.

Determination of mechanical equivalent of heat (J)

Procedure

1. Weigh the complete calorimeter (empty) along with stirrer using a suitable weighing balance. This will give the mass of complete calorimeter (M).
2. Fill the inner container of calorimeter to about 2/3rd level with water and ensure that on closing the top lid coil is completely immersed in water.

3. Weigh the complete calorimeter again. This will give the mass of complete calorimeter filled with water (M+m).
4. Complete the experimental setup as shown in diagram 1.
5. Insert the thermometer inside the smaller inner container of calorimeter through a hole in the top lid using rubber stopper with a hole. See to it that bulb of thermometer is well immersed in water inside the container.
6. Start the flow of current by inserting plug in the key and adjust the rheostat so as to obtain a current of 1-2 A with a potential difference of about 6V. Note the initial temperature of water (1) , ammeter and voltmeter readings. Start the stopwatch.
7. The temperature of water inside calorimeter will start rising. Continuously stir the water with the stirrer. Adjust any variation in ammeter and voltmeter reading to keep their value steady using rheostat.
8. When the temperature of water has risen by about 10°C, stop the stopwatch and remove the plug from key to stop flow of electric current. Note the final temperature of water (2). Also note the stopwatch reading to obtain total time of flow of electric current through the coil.
9. Allow the temperature of water to fall and note the decreased temperature (3) during the same time duration as that noted in step 9, above. This is used to correct the value of 2 for radiation error.