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Instructions on environment protection

- At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances.
- It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.
- The materials are recyclable as mentioned in its marking.
- By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.
- Please inquire at the community administration for the authorized disposal location.

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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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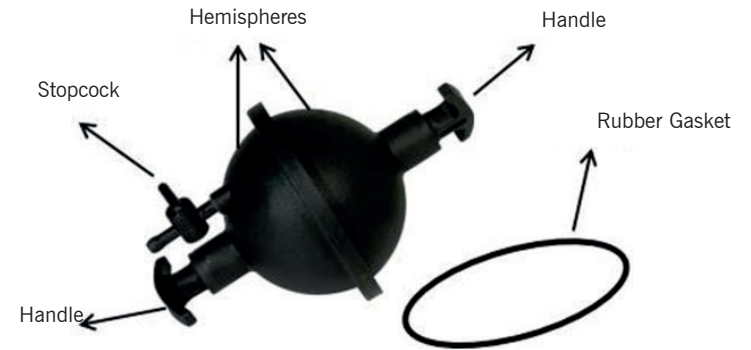
4. Procedure

- (1) Place the rubber gasket between the two hemispheres.
- (2) Connect the vacuum pump with the nozzle of stopcock via rubber tubing provided and turn the knob of the stopcock toward the nozzle direction.
- (3) Now start to evacuate the apparatus using the vacuum pump.
- (4) After evacuating the apparatus for one-minute stop further evacuating it and turn the knob of the stopcock at perpendicular to the nozzle direction.
- (5) Now pull the hemispheres; if these are easily removed then start evacuating again.
- (6) When the apparatus is proper evacuated then it will not separate out even on applying a big force on hemispheres.

5. Trouble shooting

Please check the following steps on facing the problem in evacuating the apparatus.

- (1) It may ensure that the rubber gasket is proper fitted inside the apparatus otherwise remove it and place it again inside the apparatus.
- (2) Check that the nozzle is proper fitted with the apparatus, there should not any air gap between nozzle and hemisphere.

INDEX OF LANGUAGES

Spanish	2-3
English	4-5

INDEX OF CONTENTS

1. Purpose	4
2. Other items required	4
3. Description	4
4. Procedure	5
5. Trouble shooting	5

1. Purpose

It has been designed to demonstrate the tremendous pressure of atmosphere on an evacuated object.
Apparatus consists of:

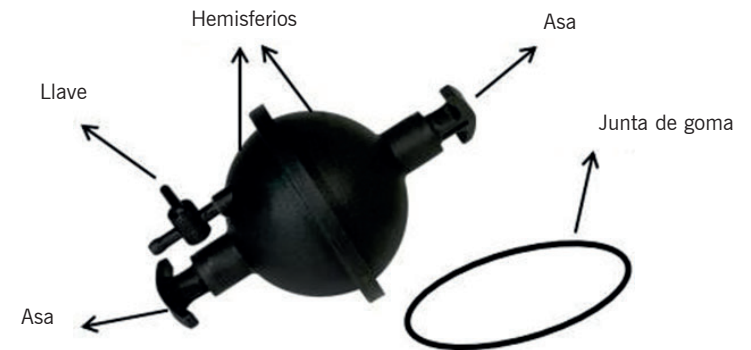
- (1) A set of Magdeburg Hemispheres
- (2) A rubber gasket

2. Other items required

- (1) Hand operated / motor driven vacuum pump

3. Description

Two plastic moulded hemispheres of about 100 mm diameter, both having especially designed handles for better gripping at their top. One of the two hemispheres has a plastic stopcock for pumping out air for generating vacuum inside the hemispheres. A rubber gasket to maintain the vacuum is provided with the apparatus. The lightweight plastic construction makes them easy to handle and convenient to use. It can be easily connected to any suitable vacuum pump for evacuation.

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