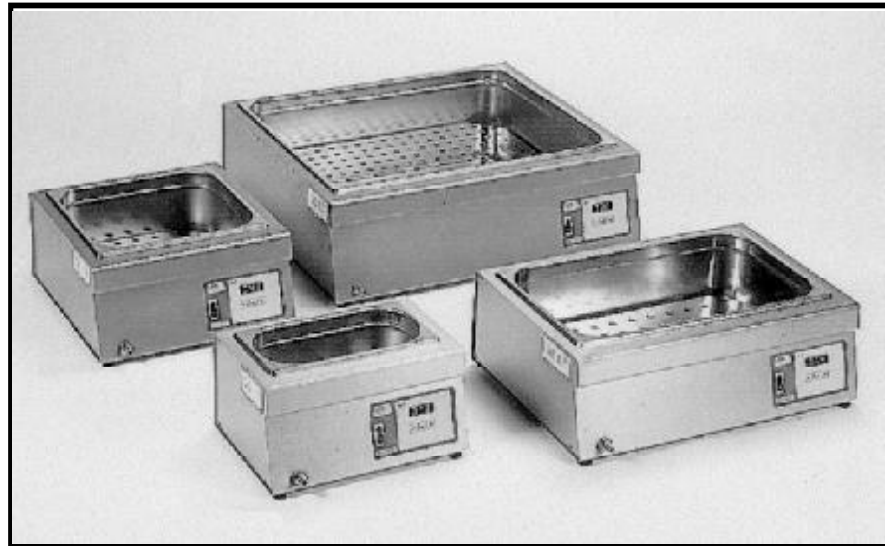
AENOR
ER
Empresa
Registrada
ER-109/1/96

PRECISDIG 6001195
PRECISDIG 6001196
PRECISDIG 6001197
PRECISDIG 6001198

PRECISDIG 6001236
PRECISDIG 6001237
PRECISDIG 6001238
PRECISDIG 6001239

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**WATER AND OIL BATHS WITH DIGITAL REGULATION
AND READING OF TEMPERATURE AND TIME**

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

SOMMAIRE

INFORMATION GENERALE
SOMMAIRE
LISTE D'EMBALLAGE
ACCESSOIRES
SPECIFICATIONS TECHNIQUES
DESCRIPTION DE L'APPAREIL
INSTALLATION
MODE OPERATOIRE
SCHEMA ELECTRIQUE
PIECES DE RECHANGE
GARANTIE
ENTRETIEN
DECLARATION DE CONFORMITE "CE"

CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
ACCESSORIES
TECHNICAL FEATURES
EQUIPMENT DESCRIPTION
INSTALLATION
OPERATION
ELECTRICAL DIAGRAM
SPARE PARTS
GUARANTEE
MAINTENANCE
"EC" CONFORMITY DECLARATION

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The standard equipment consist of the following components:

Pièce / Piece	Code/ Code	Qté / Qty.
Precisdig	6001195	1
Precisdig	6001196	1
Precisdig	6001197	1
Precisdig	6001198	1
Precisdig	6001236	1
Precisdig	6001237	1
Precisdig	6001238	1
Precisdig	6001239	1
Guide de l'utilisateur Instruction manual	80037	1

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RACKS ADAPTABLE TO THE SUPPORT-FRAME

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SPECIFICATIONS TECHNIQUES**TECHNICAL FEATURES**

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Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Code / Code	6001195	6001196	6001197	6001198	6001236	6001237	6001238	6001239	
Température maxi. °C / Maximum temperature °C	100				200				
Capacité litres / Capacity litres	5	12	20	45	5	12	20	45	
Dimens. utiles cm. Tank dimensio- ns cm	Haut. Depth	15							
	Larg. Length	30	30	48	62	30	30	48	62
	Prof. Width	14	30	30	50	14	30	30	50
Dim. extér. cm Overall dimensio- ns cm	Haut. Depth	20							
	Larg. Length	37	40	58	69	37	40	58	69
	Prof. Width	27	42	42	62	27	42	42	62
Consom. W. / Consump. W.	600	1100	1500	2400	1100	1400	2100	3800	
Poids Kg. / Weight Kg.	6	8	10	18	6	8	10	18	

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- * Remember that these are static baths.
- * Depending on the kind of fluid, the homogeneity can be altered.

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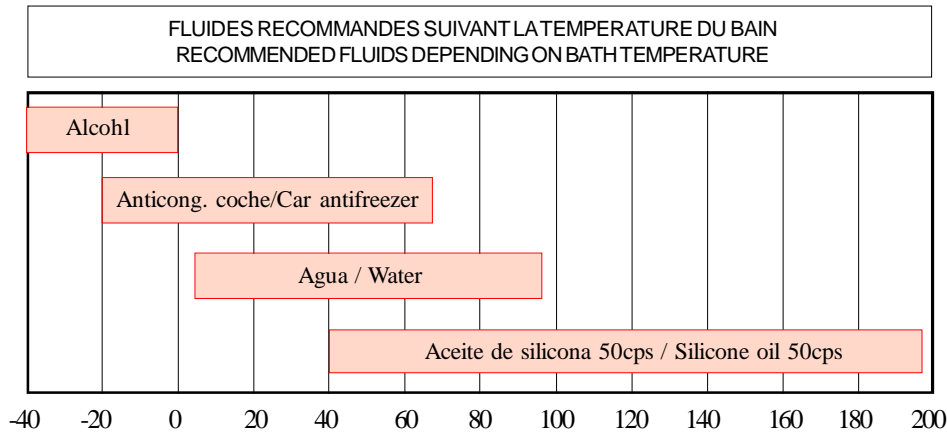
EQUIPMENT DESCRIPTION

PRECISDIG baths with adjustable temperature from ambient +5°C up to 100°C or up to 200°C, depending on the model. To use the baths fill the tank with liquid until the level covers at least the heater element. The liquid or mixture to be employed must be adequated to the model of bath and the working temperature.

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As a guidance, here are some liquids (to use these liquids, follow the manufacturer's instructions):

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INSTALLATION

Place the bath on a flat, horizontal and level surface. Bear in mind the weight of the liquid when choosing the support. Provide for bath drainage.

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Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the apparatus.

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Do not use the apparatus if it is NOT earthed.

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If you have to change the plug bear in mind the following:

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Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.

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1. Main switch with indicator lamp.
2. Temperature mode indicator.
3. Time mode indicator.
4. Temperature/Time display.

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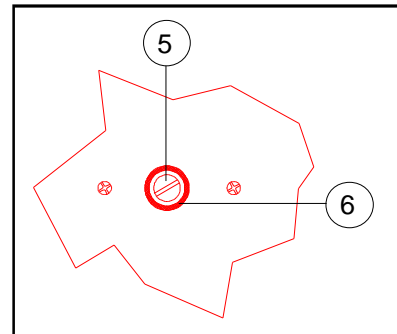
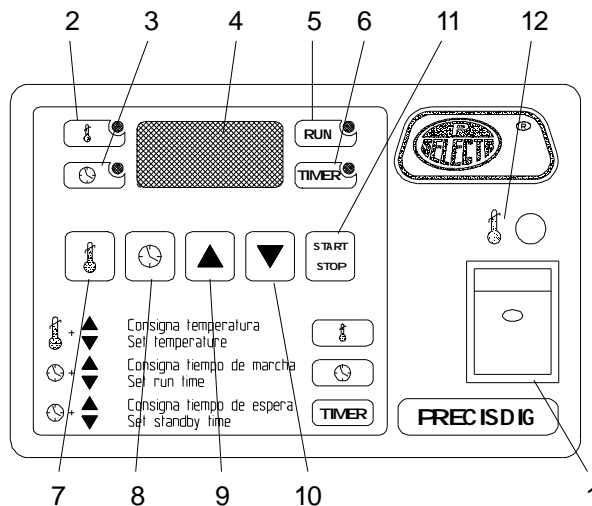
5. Starting up indicator lamp.
6. Waiting time indicator lamp.
7. Temperature selector push-button (Set of temperature from ambient +5°C up to 100°C/200°C depending on model).
8. Time selector push-button.
 - Set of starting up time: Working period between 1' and 9h 59', or up to 99.9h. once the set temperature has been reached.
 - Set of waiting time: Period of time until the starting up, between 1 and 24 hours. Decimal fraction of hour (1.5 = 1hour 30 minutes).
9. Parameter increase push-button.
10. Parameter decrease push-button.
11. Start/Stop push-button.
12. Safety thermostat indicator lamp.

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REAR PART

5. Safety thermostat regulation knob.
6. Safety thermostat reset knob (grey).

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When pressing the main switch (1), placed in the lower part of the frontal, the display lights up (4).
 Pressing the push-button (7) and the (9) or (10) we preset the desired temperature.
 During this operation the indicator (2) lights up.
 Pressing the push-button (11) the bath will start and the indicator (5) will light up.
 To stop, press the push-button again (11) and the indicator will be turned off (5).

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2. TEMPORIZED MODE AT THE DISCONNECTION:

When pressing the main switch (1) the display lights up (4).

Pressing simultaneously the push-button (7) and the (9) or (10) we preset the desired temperature.

During this operation the indicator (2) lights up.

Pressing simultaneously the push-button (8) and the (9) or (10) we preset the disconnection time.

During this operation the indicator (3) lights up.

The disconnection time begins when the bath has arrived to the preset temperature.

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In this operation mode, the indicator (5) remains intermittent until the bath reaches to the preset temperature, once reached, the indicator (5) remains lighted up.

When the preset time has lapsed, the bath stops automatically and the indicator (5) will go off.

You can stop it manually pressing the push-button (11) again.

3. TEMPORIZED MODE AT THE STARTING UP:

When pressing the main switch (1) the display lights up (4).

Pressing simultaneously the push-button (7) and the (9) or (10) we preset the desired temperature.

During this operation the indicator (2) lights up.

Pressing simultaneously the push-button (8) twice and the (9) or (10) we preset the waiting time until the starting up. During this operation the indicator (6) will light up.

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In this operation mode, the indicator (6) remains lit up during the waiting time, this over, the indicator lamp (6) goes off and the bath starts automatically.

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4. TEMPORIZED WAY AT THE STARTING UP AND AT THE DISCONNECTION:

When pressing the main switch (1) the display lights up (4).

Simultaneously pressing the push-button (7) and the (9) or (10) we preset the desired temperature.

During this operation the indicator (2) lights up.

Simultaneously pressing the push-button (8) twice and the (9) or (10) we preset the waiting time until the starting up. During this operation the indicator lamp (6) will light up.

Simultaneously pressing the push-button (8) and the (9) or (10) we preset the disconnection time.

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During this operation the indicator (3) lights up. The disconnection time begins when the bath has reached to the preset temperature.

Pressing the push-button (11) the bath will start and the indicator (5) will light up.

In this operation mode, the indicator lamp (5) remains intermittent until the bath reaches to the preset temperature, once reached, the indicator (5) remains lighted up.

Lapsed the preset time, the bath stops automatically and the indicator (5) will go off.

You can stop it manually by pressing the push-button (11) again.

Pressing the push-button (11) the bath will start and the indicators (5) and (6) will light up.

In this operation mode, the indicator lamp (6) remains lit up during the waiting time, once over, the indicator lamp (6) goes off and the bath starts automatically.

ADJUSTMENT FOR THE SAFETY THERMOSTAT:

The low limit of working temperature for the safety thermostat is nearly 50°C. The desired temperature of the safety thermostat must be at least 5°C above the working temperature. The adjusting process of the safety thermostat is long. For that reason, the safety thermostat is useful when the users work habitually at the same temperature.

1. Turn the safety thermostat screw (5) clockwise to maximum.
2. Select with the temperature selector push-button (7) the desired temperature of the safety thermostat.
3. Turn on the bath and leave it to stabilize.
4. When the bath is stabilized, slowly turn the safety thermostat screw (5) anticlockwise until it clicks and the red indicator lamp (15) lights up.
5. Wait until the bath cools to reset the safety thermostat with (6) button.
6. Select the working temperature of the bath with temperature selector push-button (7).

MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

The control elements and heating elements must only be manipulated by authorized personnel.

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The distilled or demineralized water has a great tendency to rust. To prolong the life of the bath and avoid rusting of the parts, dissolve 1 gr of CO_3Na_2 (Carbonate of Soda) per litre of water used in the bath.

CLEANING:

If normal tapwater is used calcareous deposits will appear in the bath and in the heating element. A heater covered with lime should not be used as it can lose heating power and deteriorate more quickly. Periodical cleaning of the inside of the tank and the heating element with a solution of 10% of 28% Chlorhydric acid or a solution of water and vinegar (Acetic acid) is advised. Rinse with clean water afterwards.

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: alcohol

Cleaning of plastic: Alcohol with cotton duster.

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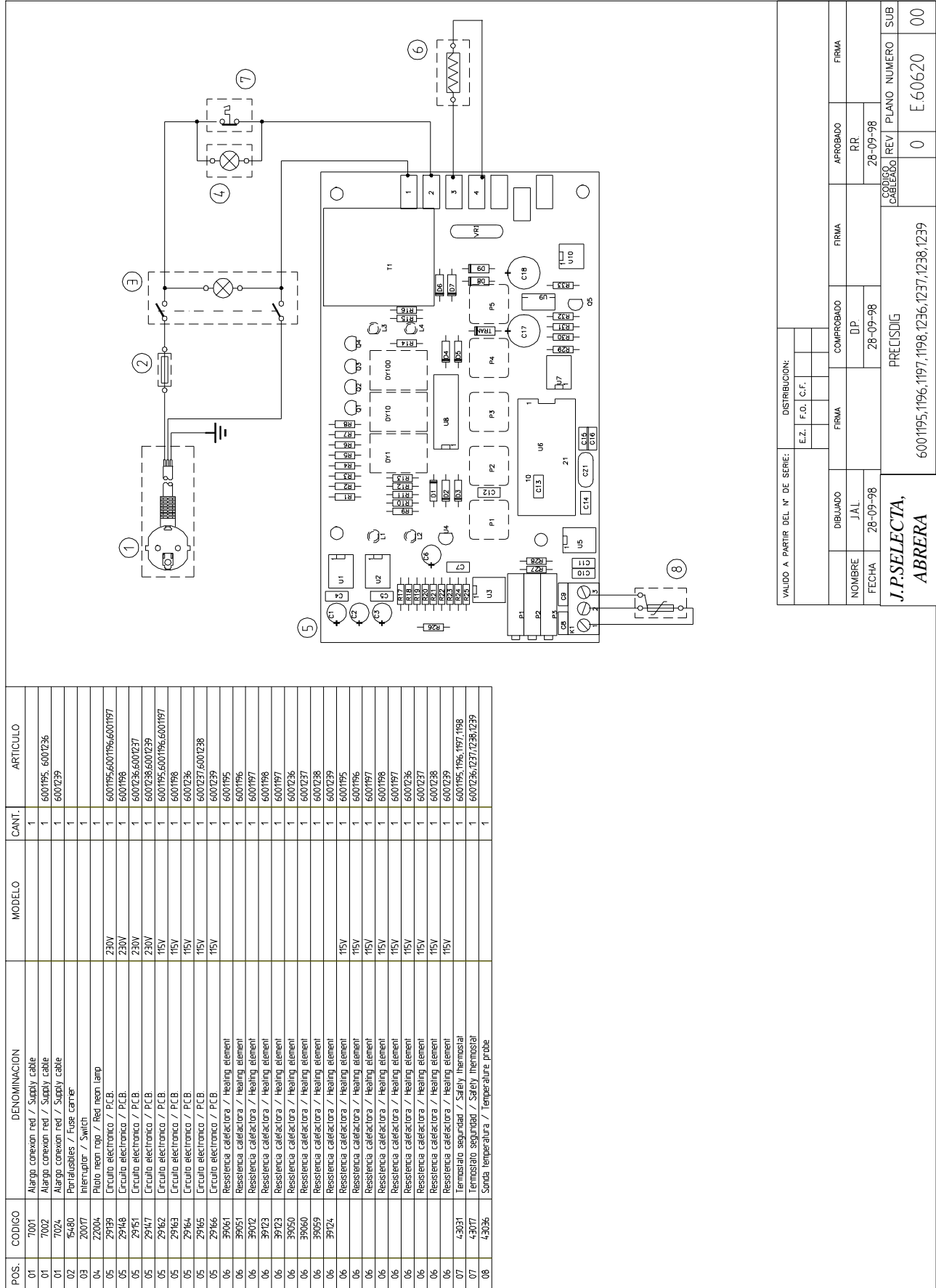
GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA, S.A..

Any manipulation of the apparatus by unathorized personnel by J.P. SELECTA, S.A. cancels the guarantee automatically.

SCHEMA ELECTRIQUE / ELECTRICAL DIAGRAM

DECLARATION DE CONFORMITE "CE" "EC" CONFORMITY DECLARATION

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PROGRAMME DE FABRICATION / MANUFACTURING PROGRAMME

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