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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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TABLE OF CONTENTS

1. Technical specifications	4
2. Operation	5
3. Incidents	5

1. TECHNICAL SPECIFICATIONS

Equipment designed to provide a hermetic and secure seal on thermoplastic bags and materials. Its robust structure and advanced functionality ensure a perfect seal, protecting the contents from any leakage or contamination. Equipped with an integrated roll holder, it facilitates material handling, optimising the sealing process and guaranteeing consistent results with every use.

Characteristics:

- Made of ABS combined with metal and with 4 rubber feet on the base.
- Easy to operate, with pressure sealing and manual cut-off
- Start-up indicator light

Code	ZZC003
Sealing width	15 mm
Max. sealing length	300 mm
Max. plastic thickness	0.4 mm
Dimensions	370x310x180 mm
Consumption	300 W
Weight	4.50 kg

2. OPERATION

- Before operating, check the different parts of the equipment.
- Plug the cable into the mains socket.
- Switch on the on/off switch.
- The green indicator light comes on. The equipment has a warm-up process, when the green indicator light turns off, the equipment is ready for use.
- Pass the paper through the tape and press the pressure clamp down.
- The equipment does not have an operating time setting. It can be adjusted by the user as required.
- After sealing is complete, press the paper cutter and slide it sideways to cut the paper.

Note: The housing is specifically designed to prevent overheating. However, to avoid burns, when using the equipment for a long period of time, pay attention to the temperature of the housing.

3. INCIDENTS

Note: If you experience any problems while using the equipment, please contact your supplier's Technical Service with a detailed description of the problem, the actions taken, and the results obtained.

- The equipment does not heat up.
Check if the switch or connector is damaged. If they are damaged or the problem persists, contact Technical Service.
- The green light does not go out for a long time.
Check if the switch is damaged. If it is damaged or the problem persists, contact Technical Service.
- Other rulings
Check that the equipment is connected under proper electrical conditions (AC 220V/60Hz). Other incidents or errors should be evaluated by a professional for repair.