


# NAHITA PARAFFIN DISPENSER

## Code ZFP013

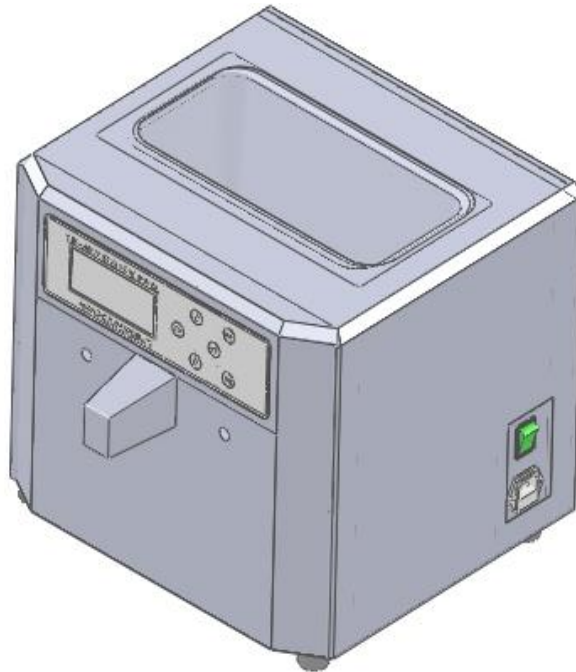

## Operation Manual

Please read this manual before using the dispenser and keep it for future consultations

# Contents

1. Use of the equipment/safety
2. Technical parameters
3. Installation
4. Operation

## Use of the equipment/safety

This equipment is used to melt and dispense paraffin.

Before heating, pay attention to the preset temperature, it must be within the melting range of the paraffin (60-80°C).

Do not press the paraffin drain switch until the paraffin is melted. In this way you avoid damaging the tap.

During heating do not open the lid of the bath.

Avoid the entry of impurities in the bath that can clog the spout.

## Technical parameters

1. Rated voltage: 220VAC, 50Hz
2. Heating temperature range: RT - 99°C
3. Programmable maximum time: 24 hours
4. Reservoir dimensions: 300 mm × 155 mm x 150 mm
5. Reservoir capacity: 5 L

## **Installation**

### *Initial inspection*

- Unpack the paraffin dispenser, remove the plastic that surrounds it and remove the poliespan protection.
- Without connecting the equipment to the power supply, make sure that it does not present any damage due to transport. If so, notify your distributor immediately so you can make the relevant claims within the deadline set by the transportation service.

### *Installation*

- Place the paraffin dispenser on a horizontal and stable table, leaving the necessary work space around it.
- Comply with all the measures required by current legislation on occupational safety and health (electrical safety devices, flammable, toxic and /or pathological products, etc.).
- Do not place the equipment in areas near heat sources, magnetic material or expose it directly to sunlight.
- The presence of flammable products (gases and liquids) in the workplace is prohibited.
- Insert the power cable into the 220V 50Hz socket provided with a ground connection and the other end to the connector of the paraffin dispenser.
- When you are not going to use the equipment for long periods of time, make sure to disconnect it from the power.

## Operation

- Turn on the equipment by pressing the “ON-OFF” switch.
- Press the OK key. The menu screen will appear with the options TIME, START, TEMPERATURE and AUTOMATIC. You can select one of them using the arrows keys and then pressing OK.
- By pressing the RESET key you can return to the menu screen.
- The paraffin dispenser has two modes of operation: normal and automatic.
- Normal operation: Select TEMPER. Press the key →. Adjust the heating temperature value by pressing ↑ or ↓. Then press OK. Set the temperatures for bath and spout.
- Select START to begin heating the bath. The set temperature (Setup) and the bath temperature will appear on the screen.
- To end the bath heating press RESET. Turn off the equipment by pressing the “ON-OFF” switch.
- Automatic operation: Select TIME. This screen shows SYS TIME, START TIME and EXIT TIME.
- Press → to adjust SYSTEM TIME. Press ↑ or ↓ to adjust the hours. Press →. Press ↑ or ↓ to adjust the minutes. Press OK.
- Select TIME again. Set the START TIME and the EXIT TIME. Press OK.
- Select AUTOMATIC. The display shows WAITING, the system time and the start time for heating.
- When the start time comes, the screen shows the set temperature (Setup) and the paraffin temperature in the reservoir.

- When the exit time comes, the screen shows again WAITING, the system time and start time for heating.
- If you are not going to use the equipment again, turn it off by pressing the “ON-OFF” switch.