MANUAL DE USO
USER'S MANUAL

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Modelo: _____

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Thank you for choosing this equipment. We sincerely wish that you enjoy your electrophoresis power supply Nahita. We highly recommend that you look after the equipment according to what is stated in this manual.

Nahita develops its products following the CE guidelines taking into account the ergonomics and safety of the user. The quality of the manufacturing materials and the correct use will allow you to enjoy this equipment for many years.

The misuse of this equipment may cause accidents, electric shocks, shorts, fires, injuries, etc. Check "Safety" to read safety guidelines.

Safety

Caution/Warning

Electrophoresis power supplies use high output voltages that are electrically isolated from earth ground through a protective impedance to minimize the risk of electrical shock to the user. The following guidelines should be observed and followed when using a power supply.

Electrophoresis power supplies have passed tests for operation at temperatures between 0° and 40°C, with non-condensing relative humidity between 0 and 95%. Operating the power supply outside these conditions is not recommended by Nahita and will void the warranty.

1. To ensure adequate cooling of the power supply, be sure that there is at least 6 cm clearance around the power supply. Do not block the fan vents at the rear of the unit.
2. Always connect the power supply to a 3-prong, grounded AC outlet, using the 3-prong AC power cord provided with the power supply.
3. Do not operate the power supply in extreme humidity (>95%) or where condensation can short the internal electrical circuits of the power supply.
4. When taking the power supply into a cold room, the unit can be operated immediately. However, when removing the power supply from the cold room, let the unit equilibrate to room temperature for a minimum of 2 hours before using it.
5. Never connect a high voltage output lead to earth ground. This defeats the floating electrical isolation of the power supply and exposes the user to potentially lethal high voltages.

Important

This instrument is intended for laboratory use only. This product conforms to the class A standards for Electromagnetic Emissions, intended for laboratory equipment applications. It is possible that

emissions from this product may interfere with some sensitive appliances when placed nearby or on the same circuit as those appliances. The user should be aware of this potential and take appropriate measures to avoid interference.

CONTENTS

1. Introduction.....	13
2. Setup and operation.....	13
3. Error alarms.....	15
4. Warranty.....	15

1 Introduction

The power supply ZFD016 has the following technical specifications and characteristic features:

Technical specifications

- Type of Output: Constant-Voltage, Current or Power
- Increment: 1V, 1mA, 1W
- Timer Range: 1min~99hr, 59min
- Display: LCD display with backlight
- Output jacks: Four sets of output jacks

Characteristic features

- Molding shell, touch keys, dual core microprocessor intelligent control;
- Indicates the preset value and the actual output value at the same time;
- It can store 10 electrophoresis methods;
- With automatic memory function;
- With standard, timing operation function;
- With constant voltage, constant current, constant power, misoperation, fault intelligent prompt functions;
- Automatic detection of no-load, over-load, short circuit, rapid resistance change, ground leak and system overheating.

2 Setup and Operation

2.1 Turn on the power switch, enter into Setting State.

EDIT[0]	
U= 100V	0v
I= 500mA	0mA
P= 300W	0w

- 2.2 The cursor will flash at “U=□□□V”. Set the value of constant voltage by pressing “▲” or “▼” button. The voltage range can be set from 10 to 600V.
- 2.3 The cursor will flash at “I = □□□mA” when you press “ENT” button. Set the value of constant current by pressing “▲” or “▼” button. The current range can be set from 1 to 500mA.
- 2.4 The cursor will flash at “P = □□□W” when you press “ENT” button. Set the value of constant power by pressing “▲” or “▼” button. The power range can be set from 1 to 300W.
- 2.5 The cursor will return and flash at “U = □□□V” when you press “ENT” button. Set the value of constant voltage again by pressing “▲” or “▼” button.
- 2.6 Power supply in Setting State now:
Press “RUN/STOP” button, it will enter into Operating State.

EDIT[0]	RUN! 00:01
U= 100V	100v
I= 500mA	350mA
P= 300W	35w

- 2.7 Power supply in Operating State now:
Press “RUN/STOP” button, it will stop output and return to Setting State.
- 2.8 In Setting State, continue to press “EDIT” button, it will enter into Editing State and display:
“SAVE [□]” → “LOAD [□]” → “T= □□:□□” → “QUIT” → “SAVE [□]”...

EDIT[0]	
U= 100V	0v
I= 500mA	0mA
P= 300W	0w

- 2.9 At “SAVE □”, press “▲” or “▼” button to select the number of memory program. The range of numbers is from 1 to 9. Then press “ENT” button to save current method into this program.

- 2.10 At "LOAD [□]", press "▲" or "▼" button to select the needed number. The range of numbers is from 1 to 9. Then press "ENT" button to load the current method from the program.
- 2.11 At "T= □□:□□", press "▲" or "▼" button to set operating time.
- 2.12 The timer range is from 1min to 99hs and 59mins. Then press "ENT" button to save it.

EDIT[0]	
U= 100V	0v
I= 500mA	0mA
P= 300W	0w

- 2.13 At "QUIT", press "ENT" button to quit Editing state, and return to Setting State.

3 Error Alarms

- 3.1 When the power supply runs with Over-load, it will stop automatically and alarm with buzzer, then display "ERROR1!". Turn off the switch immediately.
- 3.2 When the power supply runs with no-load, it will stop automatically and alarm with buzzer, then display "ERROR 2!". Turn off the switch immediately.

4 Warranty

The Power Supply is warranted for 1 year against defects in materials and workmanship. If any defects should occur during this warranty period, Nahita will replace the defective parts without charge. However, the following defects are specifically excluded:

1. Defects caused by improper operation.
2. Repairs or modifications done by anyone other than Nahita Technical Service.
3. Use with cables or connectors not specified by Nahita for this power supply.
4. Deliberate or accidental misuse.
5. Damage caused by disasters.

For inquiries or requests for repair services, contact Nahita Technical Service.

Warranty Information

Model: _____

Serial Number: _____

Date of Delivery: _____

Warranty Period: _____