APARATO LEY DE BOYLE
BOYLE'S LAW APPARATUS
REF. QJ009

INDEX OF LANGUAGES- ÍNDICE DE IDIOMAS

ESPAÑOL 2
ENGLISH..... 3

APARATO LEY DE BOYLE

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BOYLE'S LAW APPARATUS

This small and economical version gives the better and accurate quantitative results for the Pressure / Volume relationship of air at constant temperature. A graduated (0 - 300 cm³) cylinder with closely fitting piston and connected to a pressure gauge by a narrow tube to minimize the dead volume is fitted on two side supports. The cylinder is made of transparent material to see the working process inside it. We can adjust the pressure inside the cylinder by rotating the knob connected to the piston via threaded shaft. A gas tap allows the cylinder to communicate with the outside air to adjust the mass of air enclosed. The pressure can be directly read out from a big size pressure gauge circular dial, graduated in 0-7 kg/cm²

Aim: To investigate the validity of Boyle and Mariotte's law for constant amount of a gas say air.

Apparatus required: Boyle-Mariotte's Law apparatus.

Boyle-Mariotte's Law: For the limiting case of an ideal gas (sufficiently low pressures, sufficiently high temperatures), the correspondence between the state variables pressure P, volume V and temperature T is described by the ideal gas law

$$PV = nRT$$

Where R is the universal gas constant.

For cases of constant quantity of substances and isothermal process control this equation changes into the following equations:

$$PV = K$$

or

$$P = K/V$$

Procedure:

- (1) Change the pressure inside the cylinder by rotating the knob connected to the piston and note the corresponding volume.
- (2) Repeat step 1 by increasing the pressure in each step and keeping the temperature constant. Note the corresponding volume in each step.
- (3) Plot a graph between reciprocal of volume (1/V) and pressure (P), which comes out to be a straight line.

| Sr. No. | Pressure (P) | Volume(V) | 1/V |
|---------|--------------|-----------|-----|
| 1. | | | |
| 2. | | | |
| 3. | | | |
| 4. | | | |
| 5. | | | |
| 6. | | | |
| 7. | | | |
| 8. | | | |